


VÅNTEC-500 Area Detector for Pharmaceutical XRD

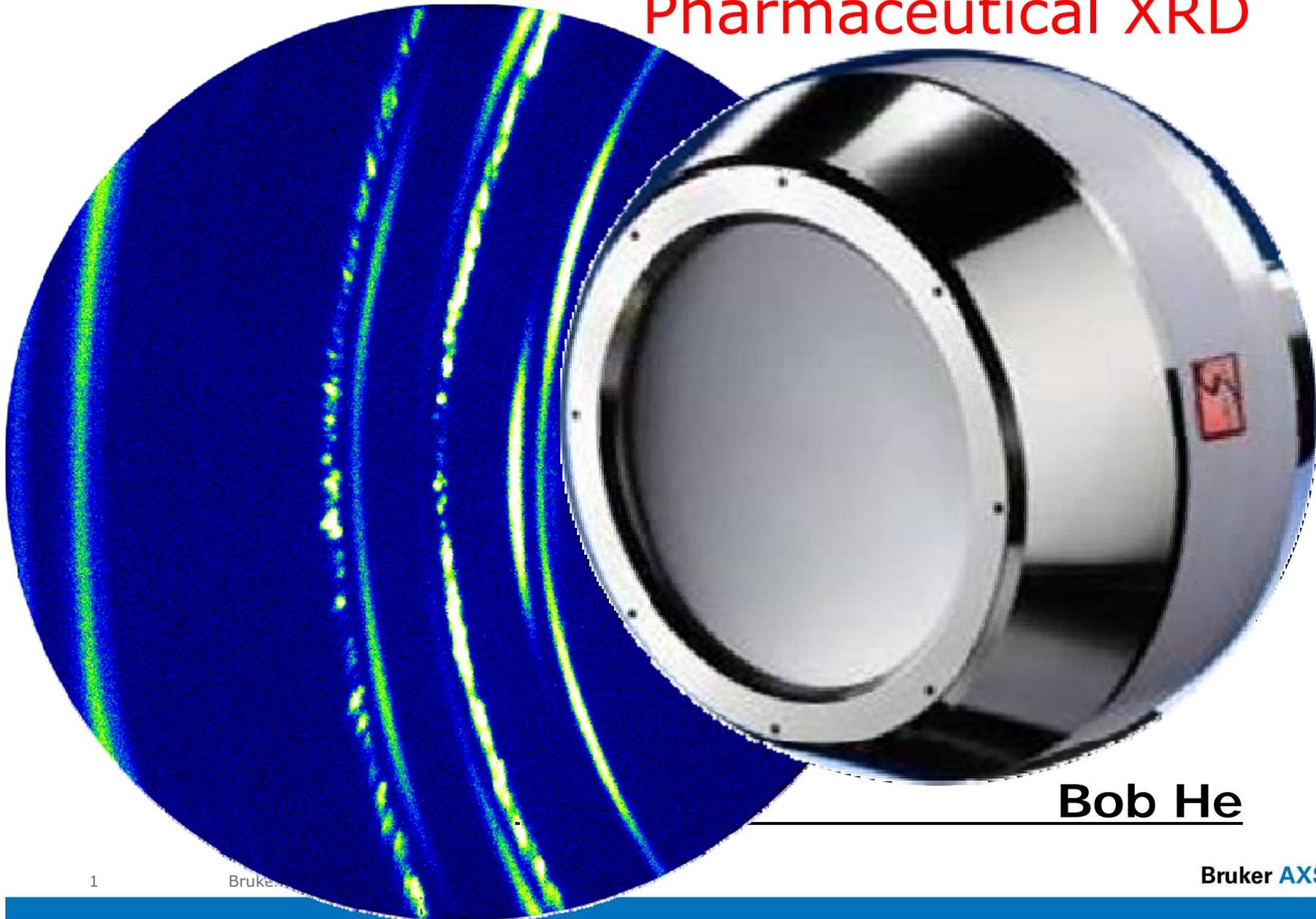

Bob He

This document was presented at PPXRD - Pharmaceutical Powder X-ray Diffraction Symposium

Sponsored by The International Centre for Diffraction Data

This presentation is provided by the International Centre for Diffraction Data in cooperation with the authors and presenters of the PPXRD symposia for the express purpose of educating the scientific community.

All copyrights for the presentation are retained by the original authors.


The ICDD has received permission from the authors to post this material on our website and make the material available for viewing. Usage is restricted for the purposes of education and scientific research.



PPXRD Website – www.icdd.com/ppxrd

ICDD Website - www.icdd.com

XRD² for Pharmaceutical: Reflection System (CS)

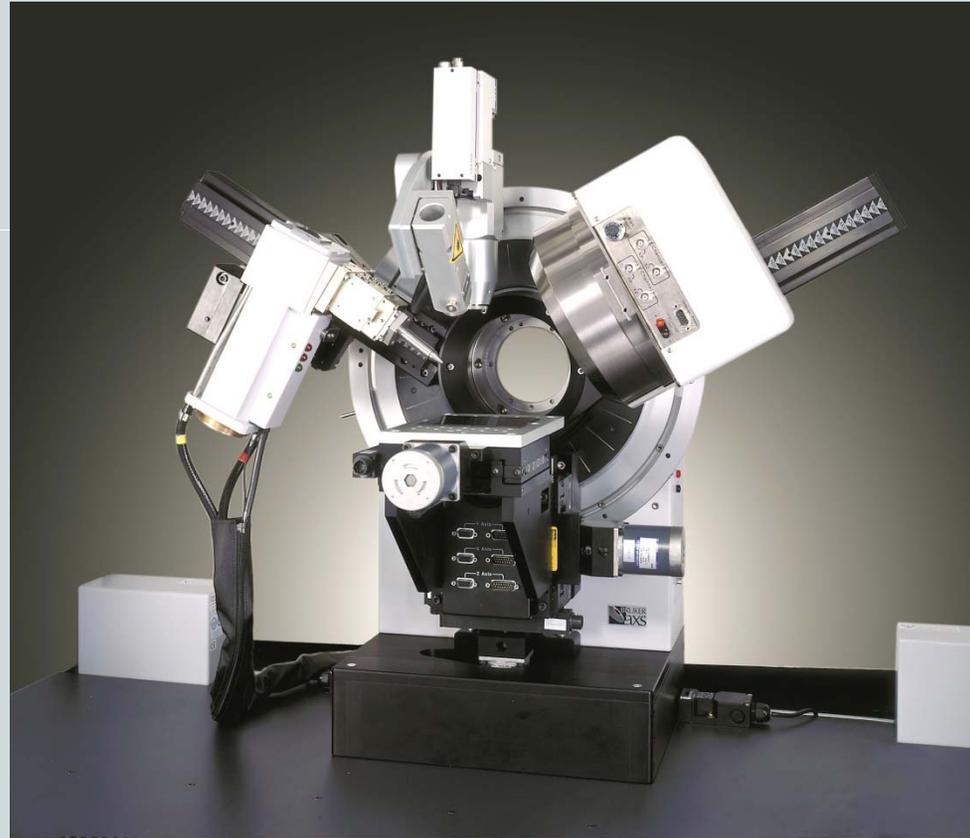

XRD² for Pharmaceutical: Reflection & Transmission (HTS)



**US Patent
#7,242,745**



XRD² : Innovation and Development

The most dramatic development in XRD² happens in three critical components and data evaluation algorithms:


- Source: required radiation energy, focal spot size and intensity.
- Optics: select wavelength, beam profile and divergence.
- Detector: collect 2D pattern with correct intensity and position.
- Data evaluation Algorithm: Diffraction Vector Approach.

X-ray Source for XRD²: Incoatec Microsource (I μ S)TM

- High brilliance
- Low energy: 30 W
- Air-cooled
- Spot size < 100 μm
- Montel mirror




$I_{\mu}S$ & VÅNTEC-2000 vs. Classic Set-up Corundum Comparison

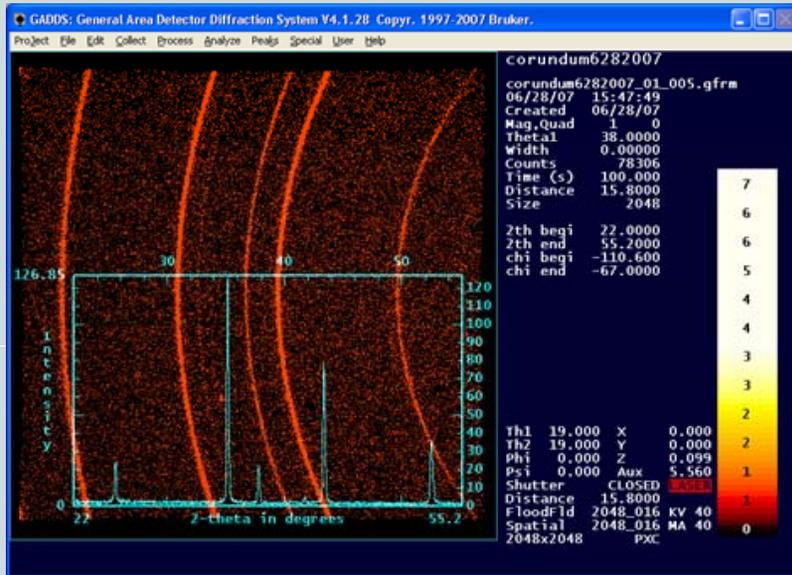


$I_{\mu}S$ & VÅNTEC-2000
45kV, 0.650mA,
0.3mm snout
total counts: 1235K

Single 40mm Göbel Mirror,
45kV, 40mA,
0.3mm collimator
total counts: 78K

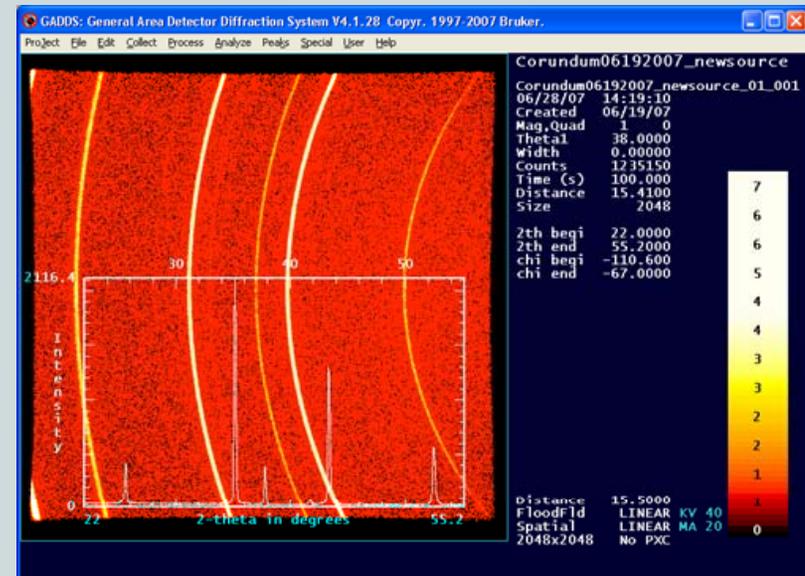

XRD²: Choice of Detectors

Sensitivity vs. Count Rate

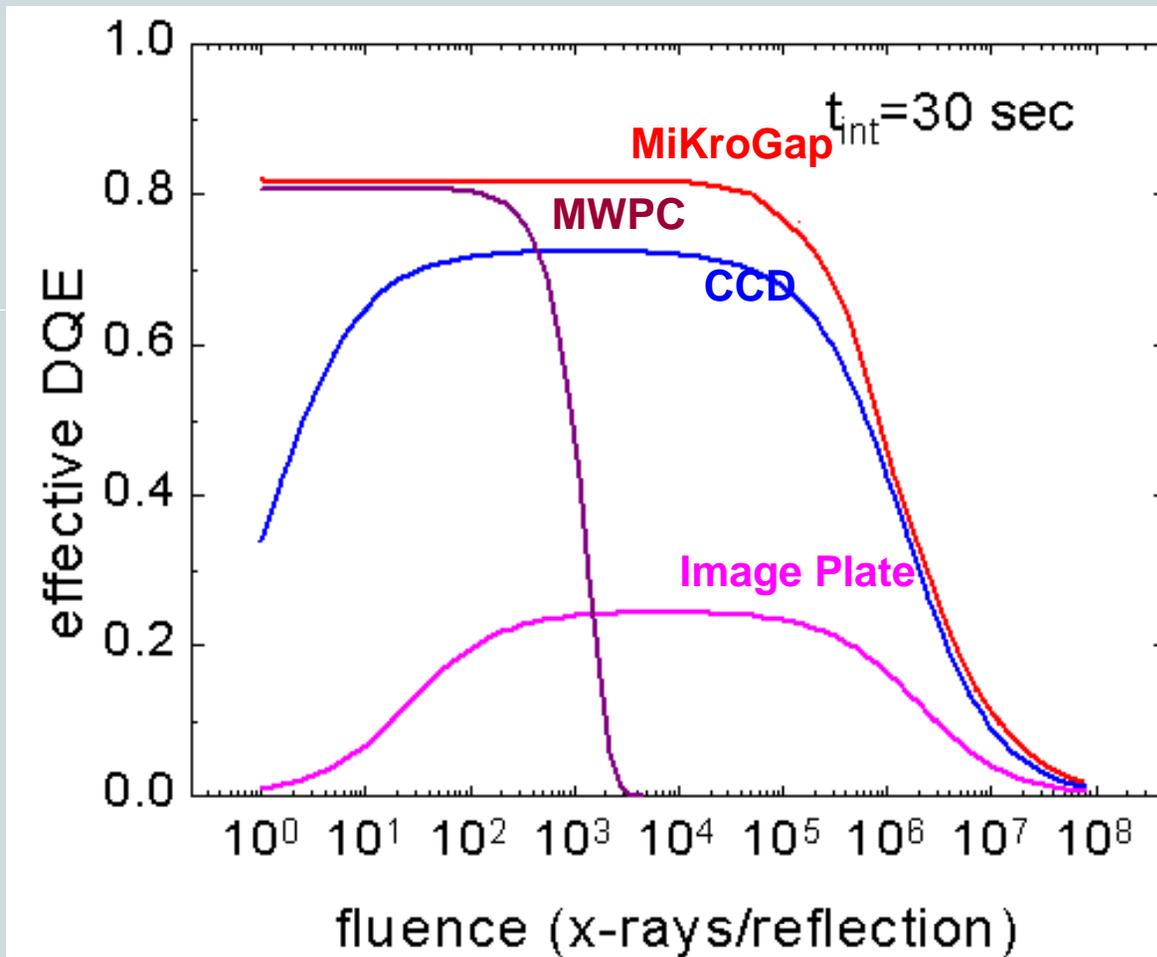

Detective Quantum Efficiency (DQE):

■ The DQE is a parameter defined as the square of the ratio of the output and input signal-to-noise ratios (SNR).

$$DQE = \left(\frac{(S/N)_{out}}{(S/N)_{in}} \right)^2$$

■ The DQE of a real detector is less than 100% because not every incident x-ray photon is detected, and because there is always some detector noise.



■ MiKroGap™ has the best overall performance.

Detector Technology from MWPC to MikroGap™



- MikroGap™ technology with resistive anode:
 - shortens drift time of ions
 - fast electrons induce charge on readout strips
- Adjusted surface resistance ($10^5 - 10^7 \Omega/\text{area}$):
 - high enough to limit discharges
 - low enough to support high count rates

US Patent US 6,340,819 B1

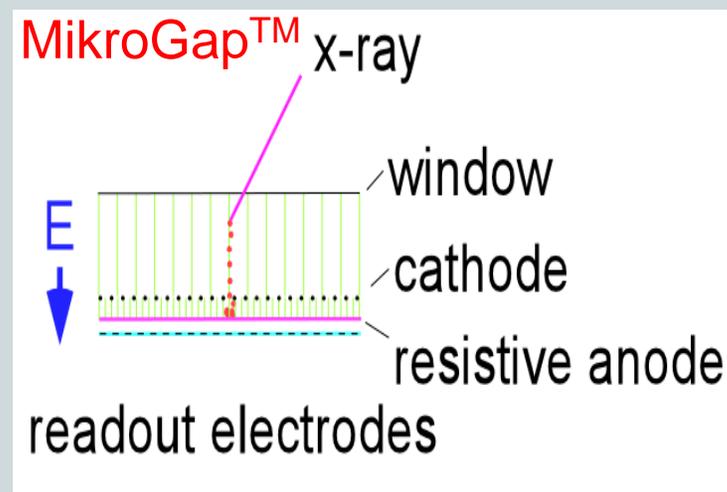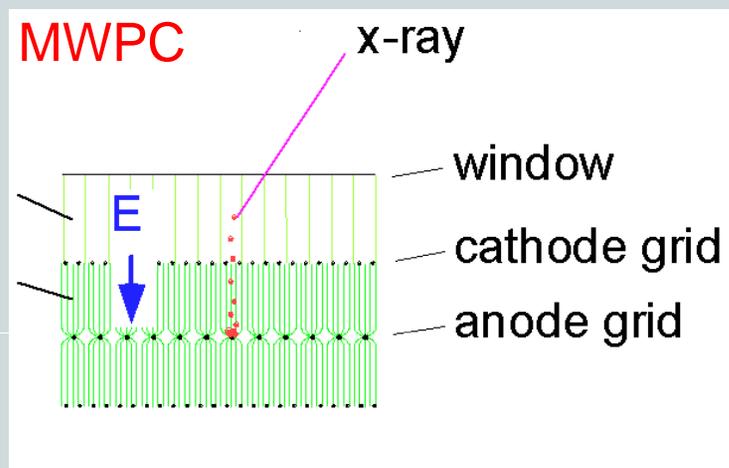

VANTEC-500 – Outperforms all previous gaseous detectors



Similar to Hi-Star (MWPC) detector:

- High sensitivity: 80% DQE (detection quantum efficiency) at 8.04 keV radiation
- Energy range: 3-15 keV (good for Cu, Co, Fe and Cr X-ray sources, not recommended for Mo)
- Energy resolution ($\Delta E/E$): 20% at 8.04 keV radiation
- Low background noise: <5 cps/global
- Readout time: real time
- No cooling
- Curved Be-window to reduce parallax

VÅNTEC-500 – Outperforms all previous gaseous detectors



Advances from MWPC:



- Tapered front end for high 2θ angle access and space for large samples and sample stages
- Doubled the spatial resolution: The FWHM of the PSF is $200\mu\text{m}$
- Two orders of magnitude higher maximum count rate:
Global count rate: 1.5Mcps
Local count rate: 250kcps per point-like reflection
- Radiation hardness: accidental intensive irradiation without permanent damage
- Maintenance-free: no re-gassing

VÅNTEC-500 – Tapered front for low and high 2θ accessibility



$$2\theta_{range} = 2 \arctan \frac{l}{D}$$

$$2\theta_{max} = \pi - \frac{m+h}{D}$$

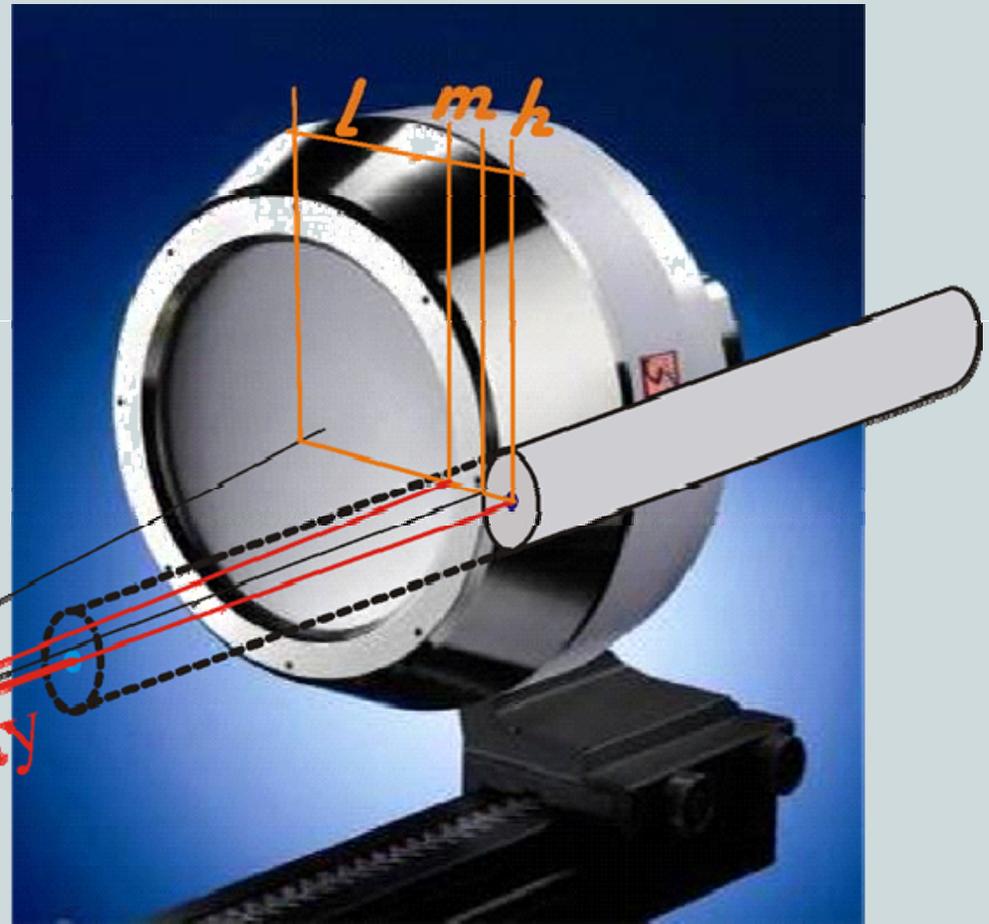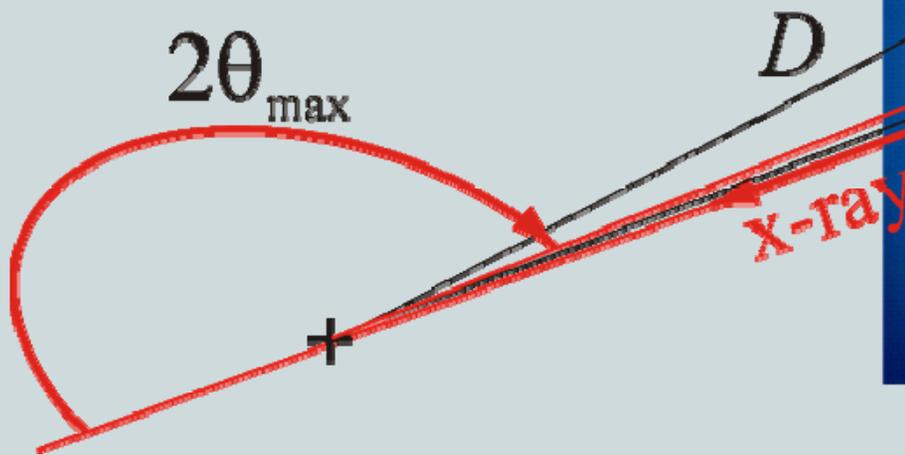

VANTEC-500 – Outperforms all previous gaseous detectors




Detector geometry:

- Be-window opening 140 mm in dia.
- Frame size:
 - 2048 x 2048 pixels
 - 1024 x 1024 pixels
 - 512 x 512 pixels
- Pixel size:
 - 68 μm x 68 μm
 - 136 μm x 136 μm
 - 272 μm x 272 μm
- Detector working distance:
5~30 cm in D8 DISCOVER enclosure
- 2θ range in a single frame:

5 cm	83°
10 cm	56°
15 cm	42°
20 cm	33°
25 cm	27°
30 cm	23°

Hi-resolution with MikroGap™ Technology

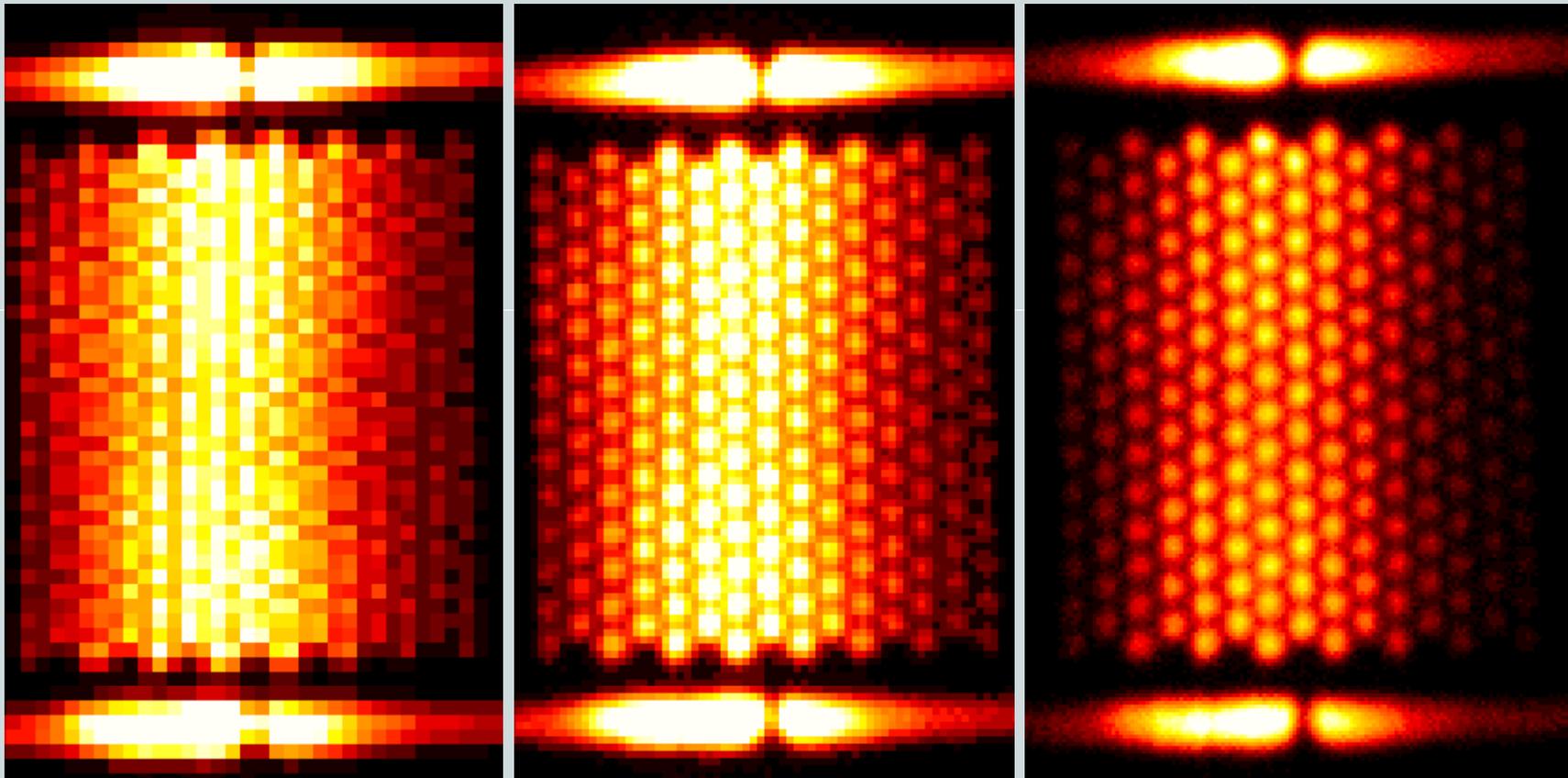

272 μm pixel (512x512) 136 μm pixel (1024x1024) 68 μm pixel (2048x2048)

- 68 μm pixel size delivers the best spatial resolution


The New D8 DISCOVER




The New D8 DISCOVER

Patented Door Mechanism

Swing door: wide opening for good accessibility

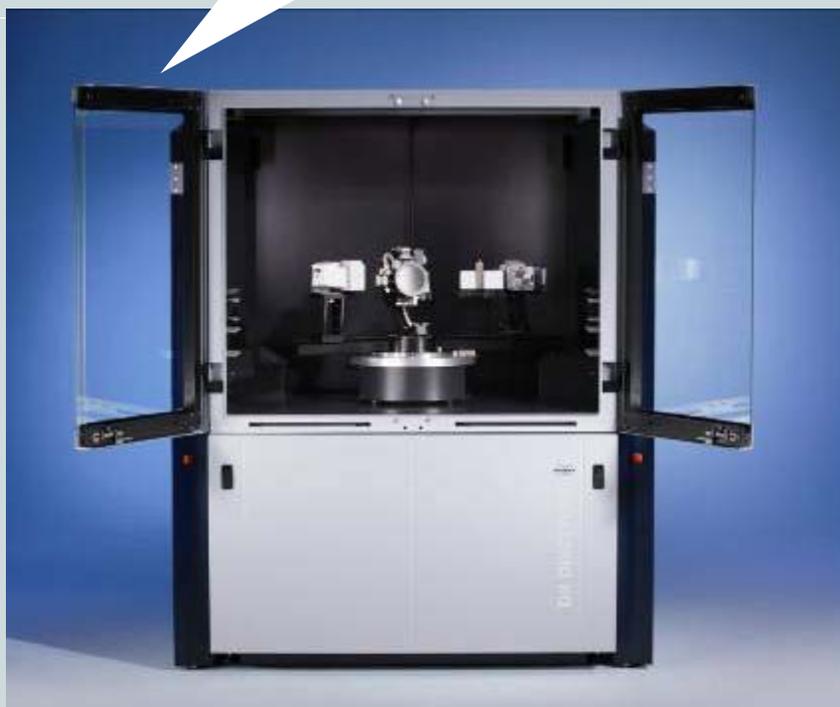

Sliding door: easy access for sample loading and configuration changes

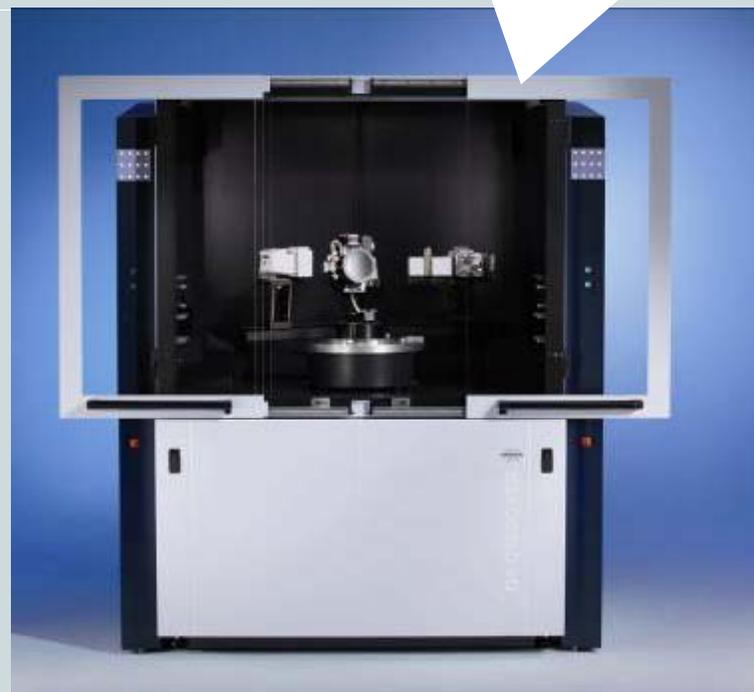

The D8 DISCOVER with DAVINCI

Tool-free mount & component recognition



DAVINCI.SNAP-LOCK
Tool-free change of optics



DAVINCI.MODE
Component recognition

SAXS




- $q_{\min} = 0.025 \text{ \AA}^{-1}$
- $d_{\max} = 25 \text{ nm}$

He beampath, SAXS beamstop,
Vantec 500

Configure: GADDS HTS

Vertical theta-theta, Reflection/Transmission

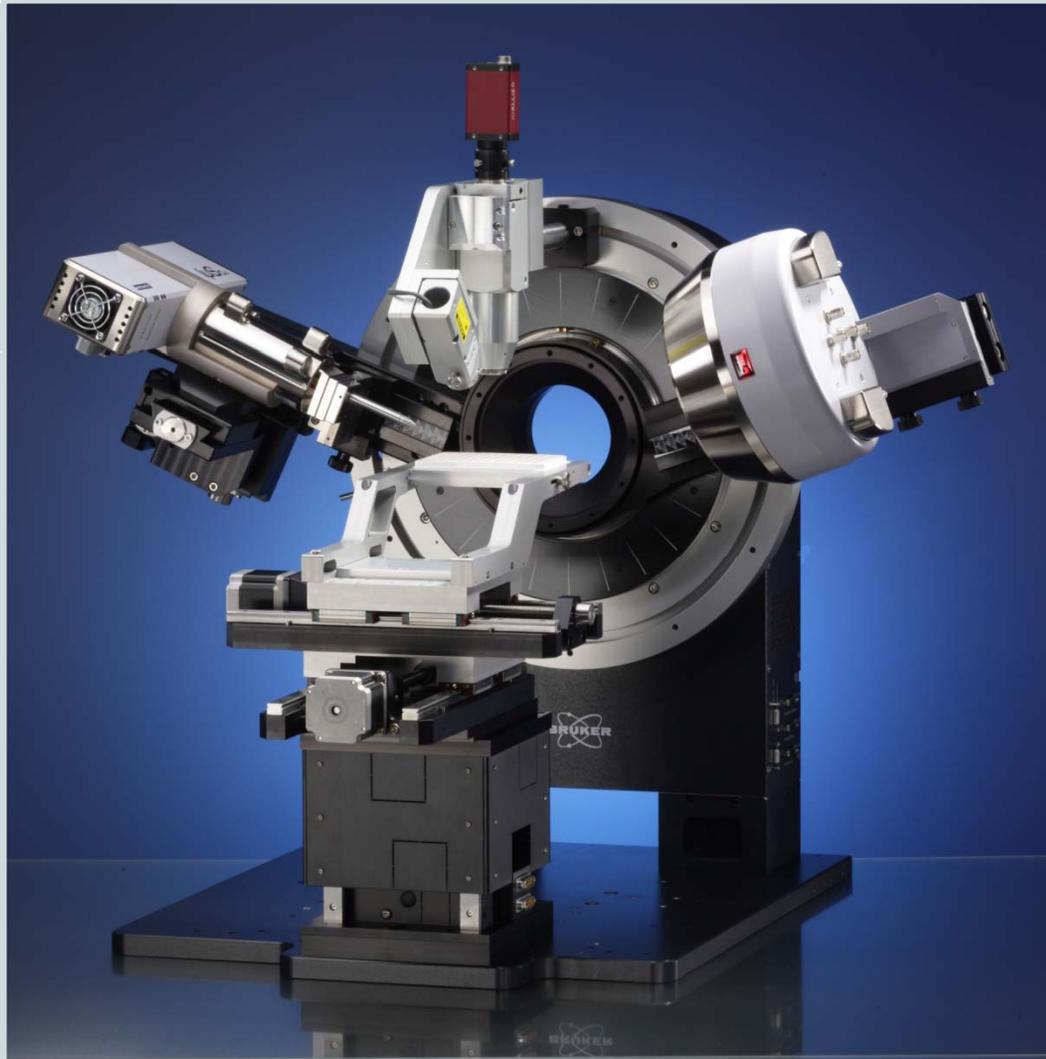

No barrier between 0D/1D/2D

Vertical theta-theta, CEC for microdiffraction/stress/texture, 0D->1D->2D



XRD²: Systems with VANTEC-500 Detector



USER DA VINCI TOOLS JOBLEST. START JOBS CONFIGURATION DBMANAGEMENT

Primary Beam Path Secondary Beam Path

1 ✓   1 ✓

Tube Cu VANTEC-500 2-D mode

2 ✓ 

POLYCAP Parallel-beam

3 ✓ 

Slit 0.6 [mm]

4 ✓ 

Soller mount empty

5 ✓ 

Slit 0.6 [mm]

6 ✓ 

Micro mask 0.5 [mm]


7 ✓ 

UBC collimator 0.5 [mm]



Sample Stage

1 ✓ 

Centric Eulerian cradle

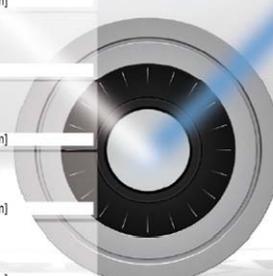

XRD² & Single Crystals

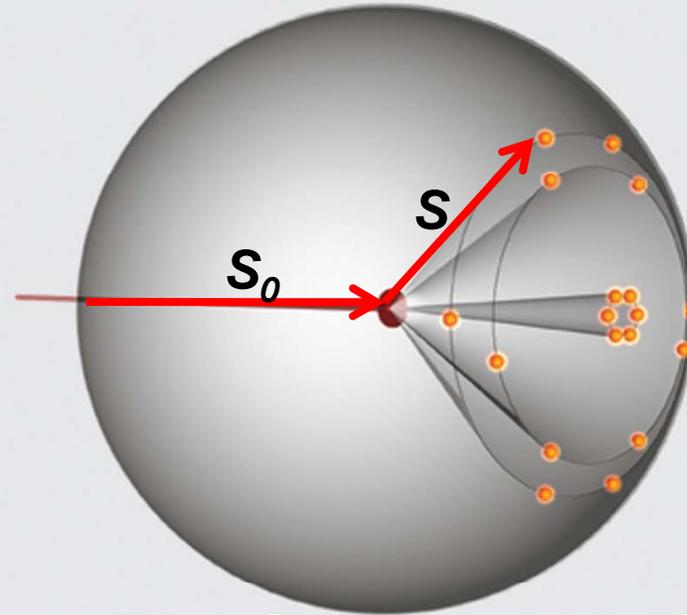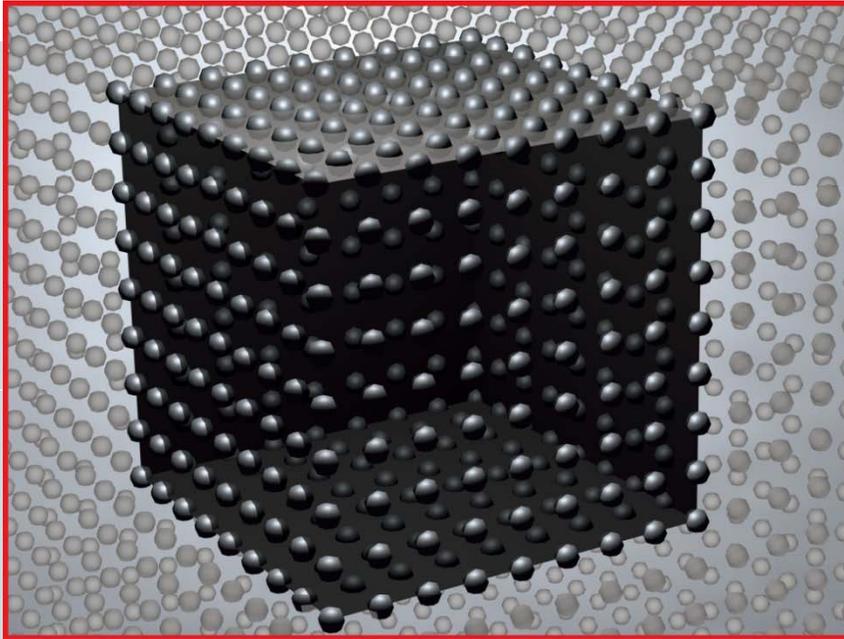


Laue equation


$$a \cdot (s - s_0) = h\lambda$$

$$b \cdot (s - s_0) = k\lambda$$

$$c \cdot (s - s_0) = l\lambda$$

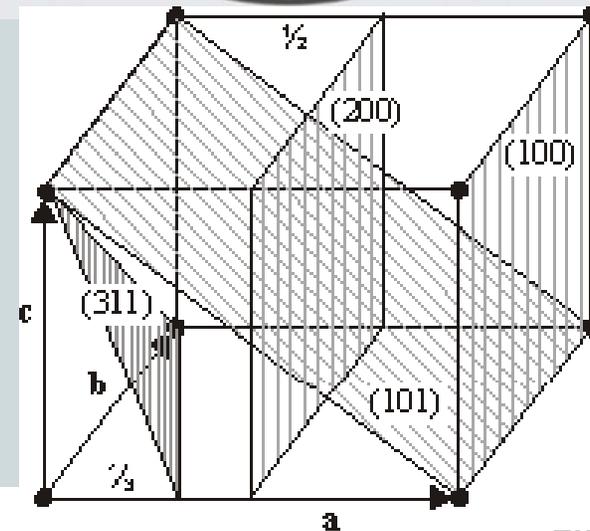


XRD² & Powders

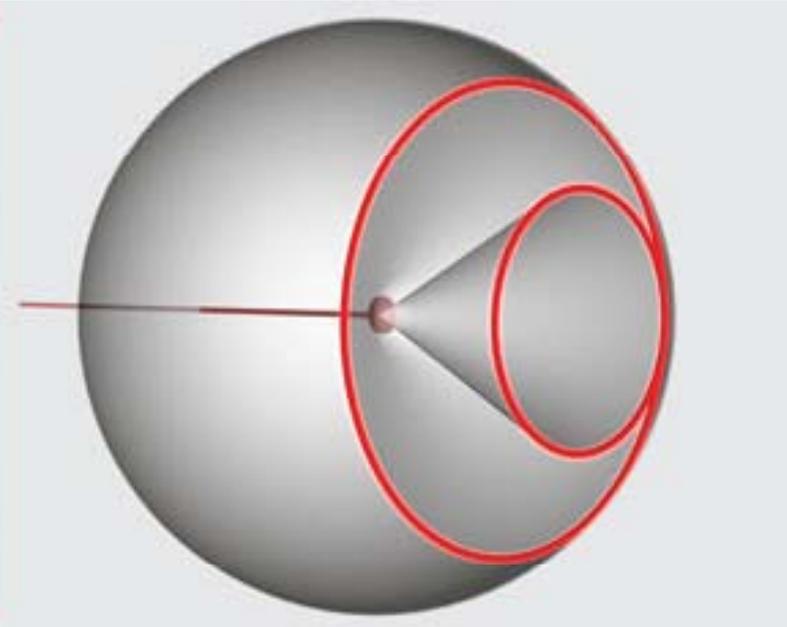


Bragg law

$$n\lambda = 2d \sin \theta$$

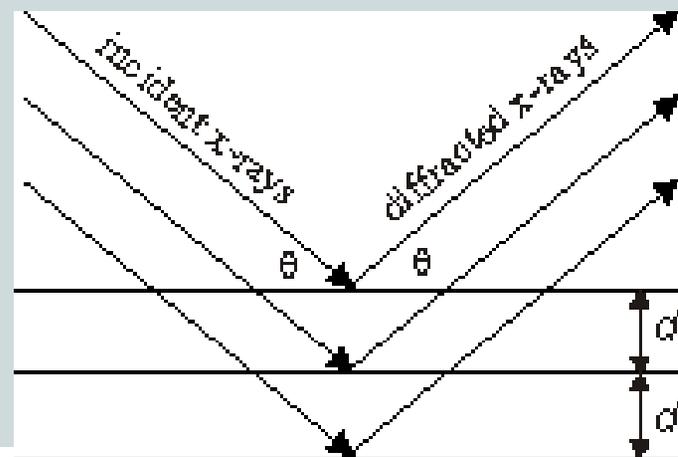


XRD²: Diffraction pattern with both γ and 2θ information

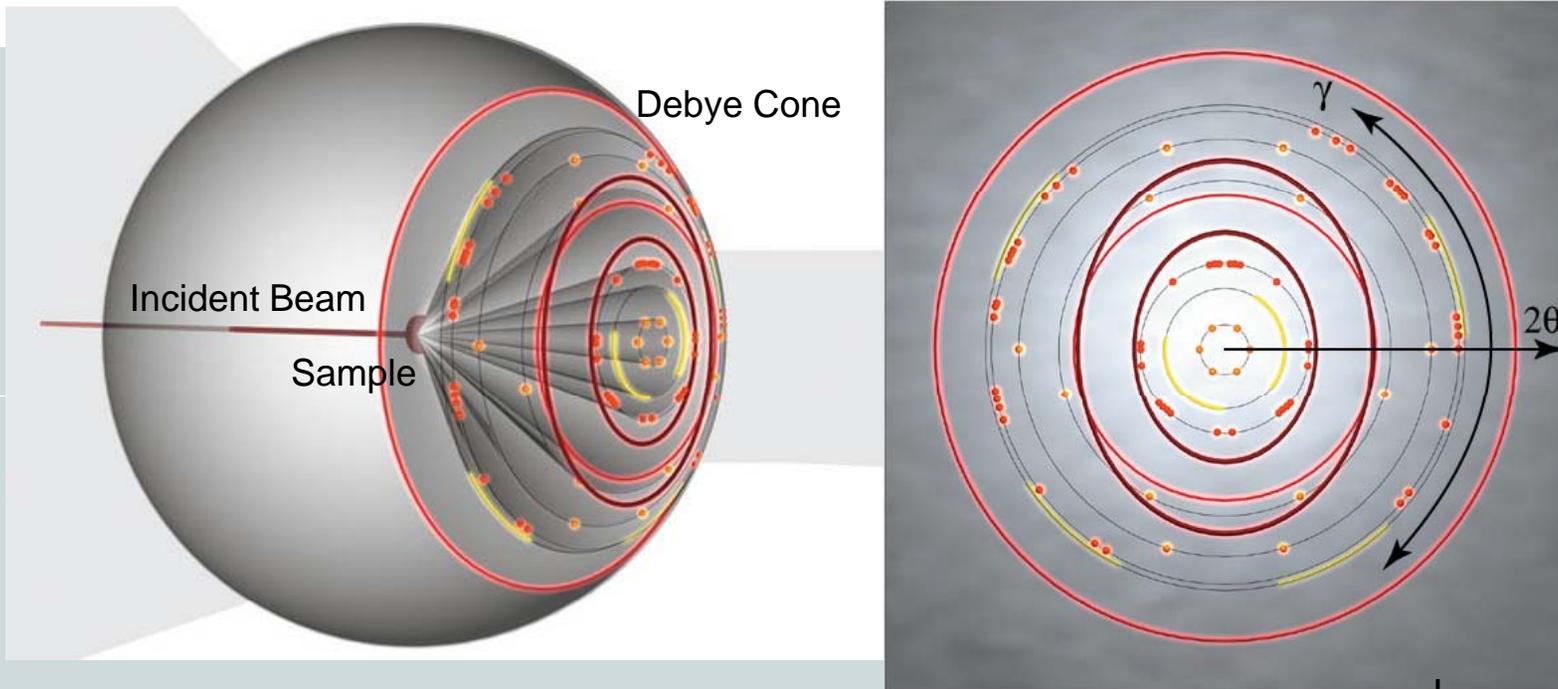

Diffraction vector with γ


$$\mathbf{H} = \frac{\mathbf{s} - \mathbf{s}_0}{\lambda} = \frac{1}{\lambda} \begin{bmatrix} \cos 2\theta - 1 \\ -\sin 2\theta \sin \gamma \\ -\sin 2\theta \cos \gamma \end{bmatrix}$$

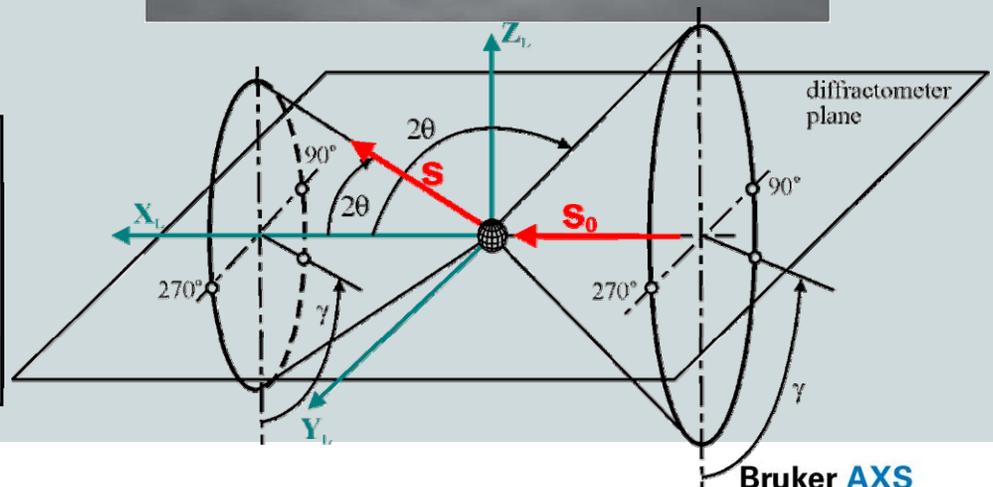


XRD²: Diffraction vector approach

Applications	Vector approaches
Phase Identification:	Polarization and absorption correction
Texture Analysis:	Orientation mapping angles; Data collection strategy (scheme)
Stress Measurement:	Fundamental equation derived by second order tensor transformation; Data collection strategy (scheme)
Crystal Size Analysis:	Equations for the effective volume calculation at both reflection and transmission modes.


XRD²: PhaseID Measurement Geometry

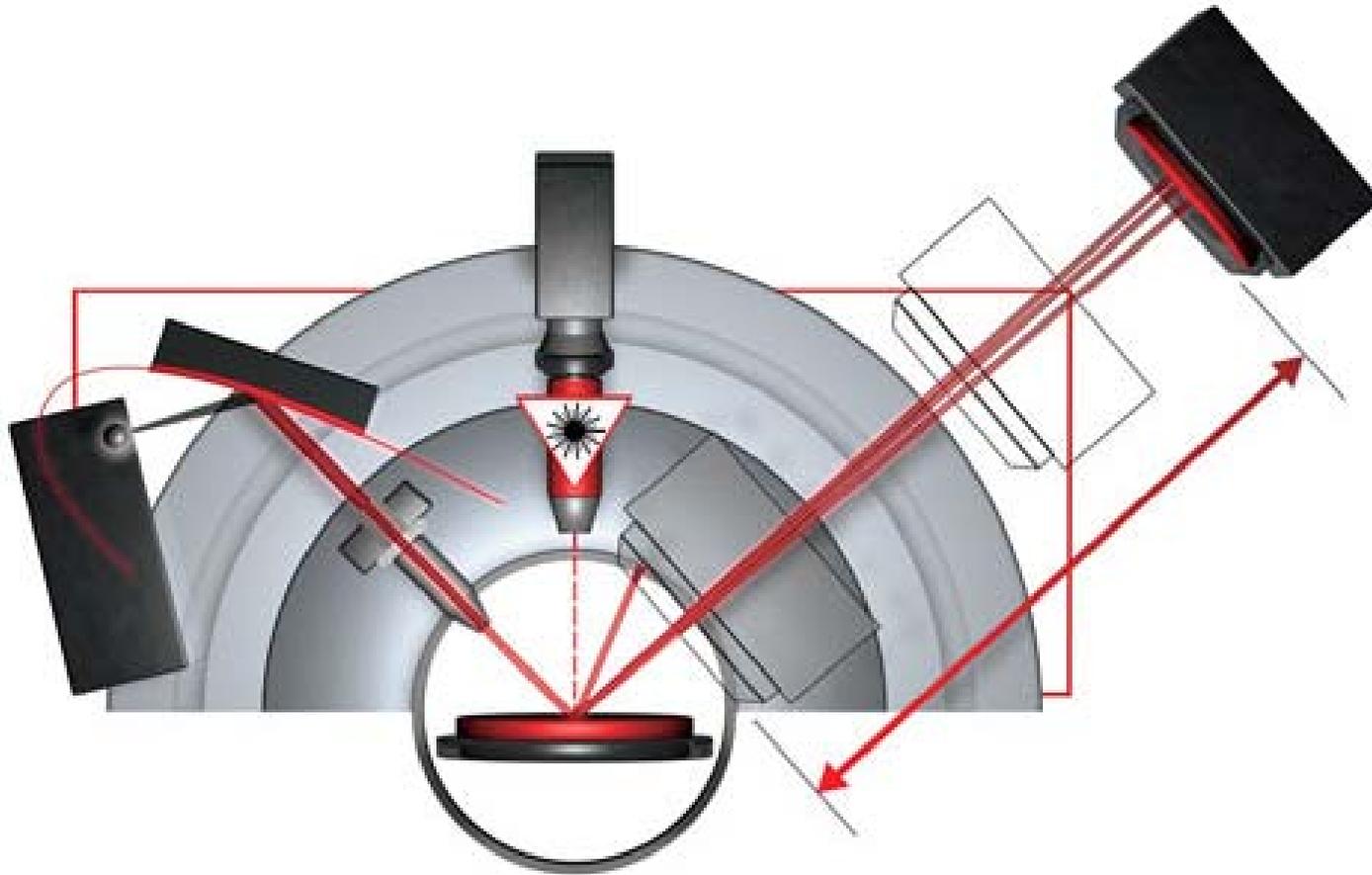



XRD²: Single Frame Covering All

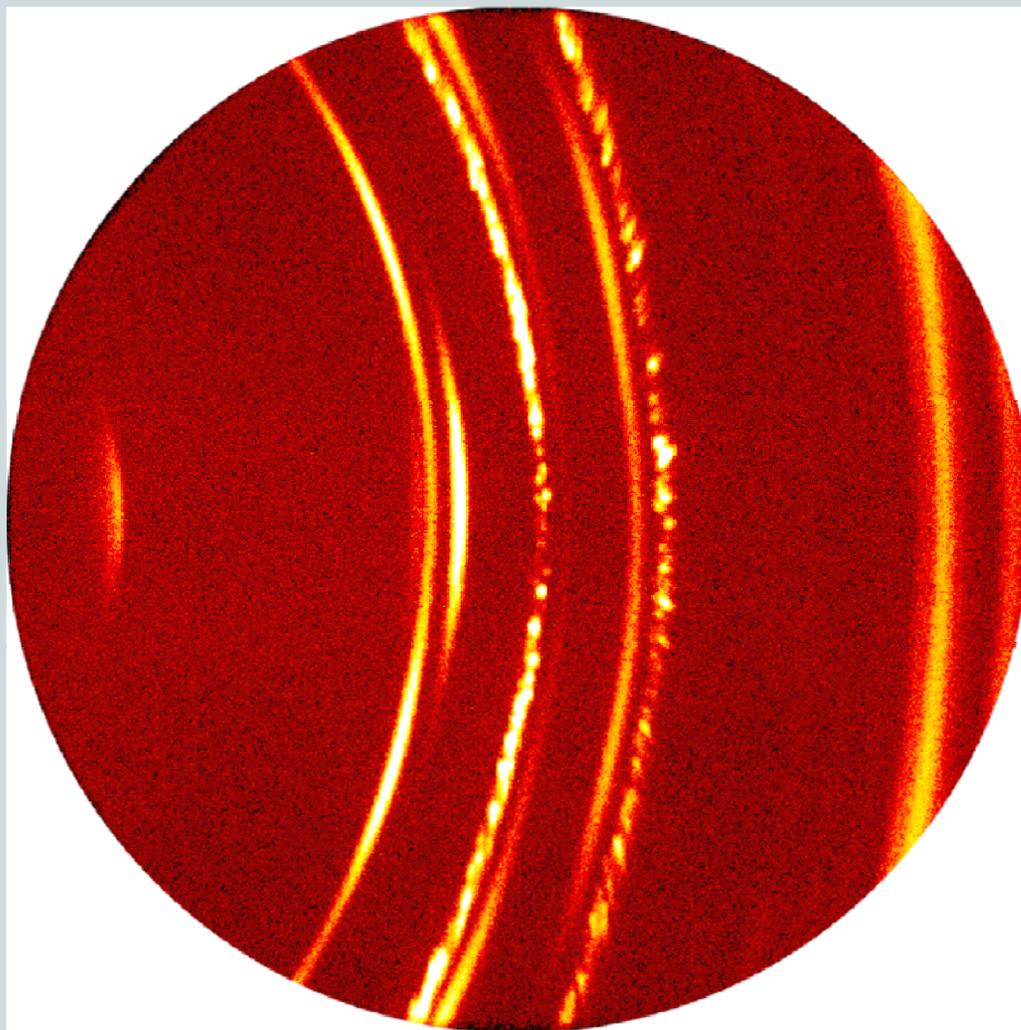

XRD²: Frame Merge and Integration

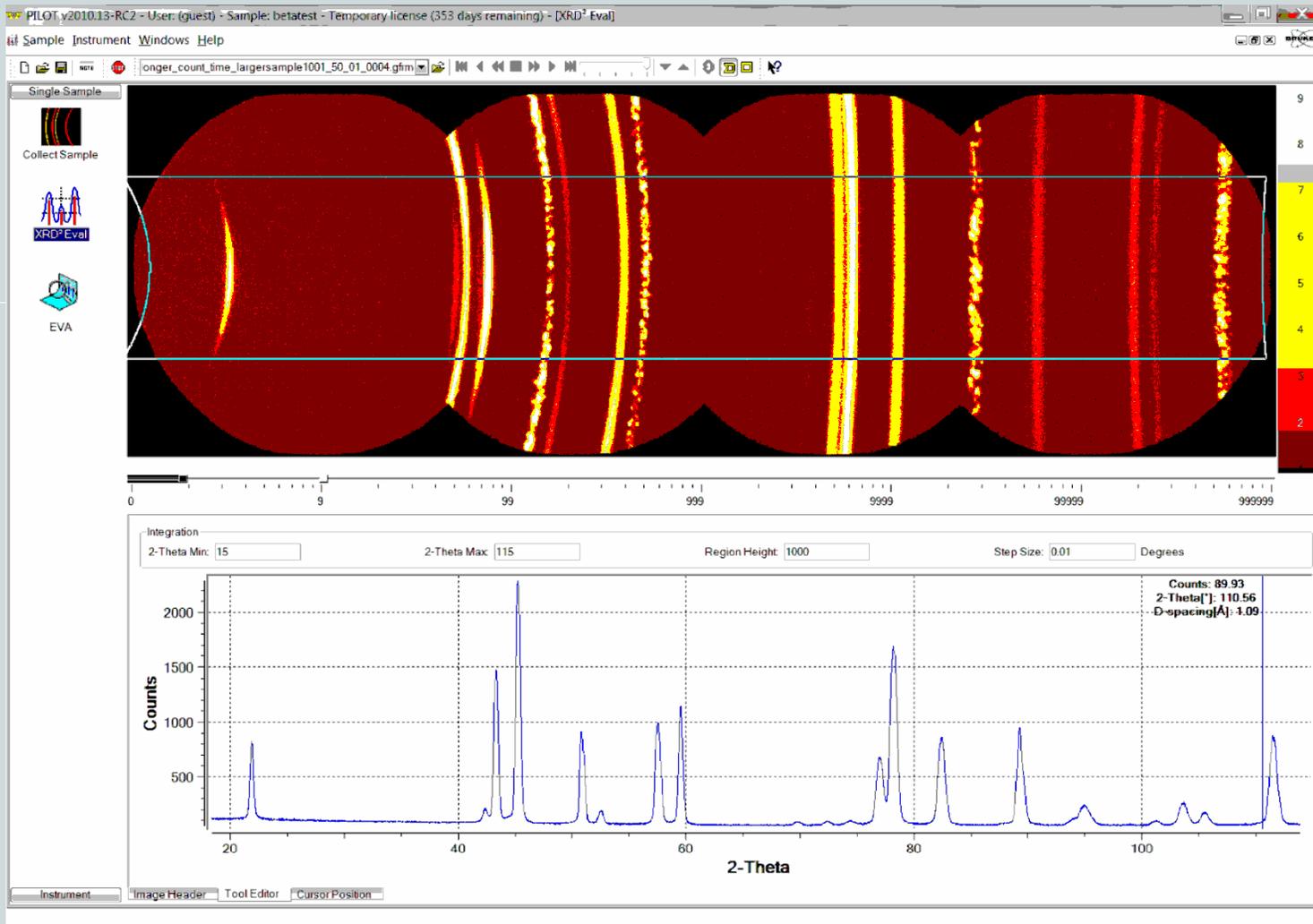

XRD²: Fundamental Equation for Texture Analysis

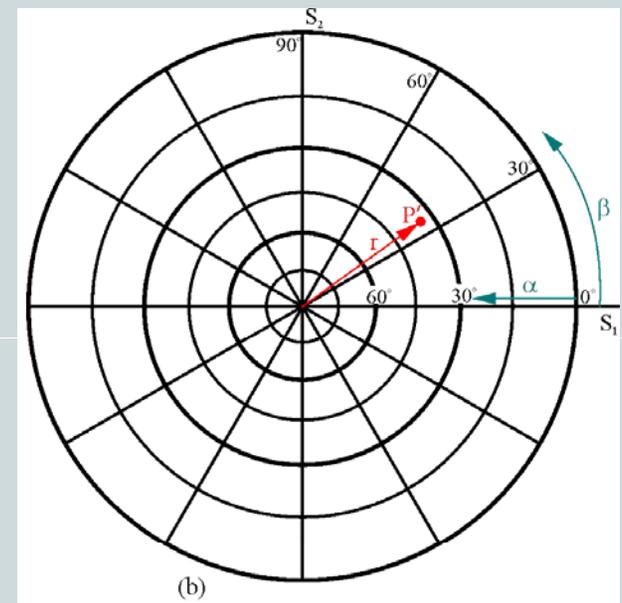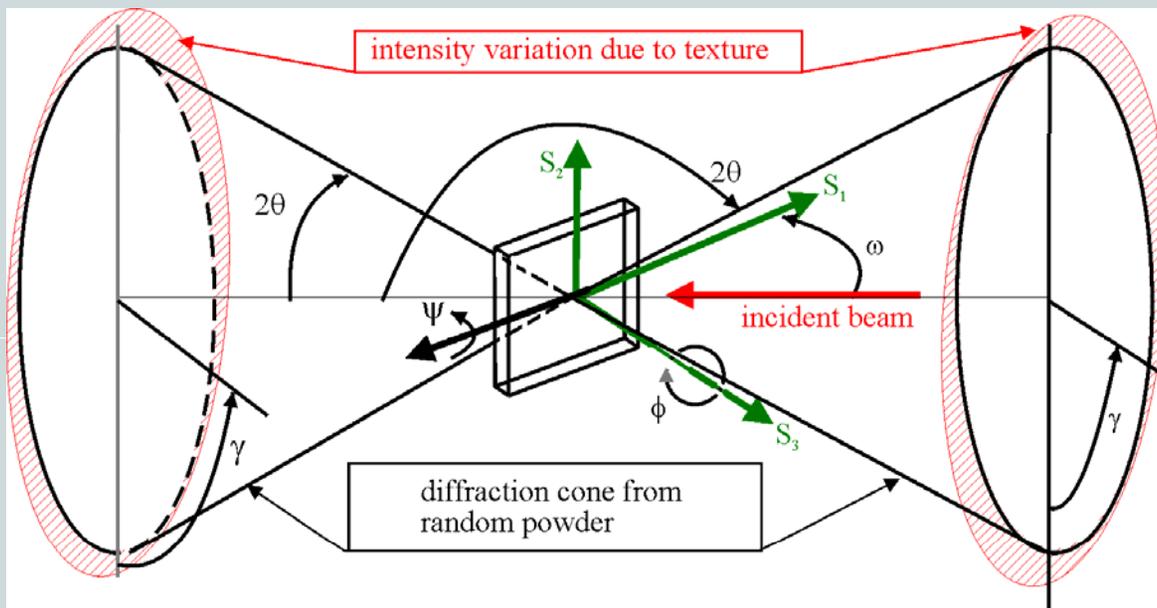


The pole figure angles (α, β) can be calculated from the unit vector components by the pole mapping equations:


$$\alpha = \sin^{-1} |h_3| = \cos^{-1} \sqrt{h_1^2 + h_2^2}$$

$$\beta = \pm \cos^{-1} \frac{h_1}{\sqrt{h_1^2 + h_2^2}} \quad \begin{cases} \beta \geq 0^\circ & \text{if } h_2 \geq 0 \\ \beta < 0^\circ & \text{if } h_2 < 0 \end{cases}$$


The D8 DISCOVER with DAVINCI VÅNTEC-500 for texture measurement

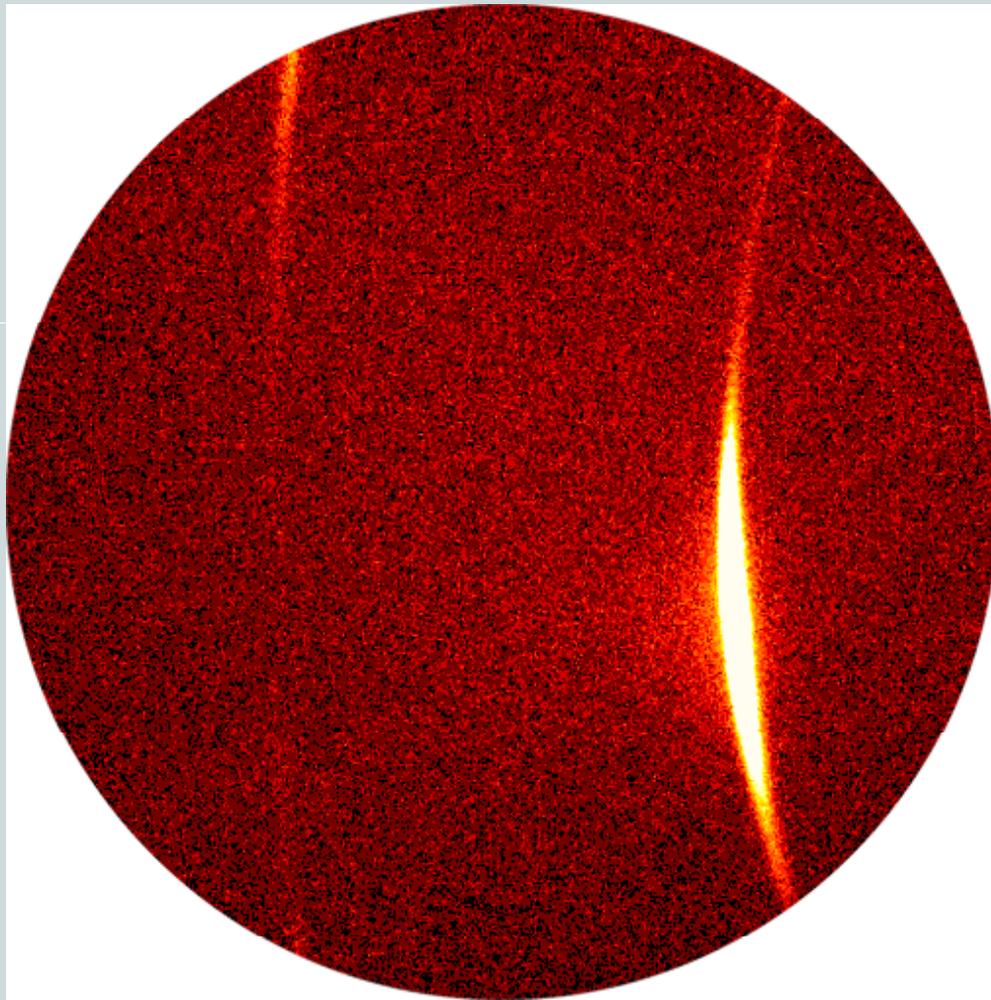


- Steel can
- (200) & (110) rings
- Intensity variation during ϕ scan

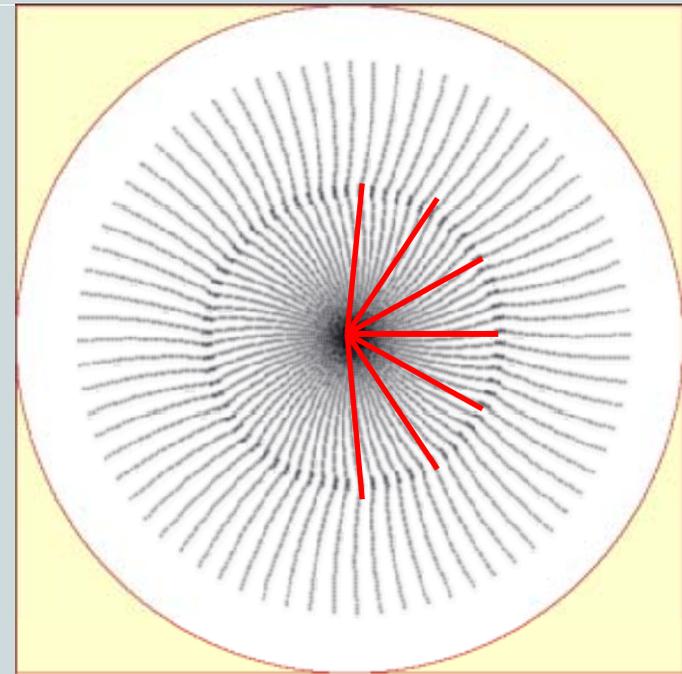


XRD²: Particle size measurement by 2θ & γ profile analysis:

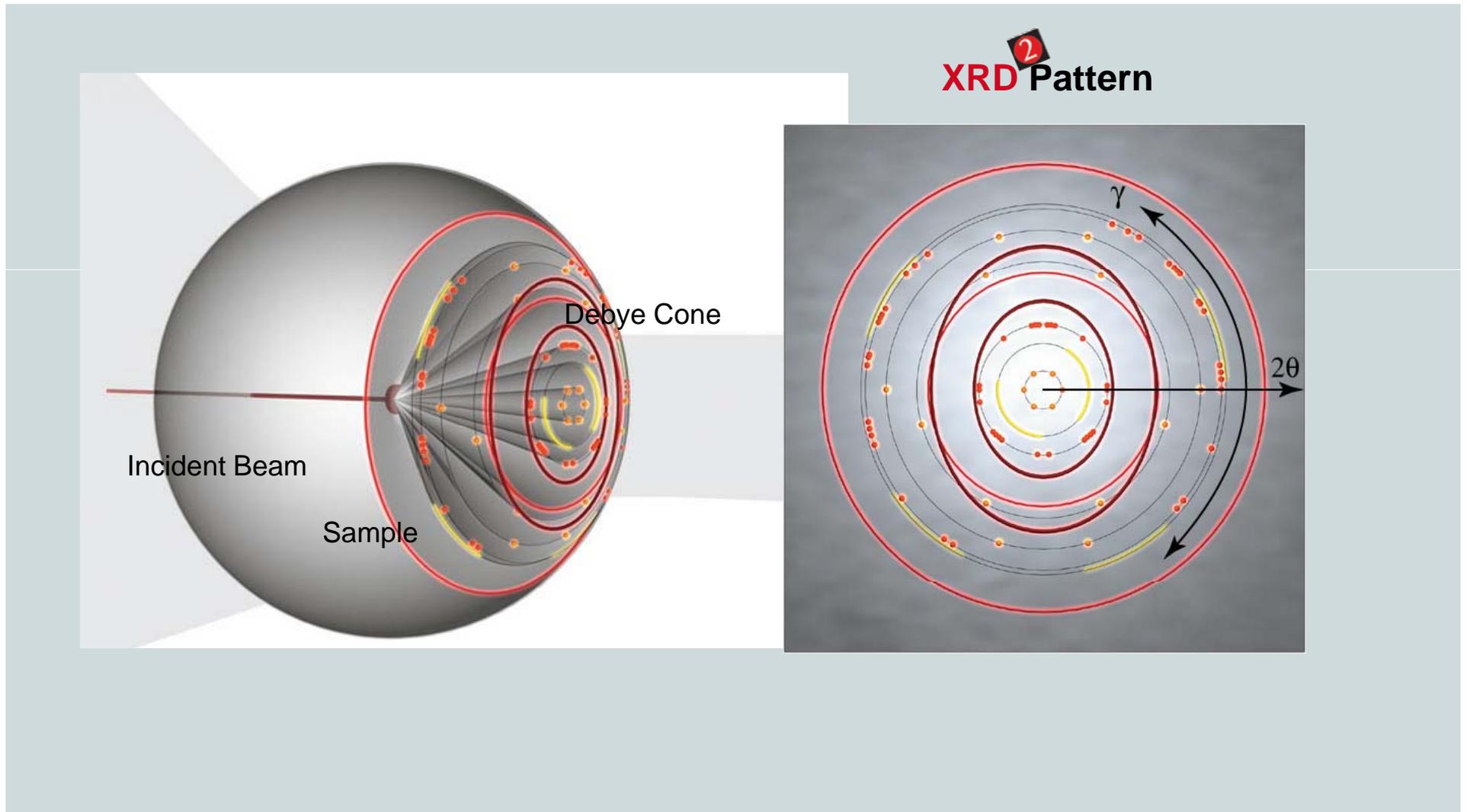

XRD²: Peak broadening-gold Nanoparticles

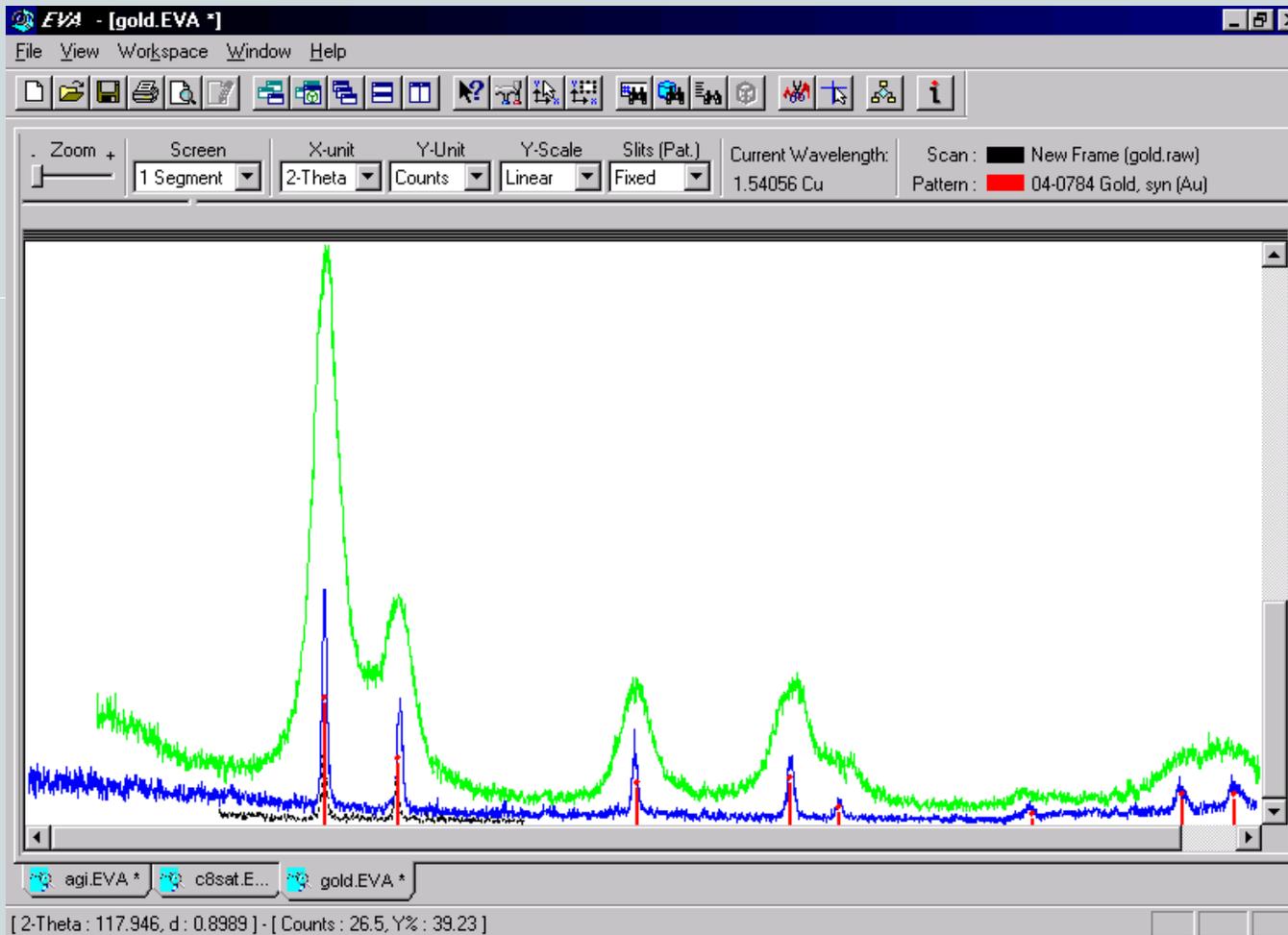

XRD²: Particle size and instrument broadening:

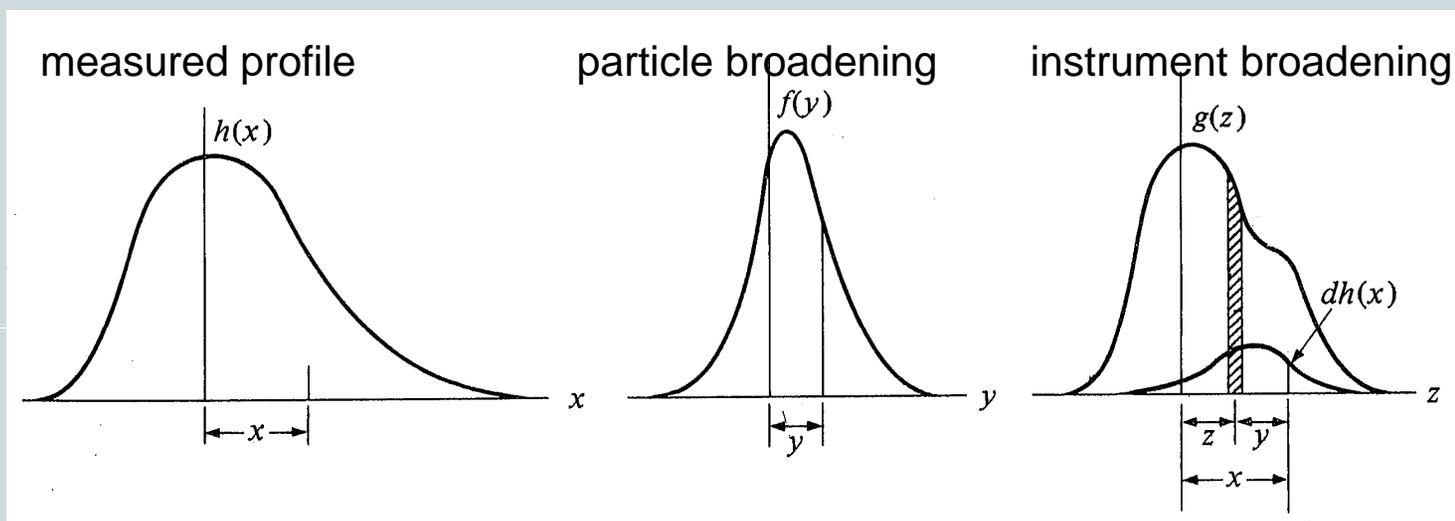

The measured profile is a convolution of the functions representing particle-size broadening and instrument broadening

$$h(x) = \frac{1}{A} \int g(z) f(x-z) dz$$

where A is the area of the $f(y)$ curve and $y=x-z$.

Ref: B. E. Warren, *X-ray Diffraction*, Dover Publications, Inc. New York, 1990.

XRD²: Particle size calculation:

Scherrer equation:

$$t = \frac{C\lambda}{B \cos \theta}$$

where λ is wavelength (Å), B is FWHM (radians) corrected for instrument broadening, θ is Bragg angle, C is a crystal shape factor from 0.9~1.

For Gaussian profiles,

$$B^2 = U^2 - S^2$$

while for Cauchy profiles,


$$B = U - S$$

where B is the corrected FWHM for crystallite size calculation by Scherrer equation, and U and S are the FWHM's of the unknown and standard peaks, respectively.


XRD²: Data Collection:

Acetaminophen powder

5 second data collection

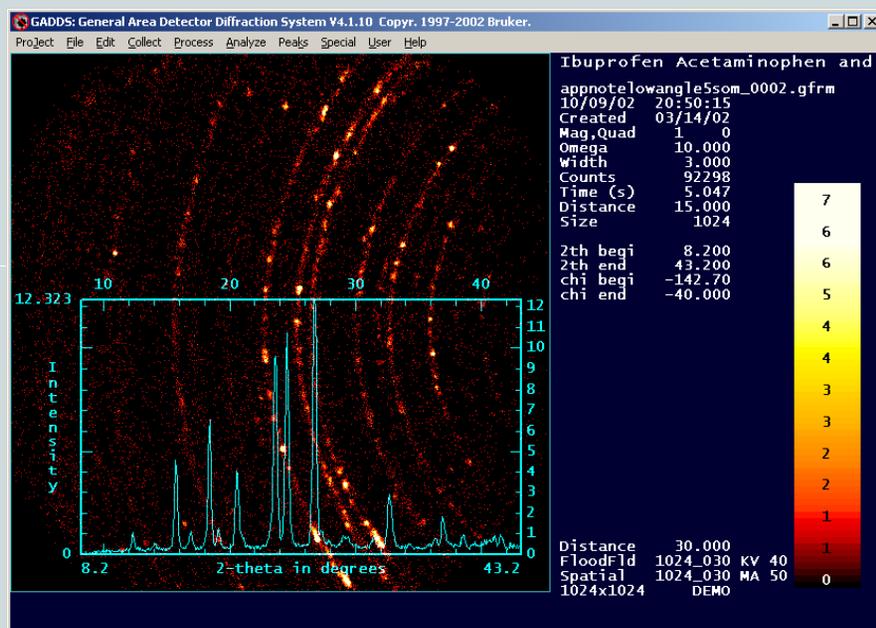

30 second data collection

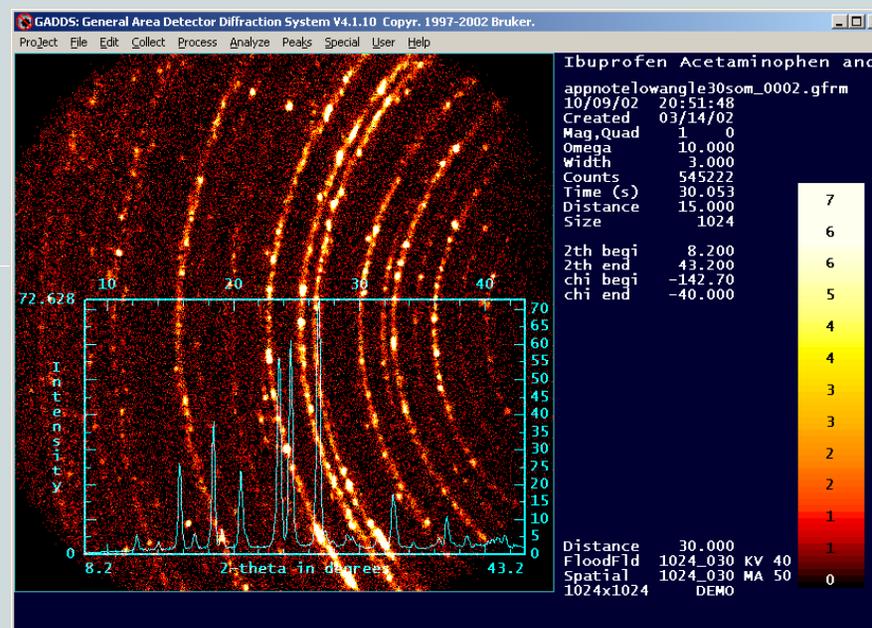

- The spotty diffraction ring is due to the large crystallites compared to the sampling volume (beam size).
- The number of spots on the ring is determined by crystallite size, instrumental window (γ -range), multiplicity of the crystal plane, and effective diffraction volume.



XRD²: Particle size measurement by γ profile analysis:

For XRD² in reflection mode, the particle size is given by

$$d = k \left\{ \frac{p_{hkl} b^2 \arcsin[\cos \theta \sin(\Delta\gamma / 2)]}{2\mu N_s} \right\}^{1/3}$$

where k is the instrumental calibration factor or can be calculated

from $k = \left(\frac{3\beta}{4\pi} \right)^{1/3}$ if the instrument broadening in 2θ direction is known.


For transmission mode with the incident beam perpendicular to the sample surface, the particle size is given by

$$d = k \left\{ \frac{p_{hkl_i} b^2 t \arcsin[\cos \theta \sin(\Delta\gamma / 2)]}{N_s} \right\}^{1/3}$$

where k is the instrumental calibration factor or


$$k = \left(\frac{3\beta}{4\pi} \right)^{1/3}$$

XRD²: Particle size measurement by γ profile analysis:

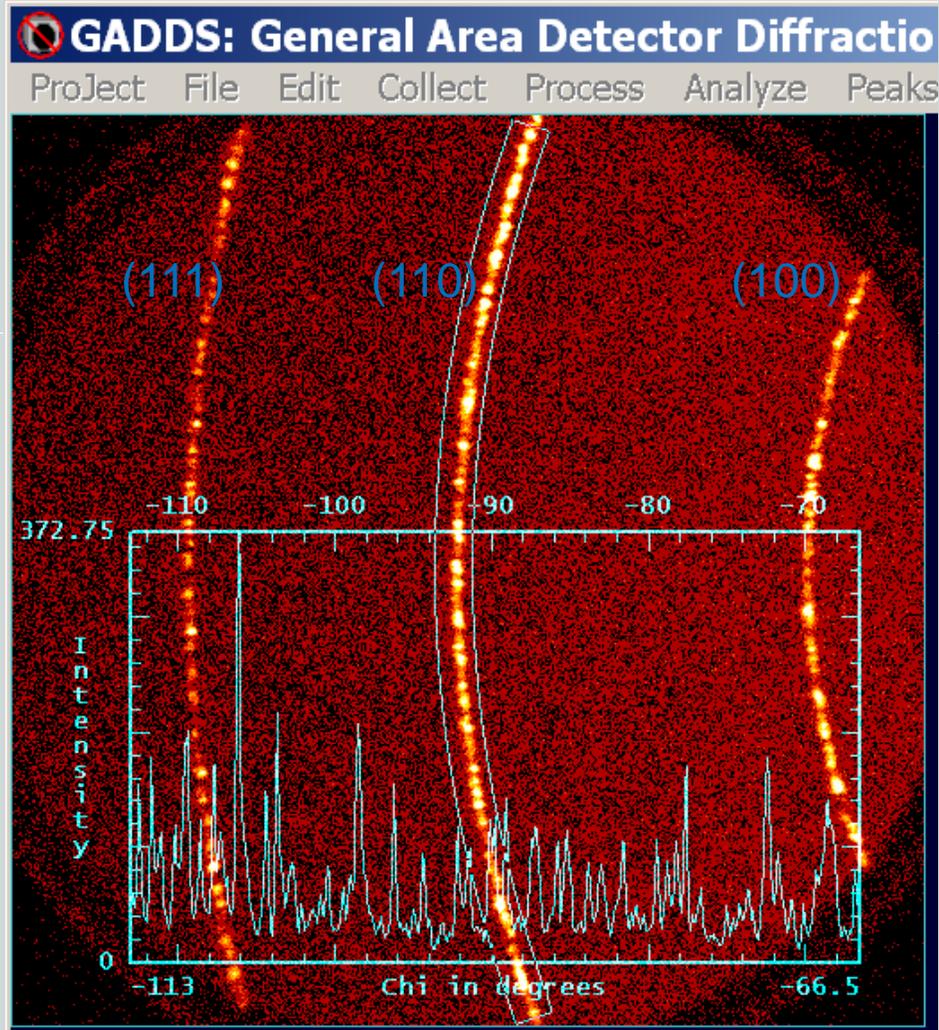

- The frame was collected from a SRM660a (LaB6) sample in transmission mode and Cu-K α x-rays.
- The 2D detector (Hi-StarTM) is set at 23.75 cm from the instrument center.
- The beam (collimator pinhole) size b is 200 μm .
- The sample thickness t is 7.0 μm , based on the calculated μ of 1138 cm^{-1} and the measured transmission of 0.45.

XRD²: Particle size measurement by γ profile analysis:

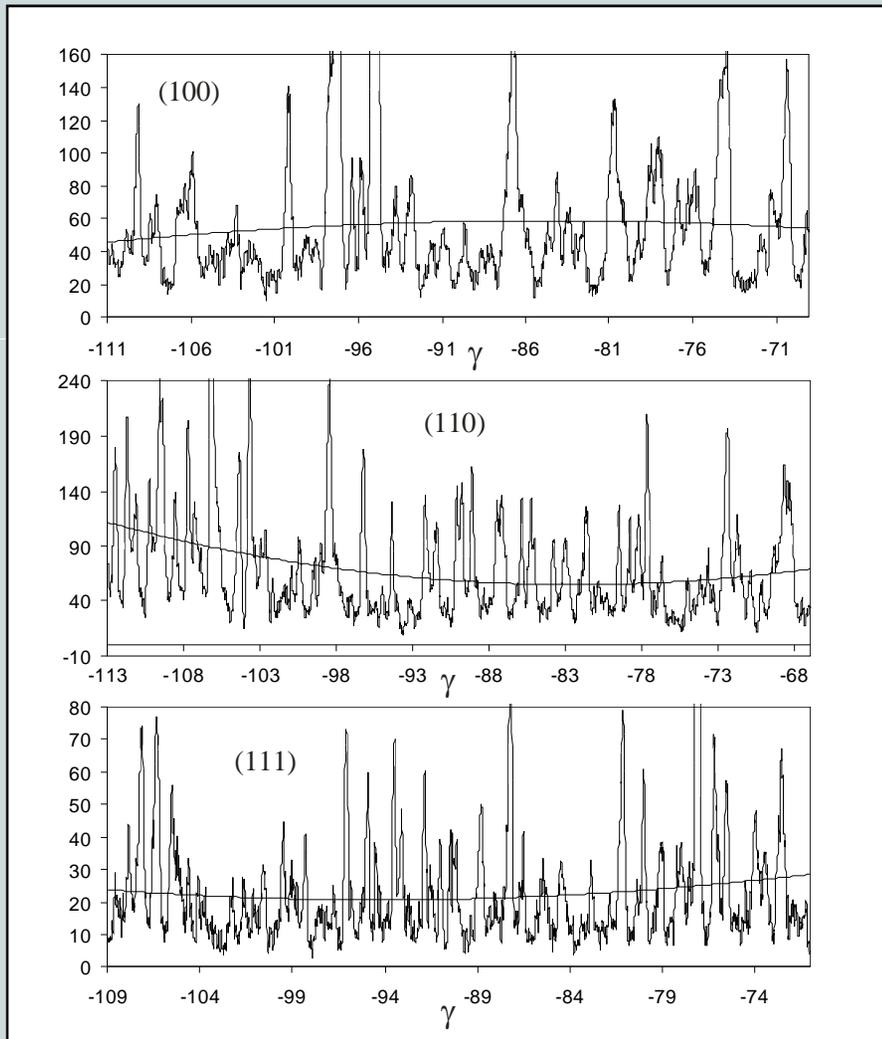

- The 2θ-integrated plots (γ -profiles) of three rings from (100), (110) and (111) planes are displayed.
- The number of crystallites is counted from the number of intersections of the γ -profile with a threshold line.
- To cancel out the effects of the overall intensity fluctuation (texture, etc.), a 2nd order polynomial trend line is fitted to each γ -profile as a threshold line.
- Every two intersections of γ -profile with the threshold line represents a crystallite.
- New analysis strategy?
- Size distribution?




XRD²: Particle size measurement by γ profile analysis:

Calibration Results:

(hkl)	P_{hkl}	2θ	$\Delta\omega$	N_s	k
(100)	6	21.36	38	23	0.1217
(110)	12	30.38	46	41	0.1106
(111)	8	37.44	42	38	0.1281

- The average scaling factor k is 0.12 in this calibration. The system can then be used to measure the crystallite size of unknown materials if the data can be collected in approximately the same condition.
- It is always necessary to calibrate the system with a known standard, preferably with a comparable sample geometry and crystallite size.
- For reflection mode, it is critical to have a standard with a comparable linear absorption coefficient so as to have similar penetration.

XRD²: Particle Size Analysis by X-ray Diffraction:

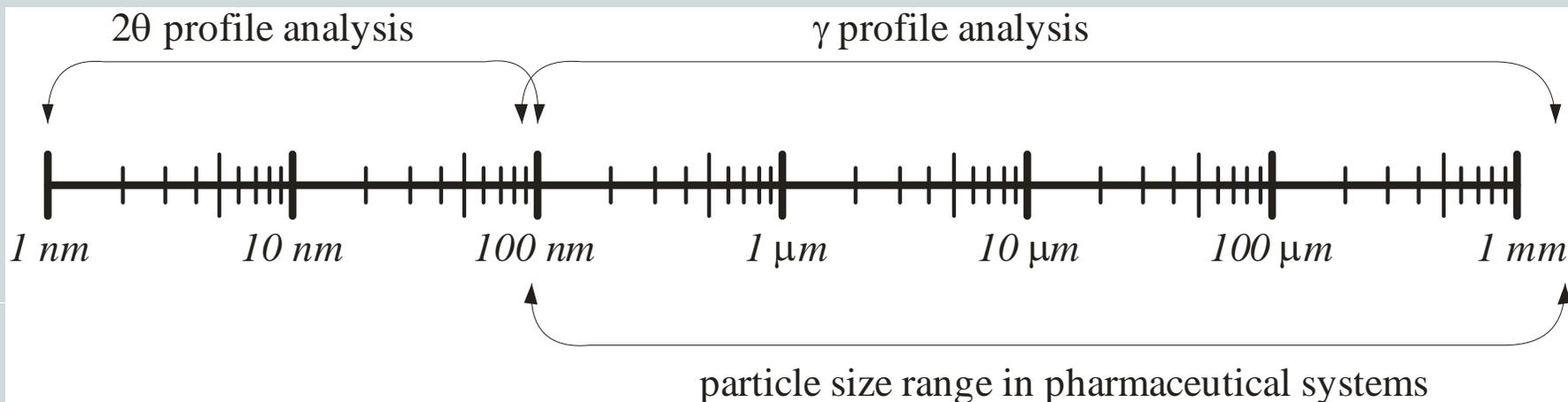


- 2θ profile analysis, including measurement from peak FWHM by Scherrer equation, or profile analysis by Stokes and Wilson, is suitable for particle size below 100 nm.
- γ profile analysis, based on sampling statistics, is suitable for particle size from sub-micrometer to a few millimeters.
- The particle size range of pharmaceutical substances is from sub-micrometer to a few millimeters.

More About XRD²



1. Introduction.
2. Geometry Conventions.
3. X-Ray Source and Optics.
4. X-Ray Detectors.
5. Goniometer and Sample Stages
6. Data Treatment.
7. Phase Identification.
8. Texture Analysis.
9. Stress Measurement.
10. Small-Angle X-Ray Scattering.
11. Combinatorial Screening.
12. Quantitative Analysis.
13. Innovation and Future Develop

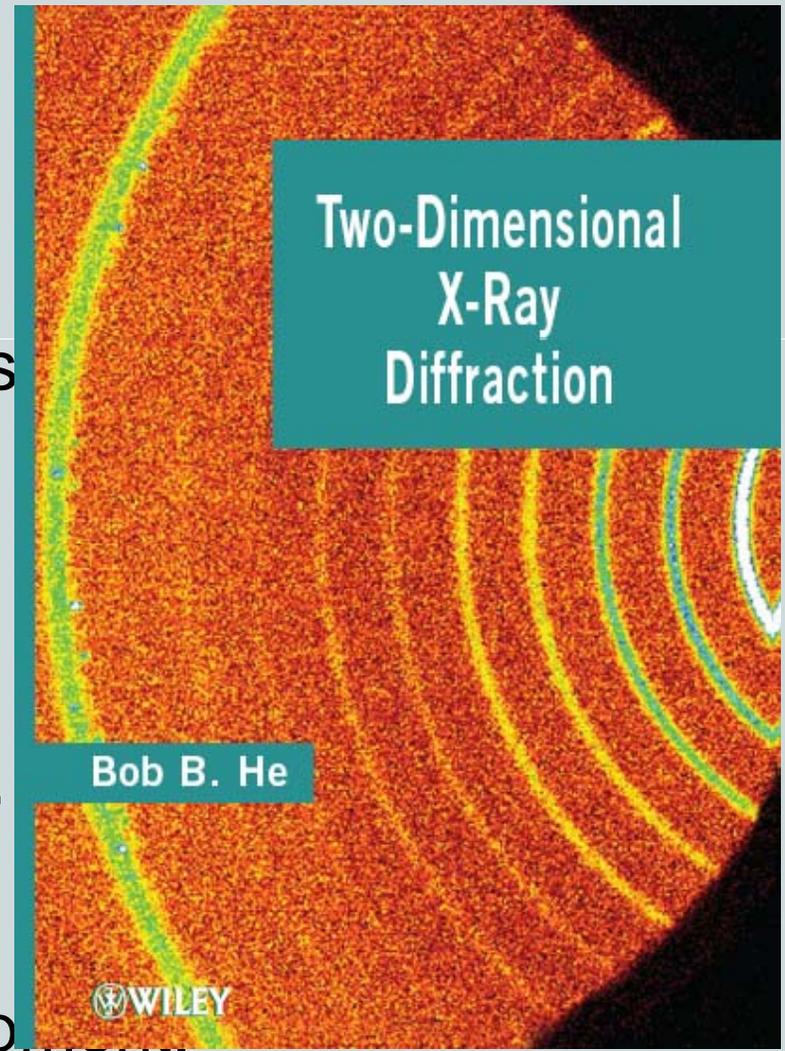



Thank You for Your Attention