

New Instrumentation For Pharmaceutical XRD/SAXS



Bob He

Director of Innovation & Business Development XRD²

Bruker AXS Inc.



This document was presented at PPXRD - Pharmaceutical Powder X-ray Diffraction Symposium

Sponsored by The International Centre for Diffraction Data

This presentation is provided by the International Centre for Diffraction Data in cooperation with the authors and presenters of the PPXRD symposia for the express purpose of educating the scientific community.

All copyrights for the presentation are retained by the original authors.

The ICDD has received permission from the authors to post this material on our website and make the material available for viewing. Usage is restricted for the purposes of education and scientific research.



PPXRD Website – www.icdd.com/ppxrd

ICDD Website - www.icdd.com

2-D XRD/SAXS

Which information can be extracted?



2-D WAXS (XRD)

- Texture
- Crystalline phases
- Strain
- Crystallinity
- Crystallite size
- Amorphous
- ...

SAXS and WAXS are **complementary** analysis methods that help better understanding samples or processes, e.g.

- Arrangement of crystals in polymers – types (spherulites, shish ...)
- Density distribution between crystal and amorphous zones
- Thermal stability and processability
- Deformation processes

2-D SAXS

- Typical size, size distribution
- Distance between particles
- Arrangement of nano-particles
- Texture of large structures
- Details about system (volume fraction degree of crystallinity)
- Porosity
- Shape and dimensions of nano-particles
- Strain
- ...

What does one want in a (home lab) XRD/SAXS system?




- Source
 - Highest flux desired to data collection time and deal with small sample volume, weak diffraction and low concentration
 - Point focus preferred with area detector
- Optics and beam path
 - Suitable to the applications
 - Easy to switch between configurations
- Detector
 - Must have high quantum efficiency
 - Very low noise, photon-counting detector preferred
 - Larger is better (to see the entire scattering pattern)
- Overall: high reliability, easy use and low maintenance

Sources

Characteristic X-ray generation

- All present day monochromatic home laboratory sources are based on characteristic radiation from a material anode
- The efficiency of this process is very low
 - Approximately 99% of the incident electron power is converted to heat, *not X-rays*
- Dissipation of this waste heat fundamentally limits the brightness of the source

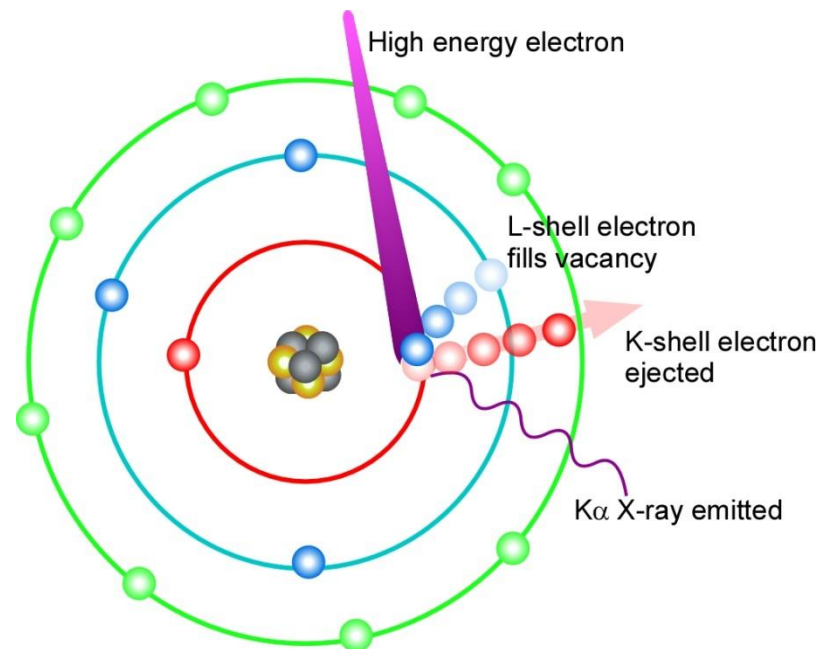


How can one make a source brighter?

- Source performance is ultimately limited by anode melting
 - There are three ways to improve performance
- Make the focus on the anode smaller
 - Allows higher heat extraction efficiency
 - ***Microfocus sources***
- Rotate the anode faster
 - Spreads the heat more efficiently
 - ***Improved rotating anode***
- Use a liquid metal anode
 - Can't melt (it is already molten!)
 - ***Metal jet sources***

How to make brighter source I: Microfocus sources

- Brightness (**B**) is proportional to power loading (*p*)
- Power loading is higher for *smaller spot focus*

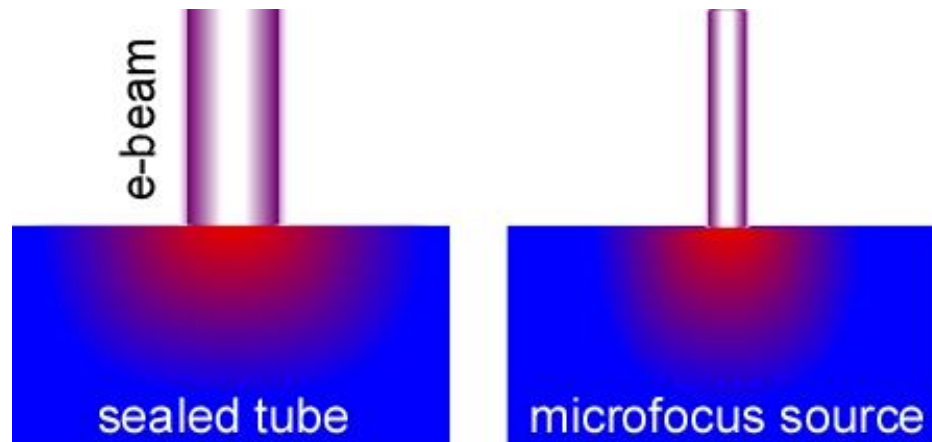

$$p_{\max} = \frac{2\kappa(T_m - T_0)}{r\sqrt{\pi \ln(2)}}$$

- | | |
|--|--|
| ▪ Large spot | ▪ Small spot |
| ▪ Quasi-one dimensional heat flow limits power loading | ▪ Two dimensional heat flow (more efficient cooling) |
| | ▪ Relative performance improved by 10 times |

I μ S microfocus source



- Intensity 3×10^{10} X-rays/mm²-sec (Cu K α)
 - 8 times higher than conventional 5.4 kW rotating anode
- Typical lifetime >5 years
 - High reliability
 - 3 year warranty
 - >300 installed
- Air-cooled
- Available in Cr, Cu, Mo, Ag




Comparison: Ibuprofen I μ S & VÅNTEC-2000 vs. Classical set-up



Sealed Tube

- 0.3 mm collimator
- Sample-Detector distance 29 cm


120 sec collection time

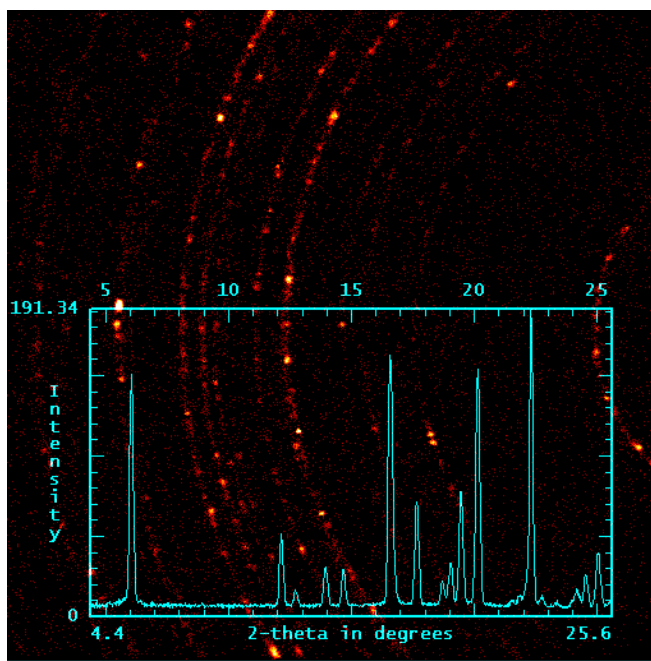

I μ S – XRD² – foc


- 2mmX2mm on sample, and 200um spot focused on detector
- small slice for integration to obtain better resolution

15 sec collection time

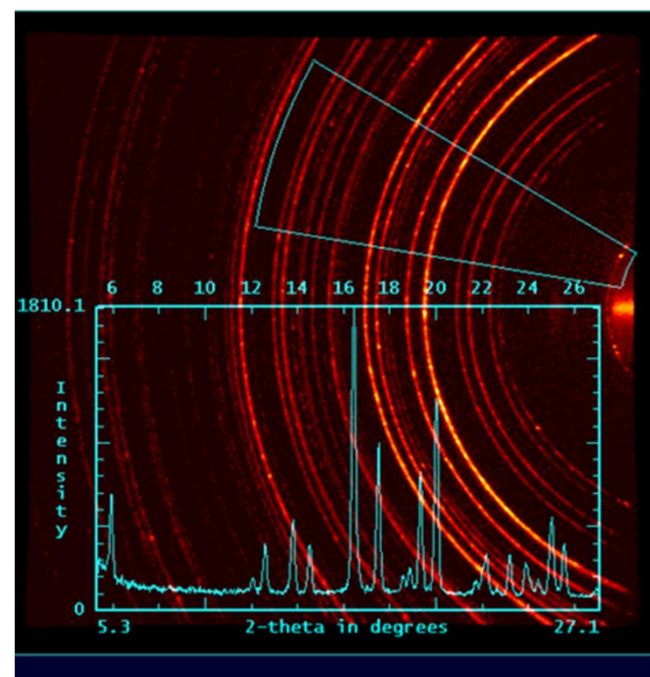

How to make a brighter source II: Rotating anode sources

- Power loading can be increased by over an order of magnitude by rotating the anode surface to spread out the heat load
 - Power load is also (modestly) increased by smaller spot
- In latest generation rotating anodes angular velocity is 10,000 rpm
- This improves performance by 50 times

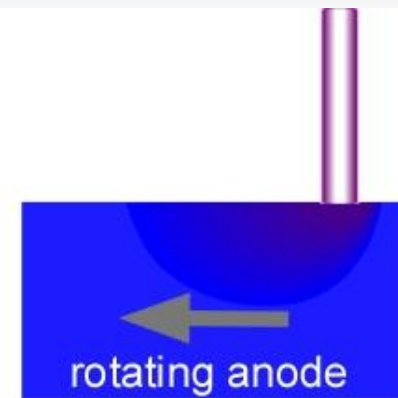


$$p_{\max} \propto \kappa (T_m - T_0) \sqrt{\frac{v}{w}}$$

w =beam width
 v =anode velocity

TXS HB High brilliance rotating anode

- Highest intensity rotating anode
 - 2×10^{11} X-rays/mm²-sec Cu K α
 - 50 times the intensity of a 5.4 kW classic RAG with multilayer optics
- Cu, Mo, Ag anodes
- Low maintenance
- Easy to align, highly stable mount
 - No alignment base
- Precrystallized, prealigned filaments
 - *No realignment required after filament changes (2X per year)*
- Precision aligned anode
 - *No realignment required after anode exchange (1X per year)*

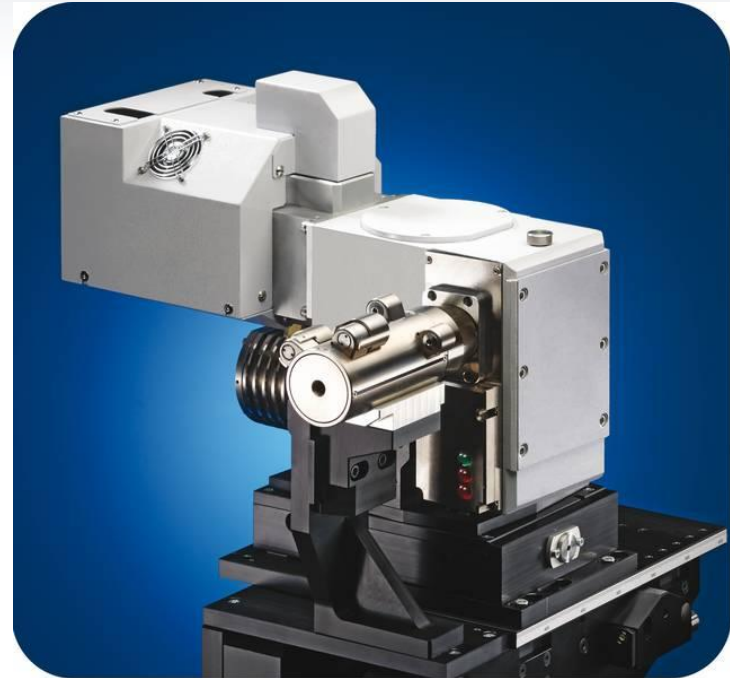

What are the limits of rotating anode performance?


- Faster rotation (ω) allows higher power loading

$$p_{\max} \propto \sqrt{\frac{\omega R}{w}}$$

- However, faster rotation also increases mechanical stress

$$\sigma_h = \frac{PR}{t} = \frac{\rho \omega^2 R^3}{2t} + \rho \omega^2 R^2$$

- State-of-the-art rotating anodes are operated close to mechanical failure limits
 - ***Little room for further improvement***


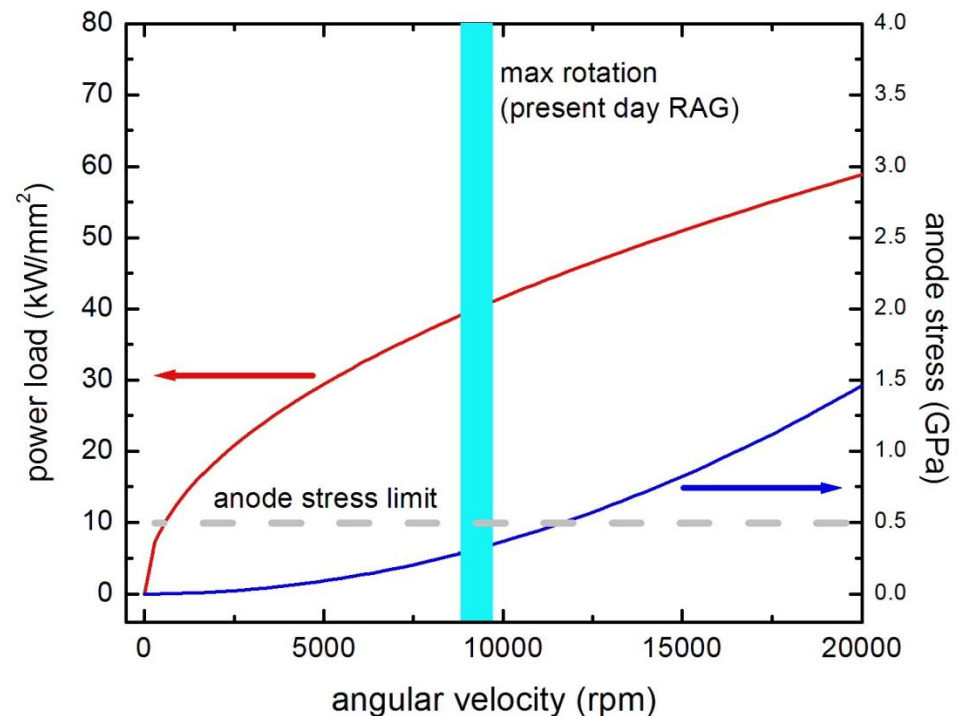


Source Comparisons

Source	Beam size (μm)	Intensity (X-rays/sec/ mm^2)	Relative Intensity
Classic 5 kW RAG	300	4×10^9	1
I μ S microfocus	110	1.3×10^{10}	8
TXS rotating anode	180	2×10^{11}	50

- Home lab sources intensities have significantly improved
 - Microfocus sources now up to 8 times the intensity of a classic 5 kW RAG
 - Latest microfocus RAGs up to 50 times the intensity of a classical 5 kW RAG
- However, beamlines still more than two orders of magnitude brighter!
- *Is it possible to produce a true synchrotron-class source in the home lab?*

NEW: Liquid metal sources

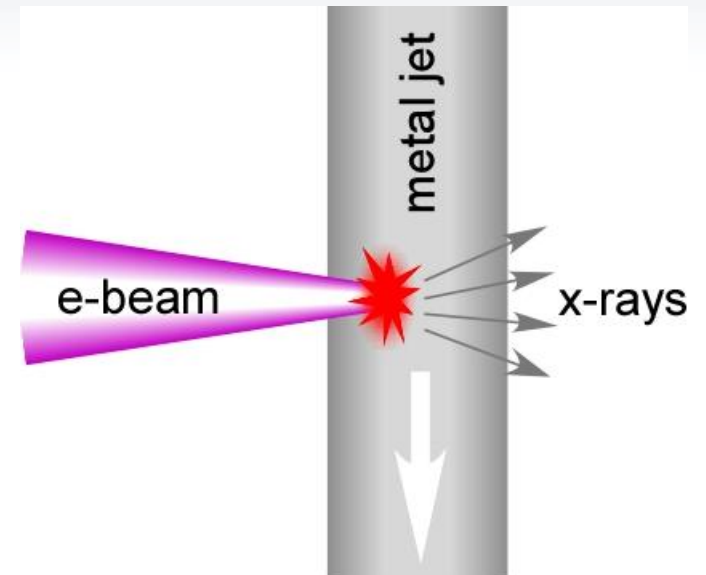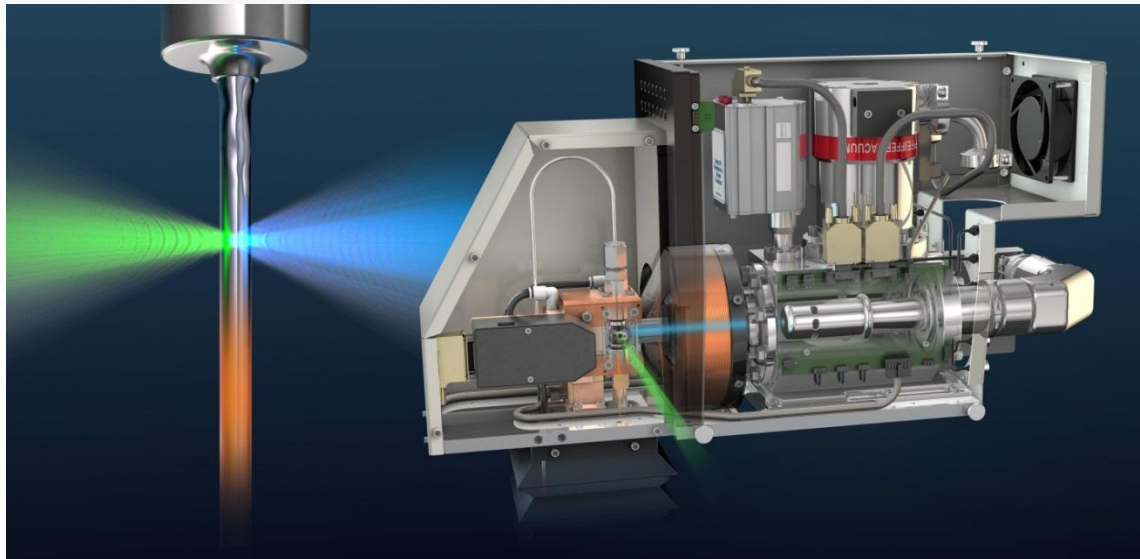


- High-speed liquid-metal-jet anode
- Anode is regenerative
- No longer limited by melting
- $>500 \text{ kW/mm}^2$ e-beam power density
 - Rotating anode limited to maximum 50 kW/mm^2

MetalJet Source Details

The path of the continuously recycled liquid metal alloy.

X-rays are emitted from the interaction point between the metal jet and the e-beam.

Electronics is housed in two 19" boxes.

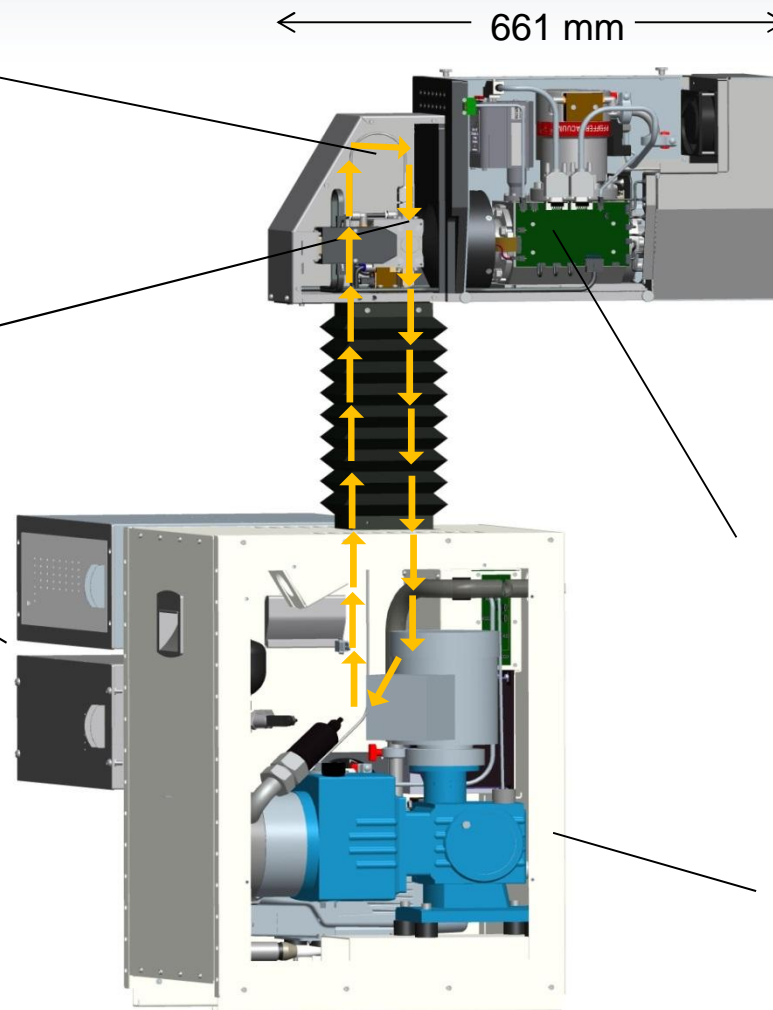


Advanced electromagnetic focusing optics together with a high brightness LaB6 cathode results in a very high quality e-beam focus.

Pumps etc. are housed in a 19" box.

Spot Size



5 μm


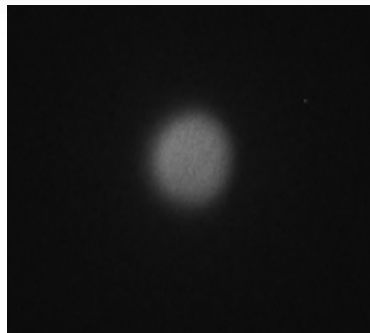


10 μm




20 μm

- Variable spot size: 5-20 μm

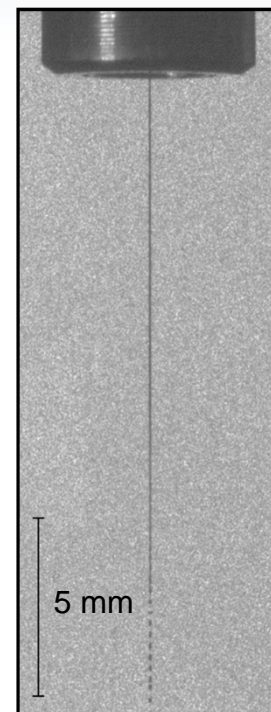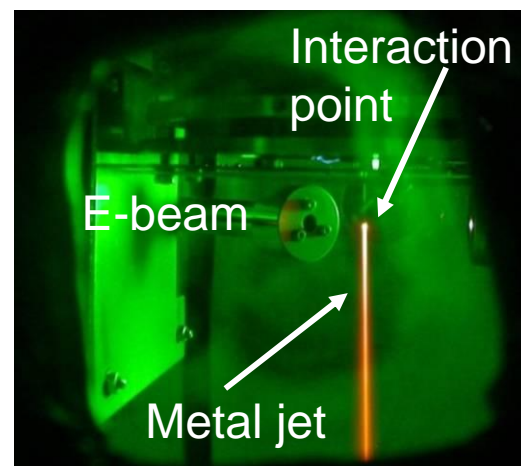

Source Spectrum and Brightness

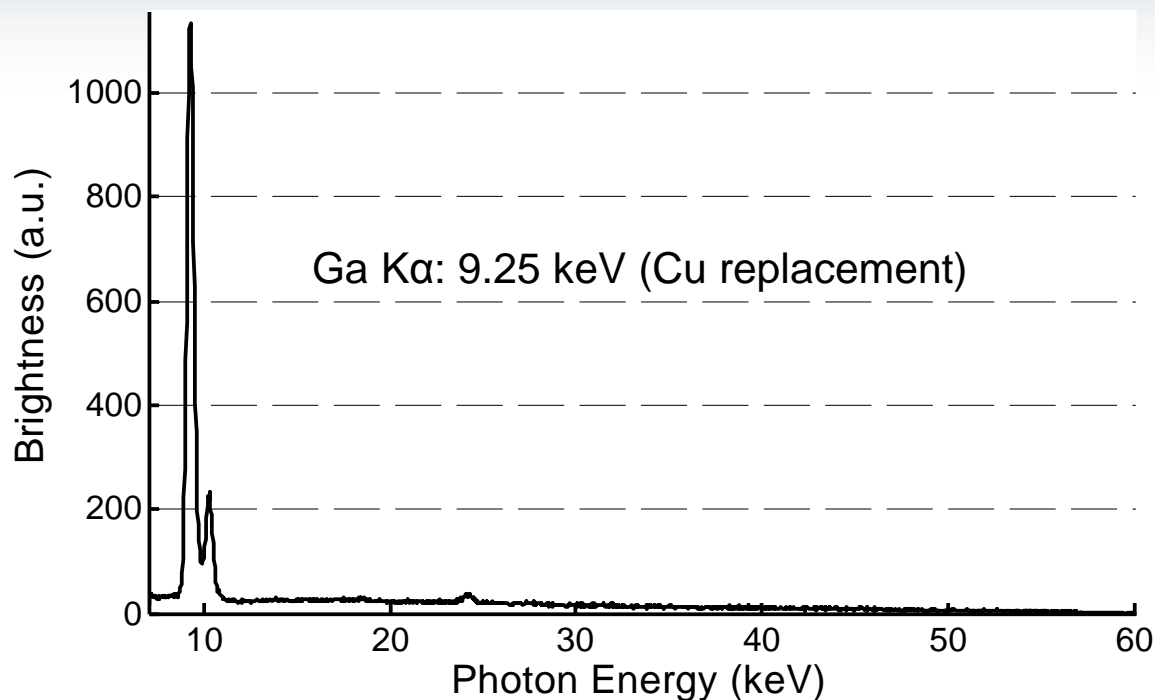


Spot size [μm , FWHM]	Voltage [kV]	Power [W]	Ga K α Brightness [Photons/(s \times mm ² \times mrad ² \times line)]
5	60	50	1.5×10^{11}
10	60	100	7.6×10^{10}
20	60	200	3.8×10^{10}

NANOSTAR with MetalJet




- Extraordinary flux, *comparable to synchrotron beamlines*
 - Up to 5×10^9 cps
- q range 0.005-0.45 \AA^{-1}
- High flexibility
 - Isotropic or anisotropic samples
- Easy maintenance
 - Comparable to standard rotating anode
- Optional WAXS attachment (IP)
- Rat tail tendon

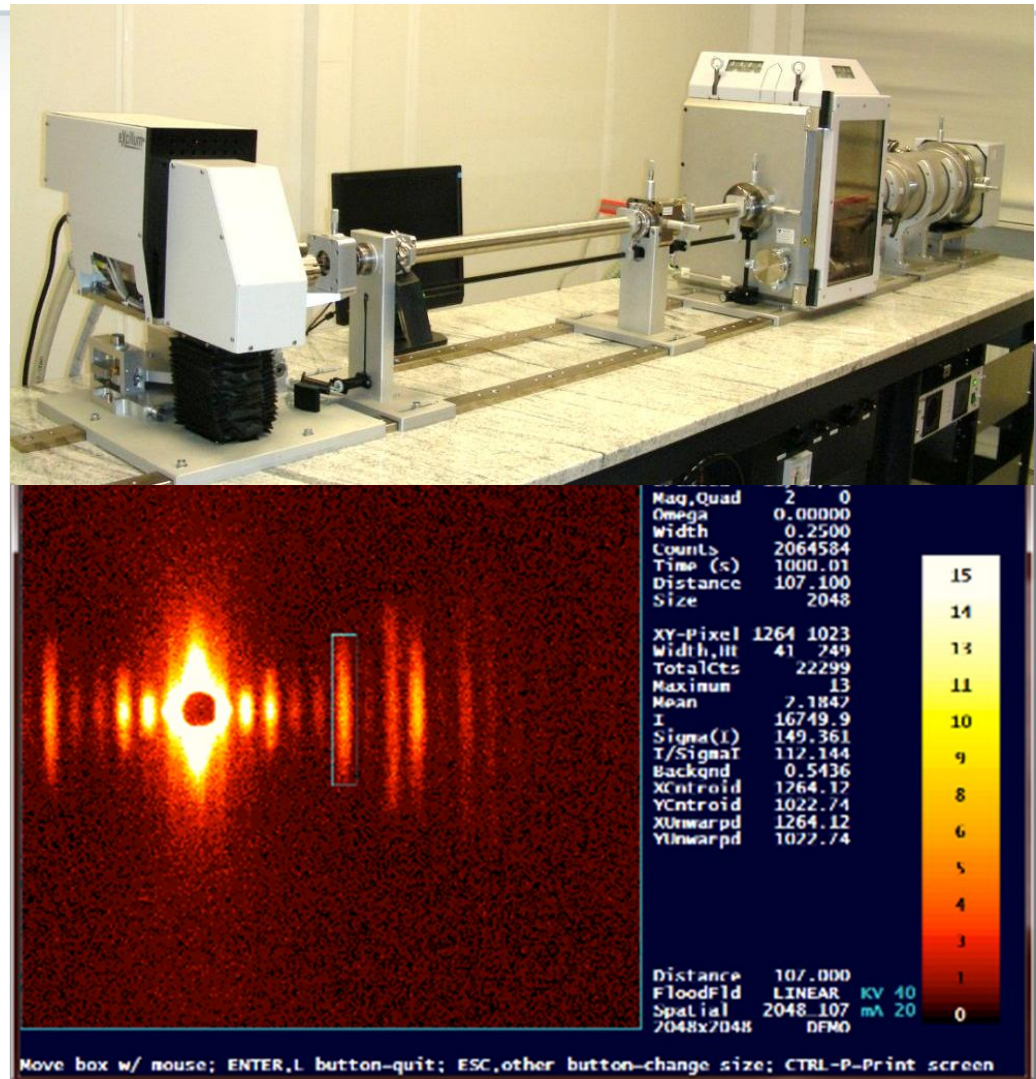

Flux comparison from glassy carbon



	Microfocus Cu-1 μ S	Rotating anode HB-TXS	MetalJet JXS-D1
Glassy Carbon			
Flux (cps)	1,50E+05	7,00E+05	4,00E+06

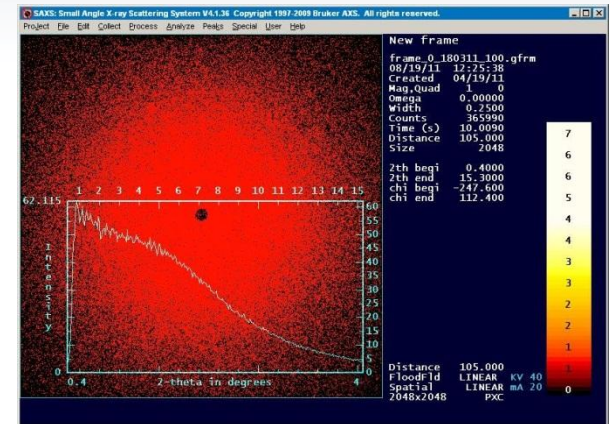


- Signal from glassy carbon 6 times higher than the brightest available rotating anode (HB-TXS)
 - **>140 times** higher intensity than classical rotating anode (5 kW)

So, is it possible to put a synchrotron beamline on a table top?



- Yes, at least the equivalent of a typical present generation bending magnet beamline


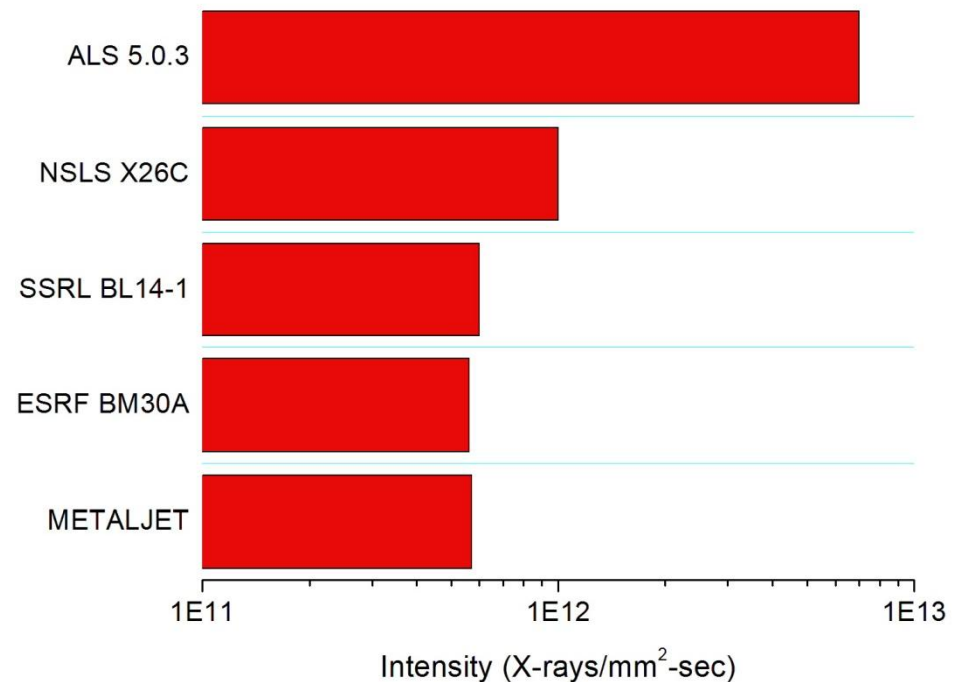


Optics and beam collimation

DAVINCI Innovations



DIFFRAC.SNAP-LOCK

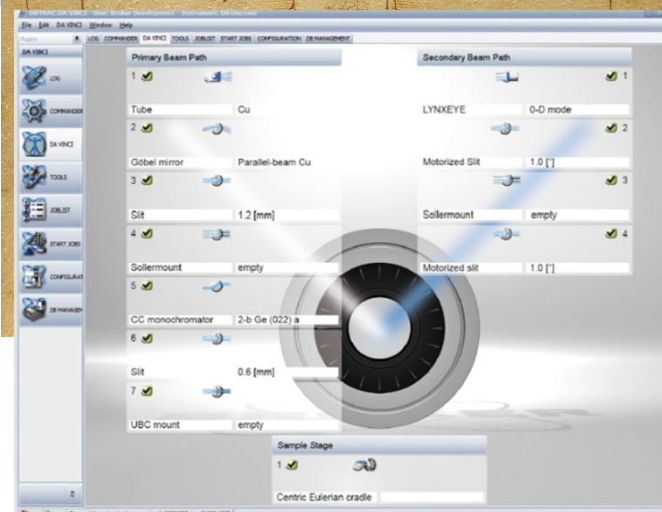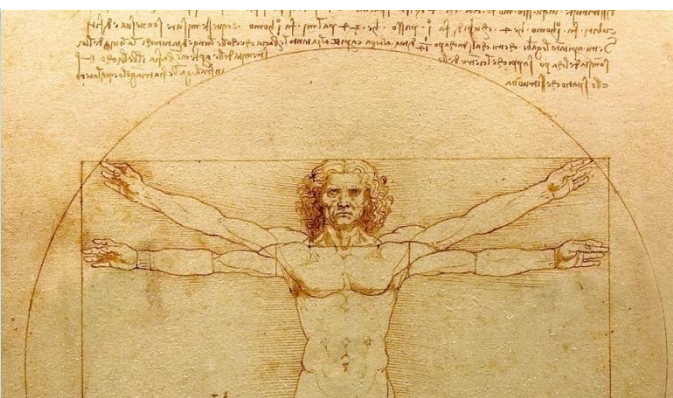


DIFFRAC."Da Vinci"
"The virtual goniometer"



DIFFRAC.MODE

No barrier between 0D/1D/2D


Vertical theta-theta, CEC for microdiffraction/stress/texture

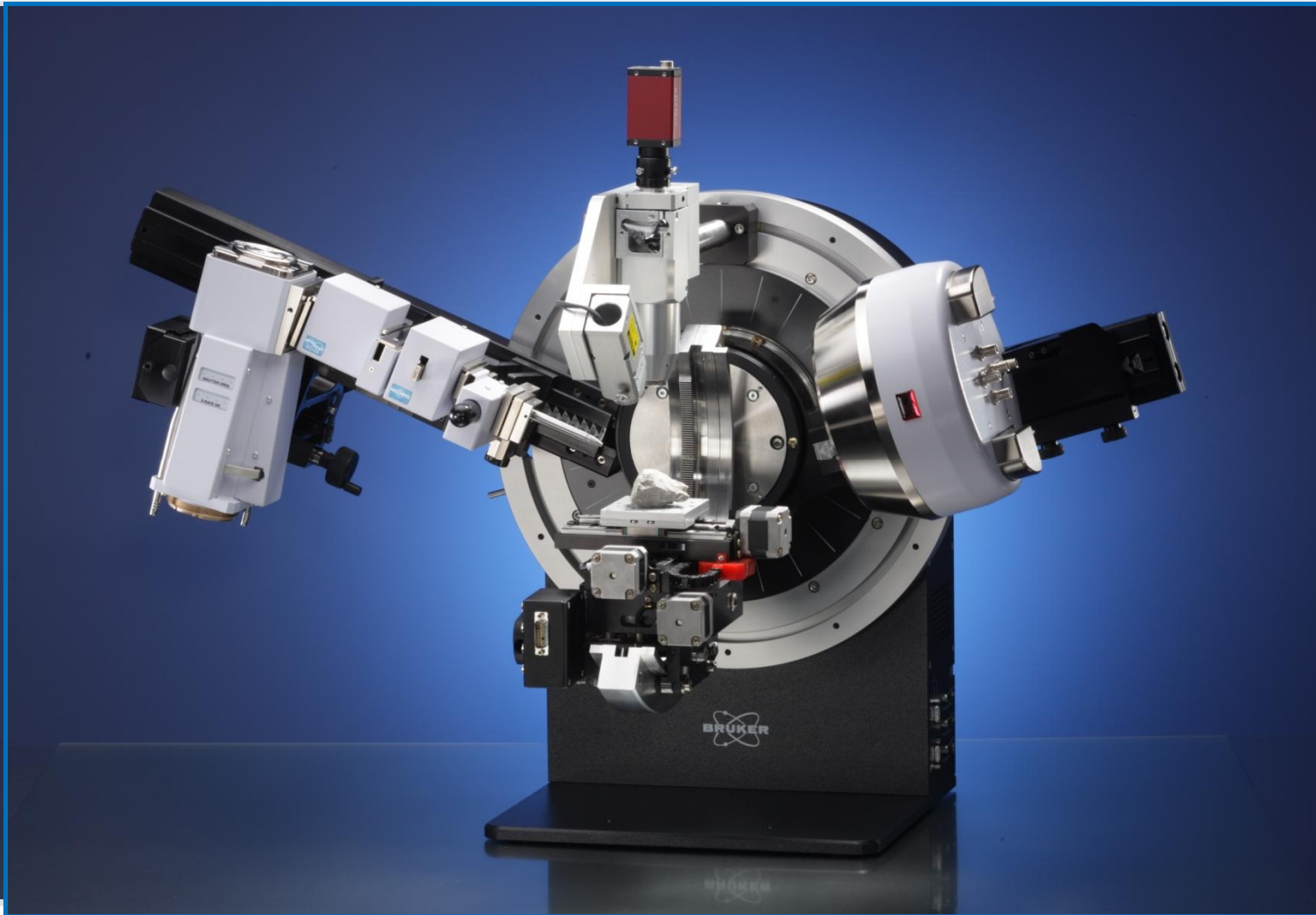


SAXS beam collimation



- Kratky camera
 - Advantages: Higher flux (for a given source), compact
 - Disadvantages: Half of scattered flux is lost, desmearing required.
- 3 pinhole
 - Advantages: uniform beam, no desmearing required, anisotropic samples
 - Disadvantages: lower flux, physically longer
- 2 pinhole
 - Apertures made of material that *does not scatter X-rays* (see Y. Li et al, App. Crys. 26, 2008, p1134)

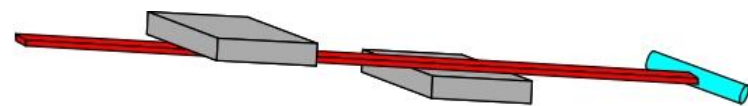

Kratky

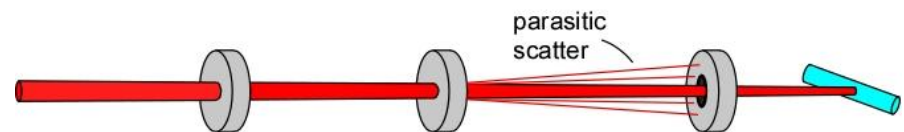


3 pinhole




Introducing the Bruker MICROpix, and MICROcalix



- Low cost
- High flux at sample
 - 1.4×10^8 cps
- High resolution
 - $q = 0.0056 - 0.45 \text{ \AA}^{-1}$
- Compact 2-D Kratky camera
 - Proven Hecus technology
- 50 W ImS source
 - Air-cooled
 - Low maintenance: 5 years tube life typical (3 years guaranteed)
- Air-cooled VANTEC detector
 - Quantum-limited sensitivity, small pixels, lowest noise
- SAXS/WAXS options

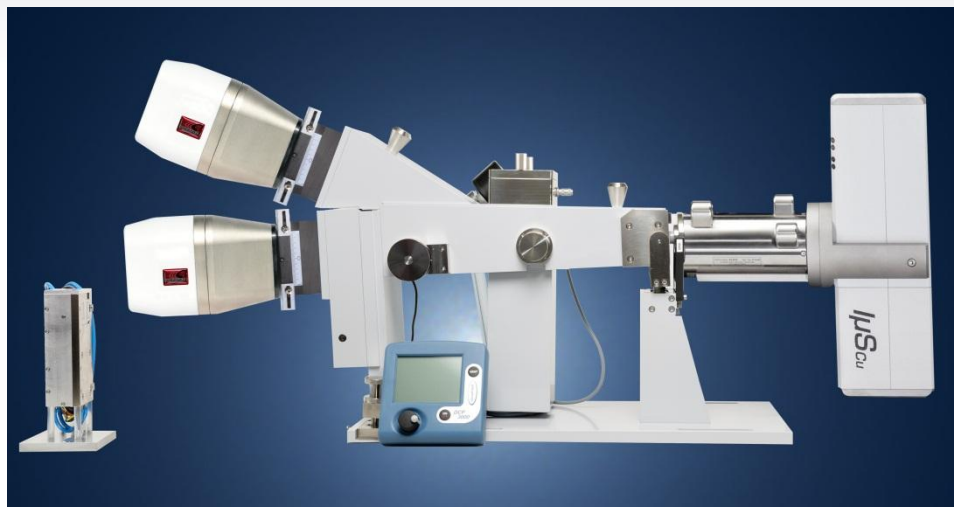

NANOSTAR

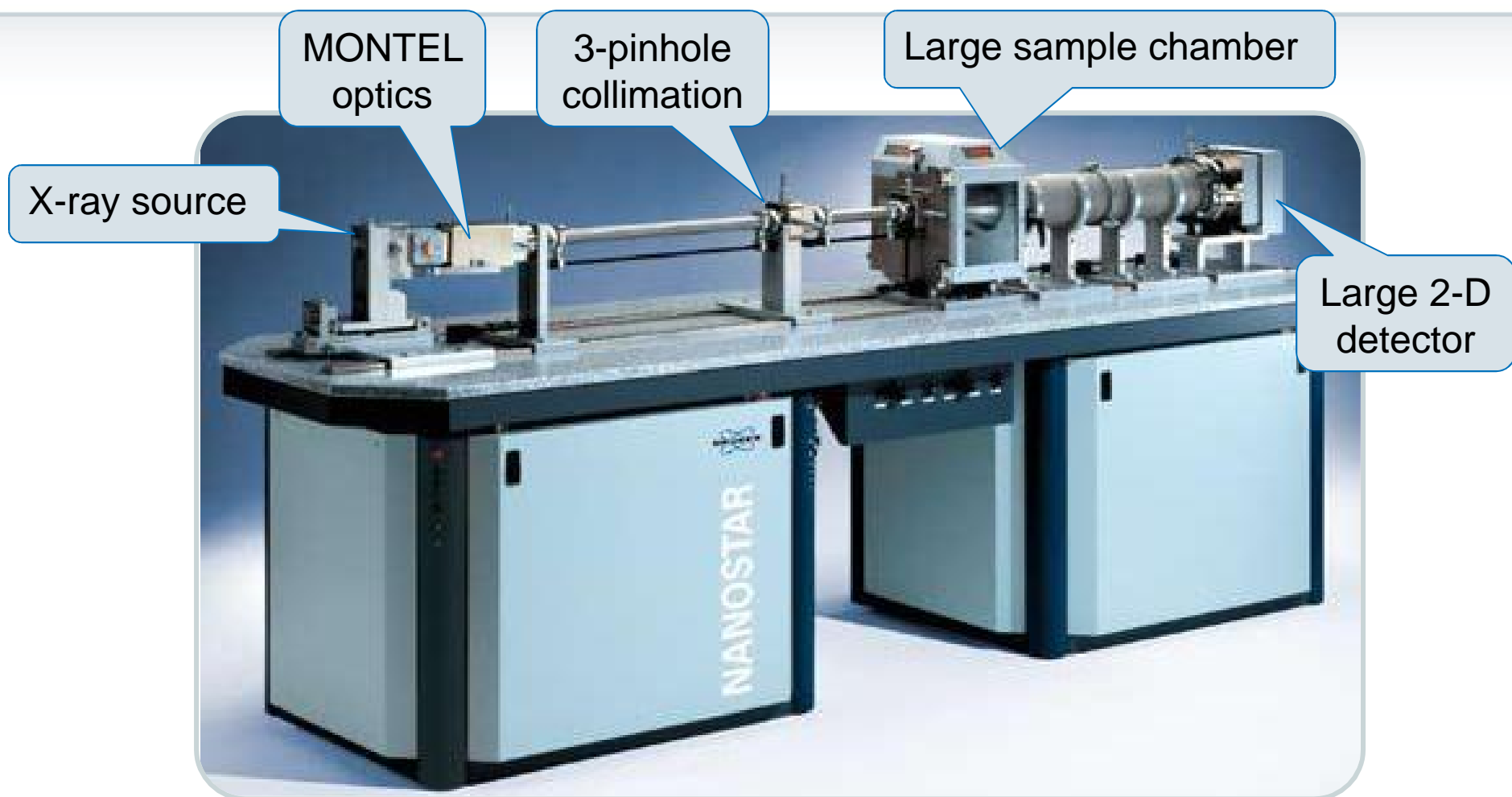

Detector


VANTEC-2000 - 2-D MIKROGAP™ X-Ray Detector




US Patent 6,340,819

- 140mm x 140mm detector window
 - Covering large Q-range
- True photon counter
 - Real time data collection and display
- No intrinsic detector noise
 - High sensitivity and low background
- High Local and Global Count Rate
- Extended Dynamic Range
 - $> 10^8$ (local max. count rate / local noise rate)
- High Spatial Resolution
 - $68 \mu\text{m} \times 68 \mu\text{m}$ pixel size
- Proven radiation-hardness
- No maintenance required

2-D SAXS/WAXS Typical Data

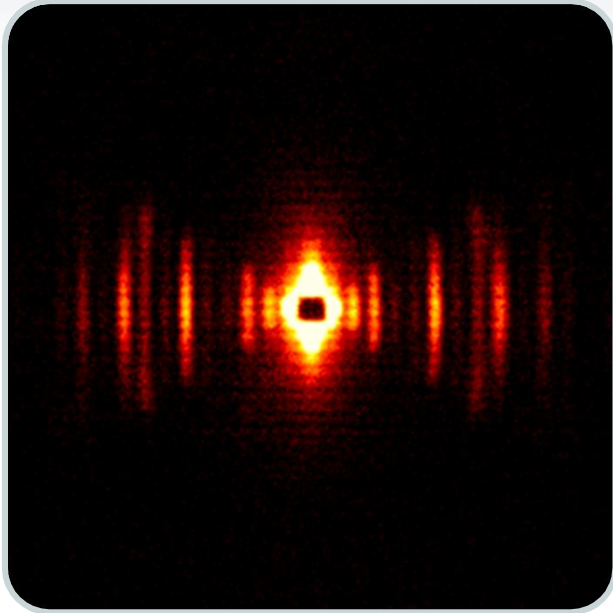

2-D SAXS pattern of collagen fibrils

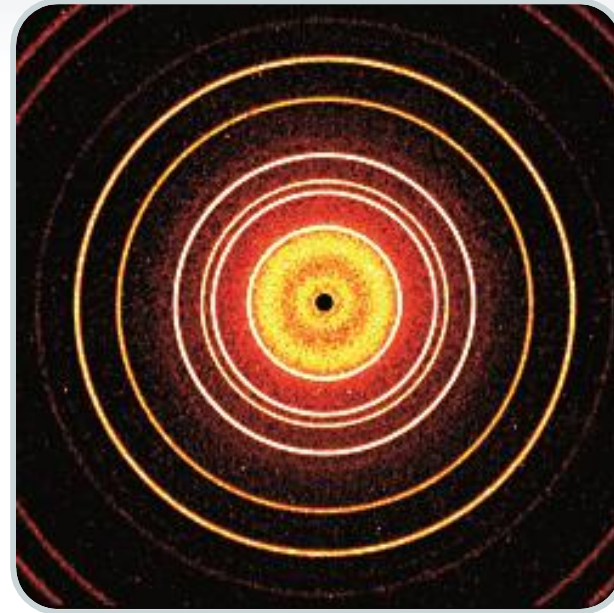

2-D WAXS pattern of Al_2O_3

- A 2-D detector is essential for measuring anisotropic data, but also enhances sensitivity for isotropic data
- Data evaluation is done on integrated 1-D data

VÅNTEC-500 – Outperforms all previous gaseous detectors.




Detector geometry:

- Be-window opening 140 mm in dia.
- Frame size:
 - 2048 x 2048 pixels
 - 1024 x 1024 pixels
 - 512 x 512 pixels
- Pixel size:
 - 68 μm x 68 μm
 - 136 μm x 136 μm
 - 272 μm x 272 μm
- Detector working distance:
5~30 cm in D8 DISCOVER enclosure
- 2θ range in a single frame:


5 cm	83°
10 cm	56°
15 cm	42°
20 cm	33°
25 cm	27°
30 cm	23°

XRD²: Single Frame Covering All

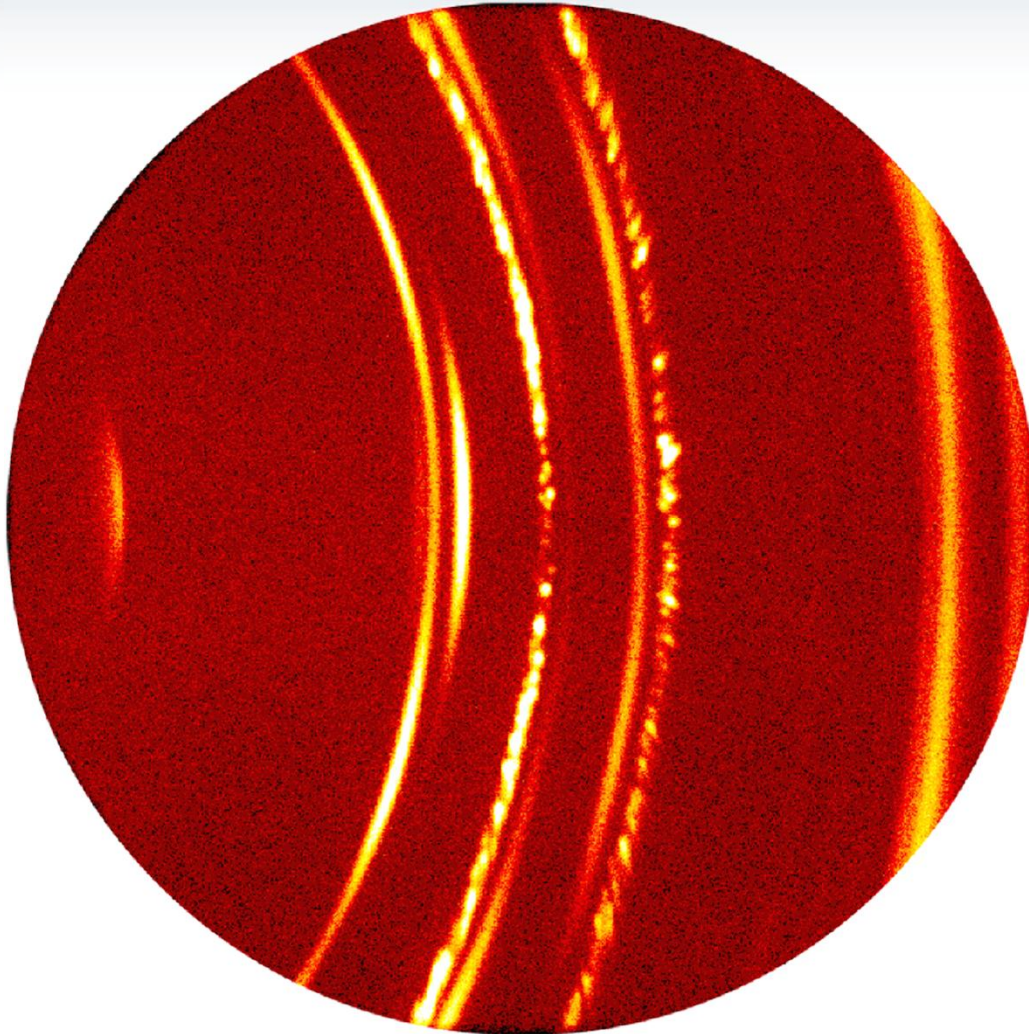


- 2θ coverage: 70° at 8 cm detector distance
- In-situ measurement for chemical reaction, phase transformation or other real-time physical changes.
- Sample with strong texture and large grain

XRD²: Frame Merge and Integration

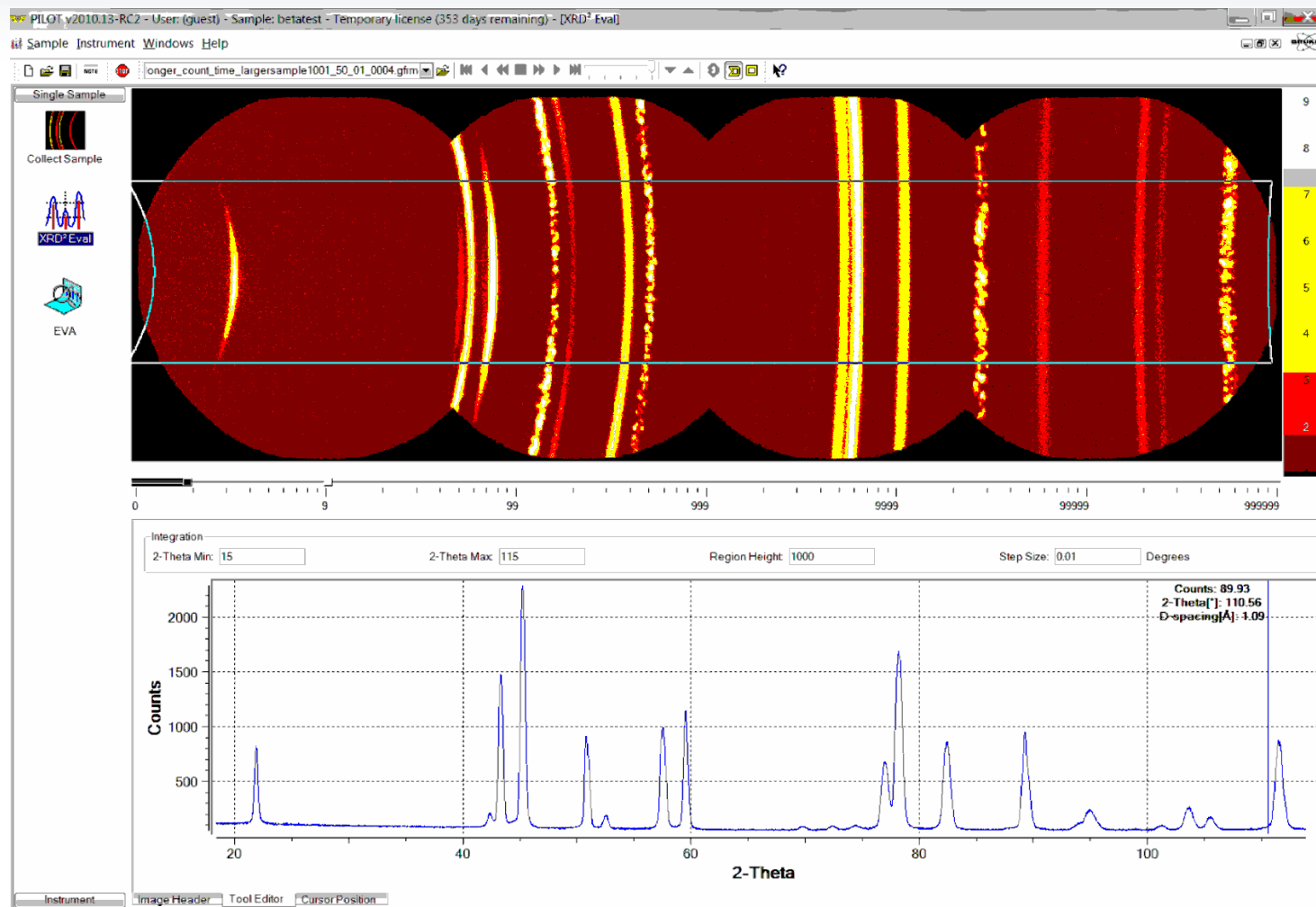


- 4 frames at 20 cm
- Merged frames 20 coverage: 100°
- Integrated profile for phase ID search/match

The D8 DISCOVER with DAVINCI VÅNTEC-500 for texture measurement

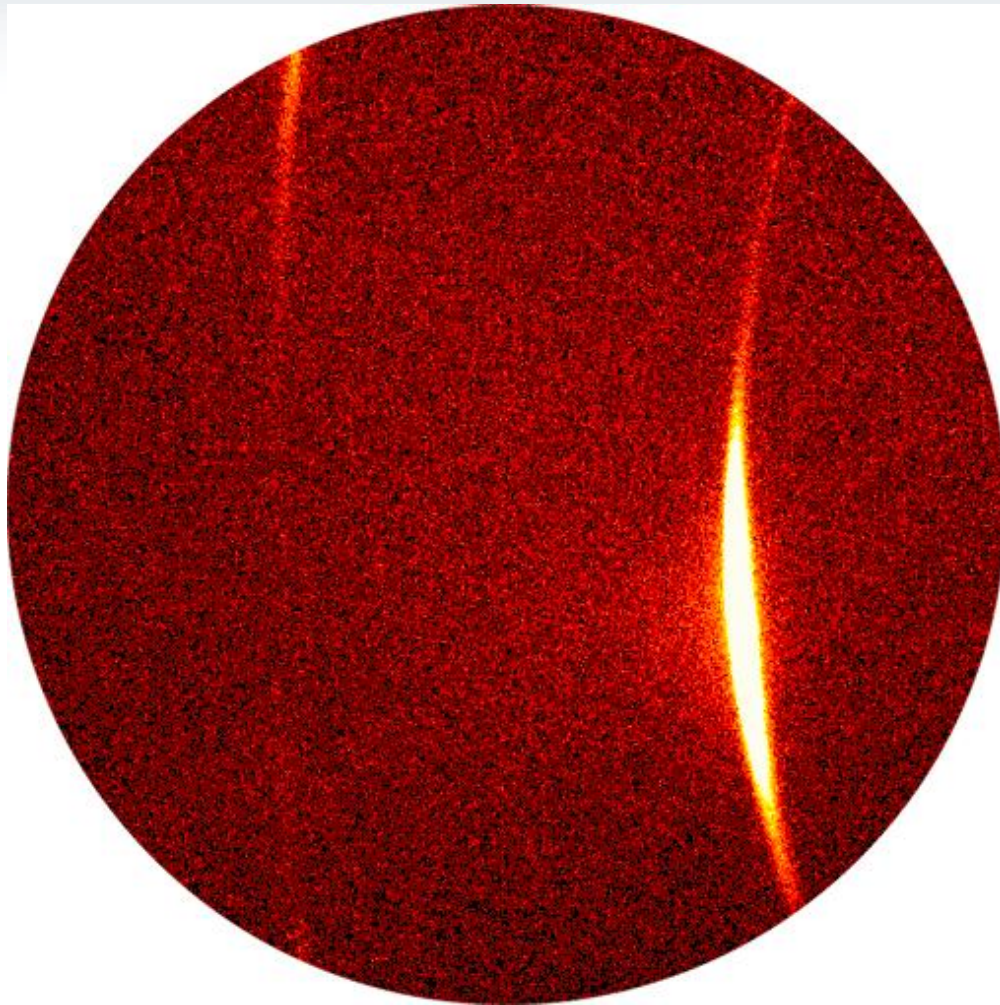

- Steel can
- (200) & (110) rings
- Intensity variation during ϕ scan

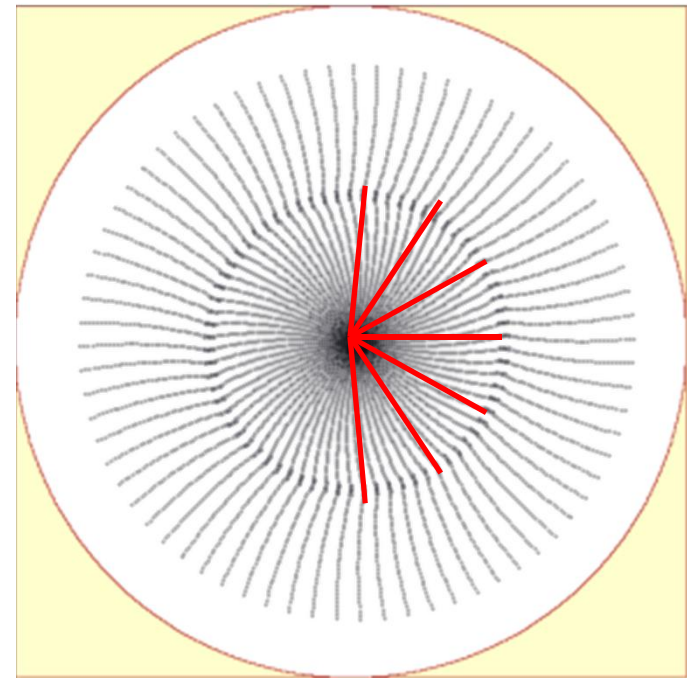

XRD²: Crystal Size by γ profile analysis: Organic glass for food & drugs




Food

Bio-preservation



In sugar glasses


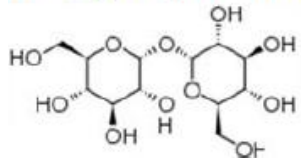


In amber





Annealing of GC crystals

LM

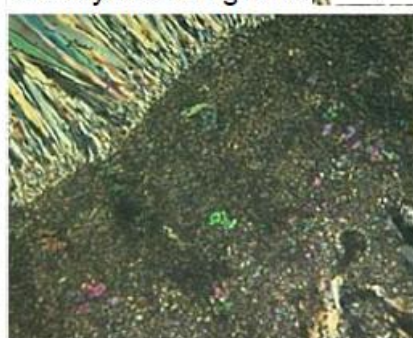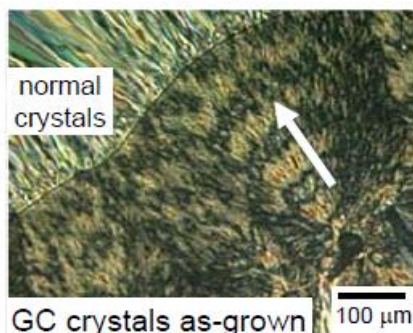


50 min/313 K

AFM

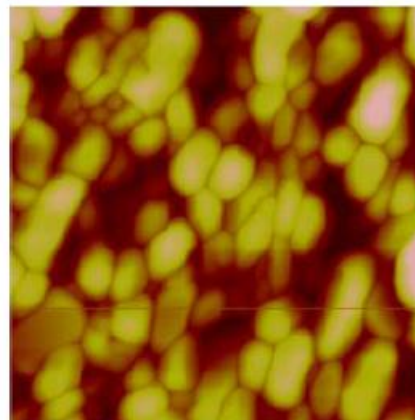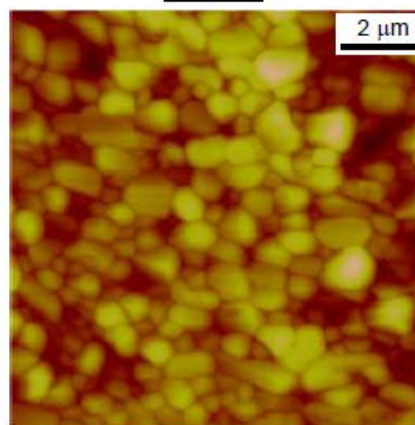


30 min/313 K


2D XRD

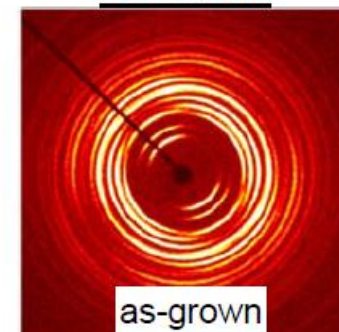

as-grown

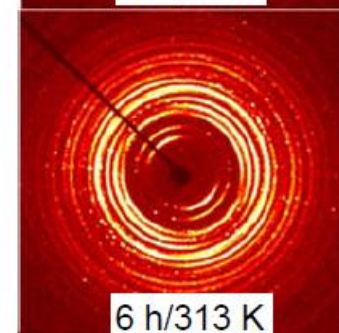


6 h/313 K

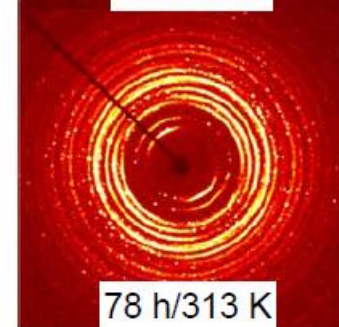

78 h/313 K


*Courtesy of Prof. Lian Yu, U. of Wisconsin - Madison

XRD²: Crystal Size by γ profile analysis: Acetaminophen powder

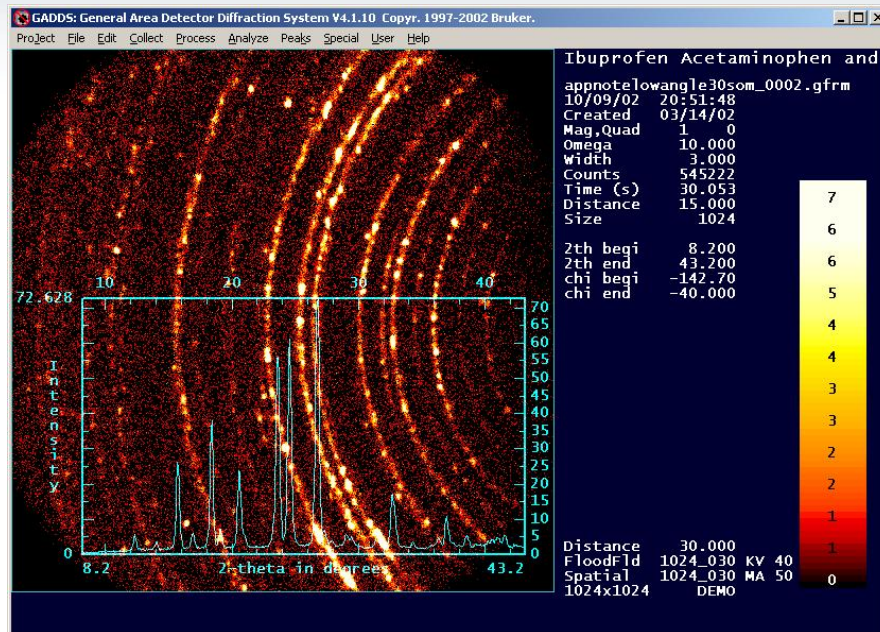


- The spotty diffraction ring is due to the large crystallites compared to the sampling volume (beam size).
- The number of spots on the ring is determined by crystallite size, instrumental window (γ -range), multiplicity of the crystal plane, and effective diffraction volume.
- The size of jelly beans and candy bin determines how many you can fill.

XRD²: Particle Size Analysis

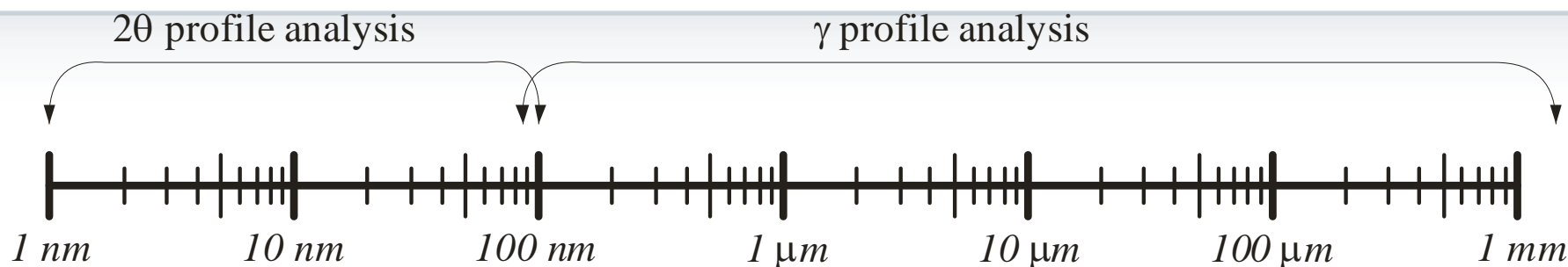


- 2θ profile analysis, suitable for <100 nm. Scherrer equation:

where B is FWHM corrected for instrument broadening,
C is 0.9~1 (crystal shape).

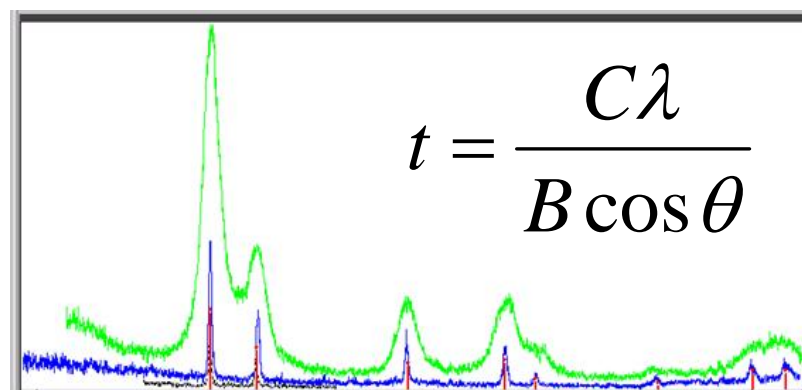

Gold nano-particle

- γ profile analysis is suitable for particle size from sub-micrometer to a few millimeters.

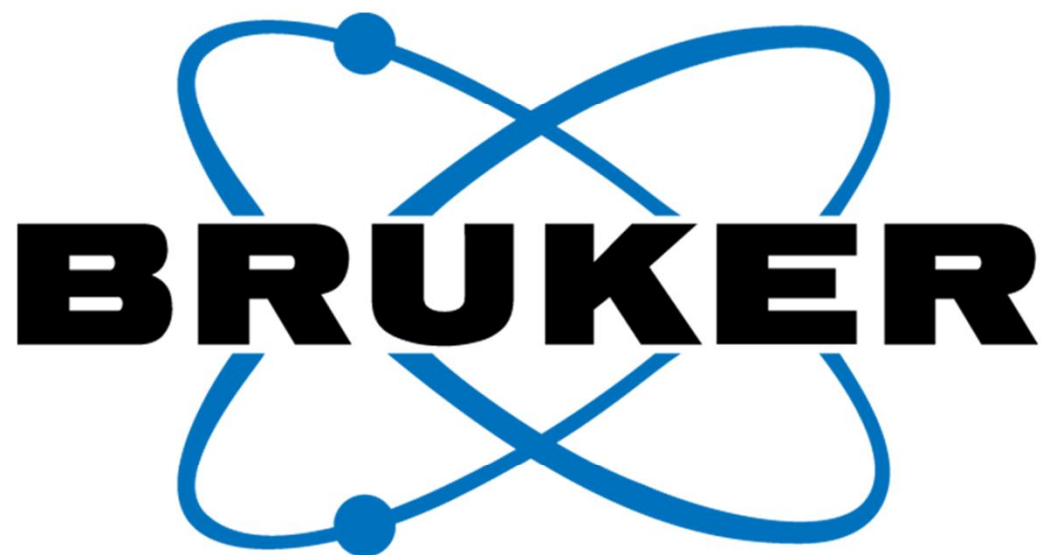

www.bruker.com

Thanks for your attention