


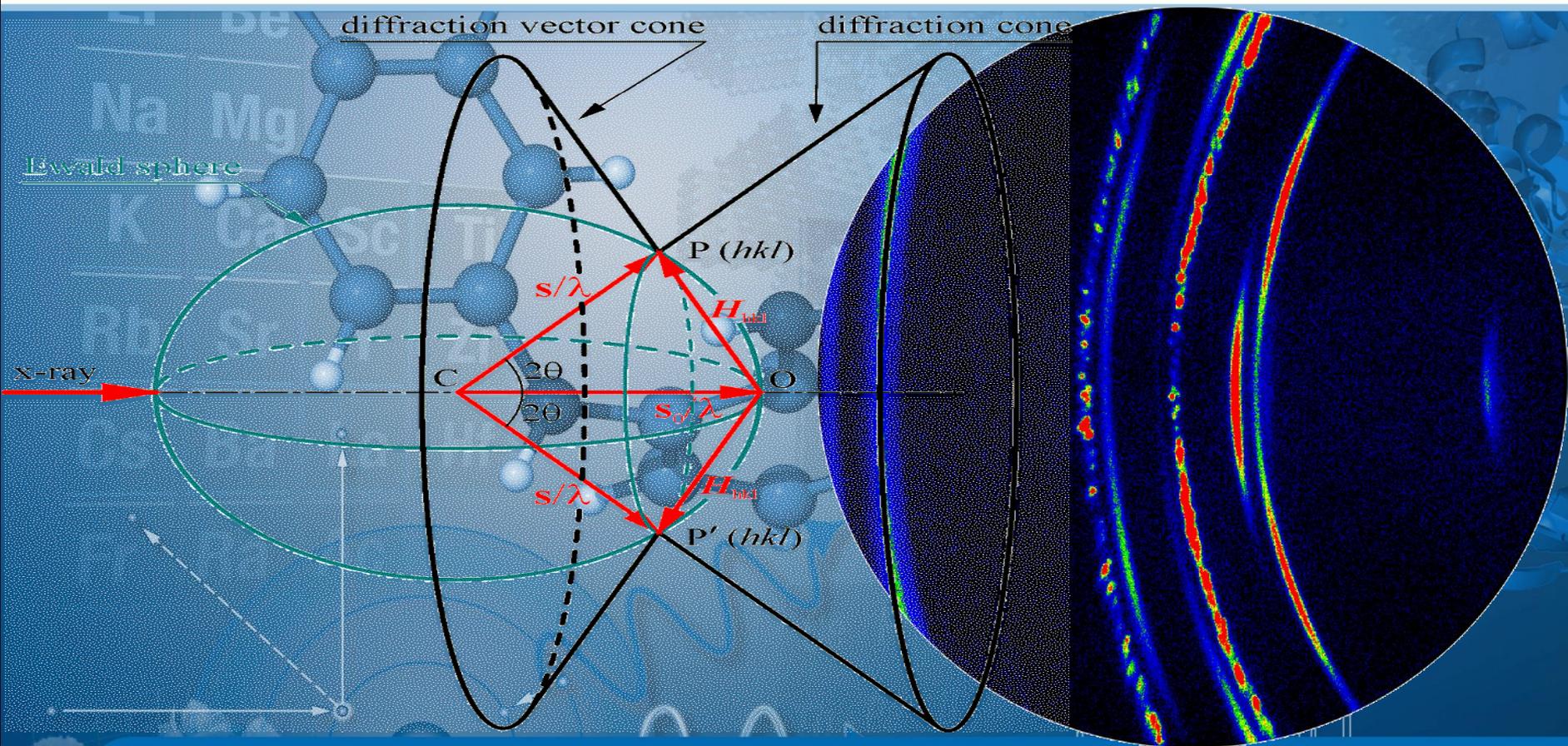

Instrumentation and Applications of XRD² for Pharmaceuticals

Bob He, Bruker AXS



This document was presented at PPXRD - Pharmaceutical Powder X-ray Diffraction Symposium

Sponsored by The International Centre for Diffraction Data

This presentation is provided by the International Centre for Diffraction Data in cooperation with the authors and presenters of the PPXRD symposia for the express purpose of educating the scientific community.

All copyrights for the presentation are retained by the original authors.

The ICDD has received permission from the authors to post this material on our website and make the material available for viewing. Usage is restricted for the purposes of education and scientific research.



PPXRD Website – www.icdd.com/ppxrd

ICDD Website - www.icdd.com

X-ray Applications for typical pharmaceutical samples



XRD & XRD ²	Single Crystal	Several Grains	Powder	Finished Product	Solutions
Qualitative Phase ID	✓⊕	✓⊕	✓Φ	✓Φ	✓Φ
Quantitative Rietveld analysis			✓		
Quantitative analysis with standards		✓	✓	✓	
X-ray movie, Non-Ambient	✓Φ	✓Φ	✓Φ	✓Φ	✓Φ
Structure solution, Indexing	✓Φ		✓		
Microdiffraction/ Mapping		✓⊕	✓⊕	✓⊕	
Shape analysis			✓Φ	✓Φ	✓Φ
HTS	✓⊕	✓⊕	✓⊕		
Grain-Size det.		✓Φ	✓Φ		
%Crystallinity		✓⊕	✓⊕	✓⊕	✓⊕


✓ - can be performed by either XRD or XRD²

Φ - better with XRD²


⊕ - accept performance and accurate results only with XRD²

Basic Concept – XRD²


Conventional X-ray Diffractometer

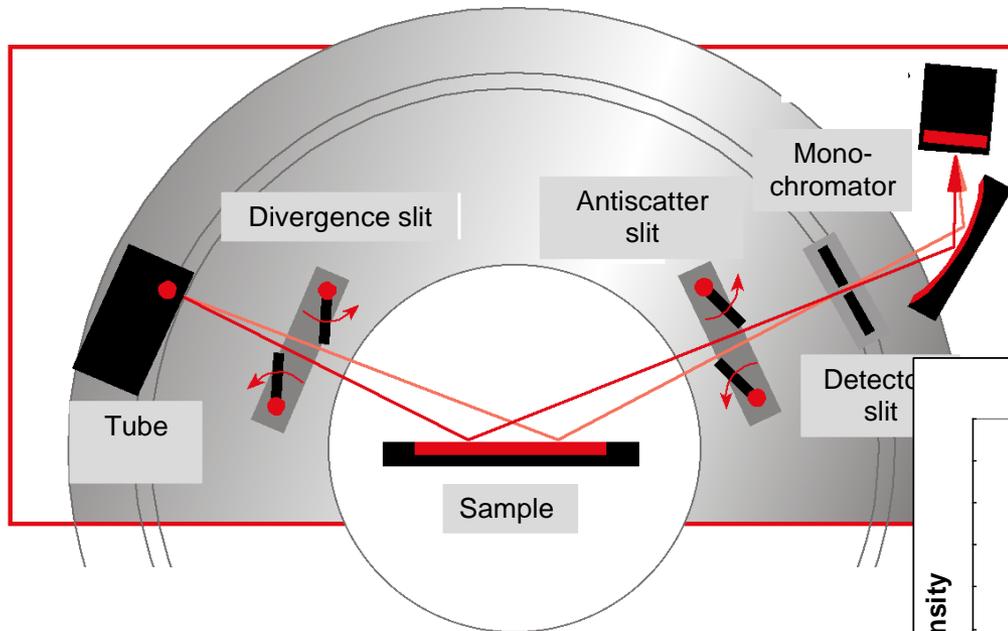



- Bragg-Brentano Geometry.
- Scanning over 2θ range to collect XRD pattern.

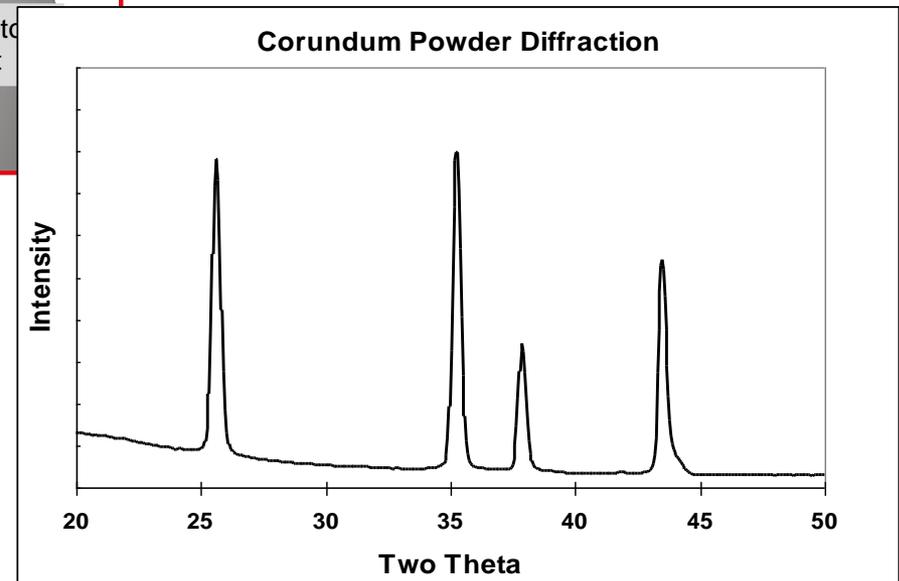

XRD²: Two-dimensional X-ray Diffraction

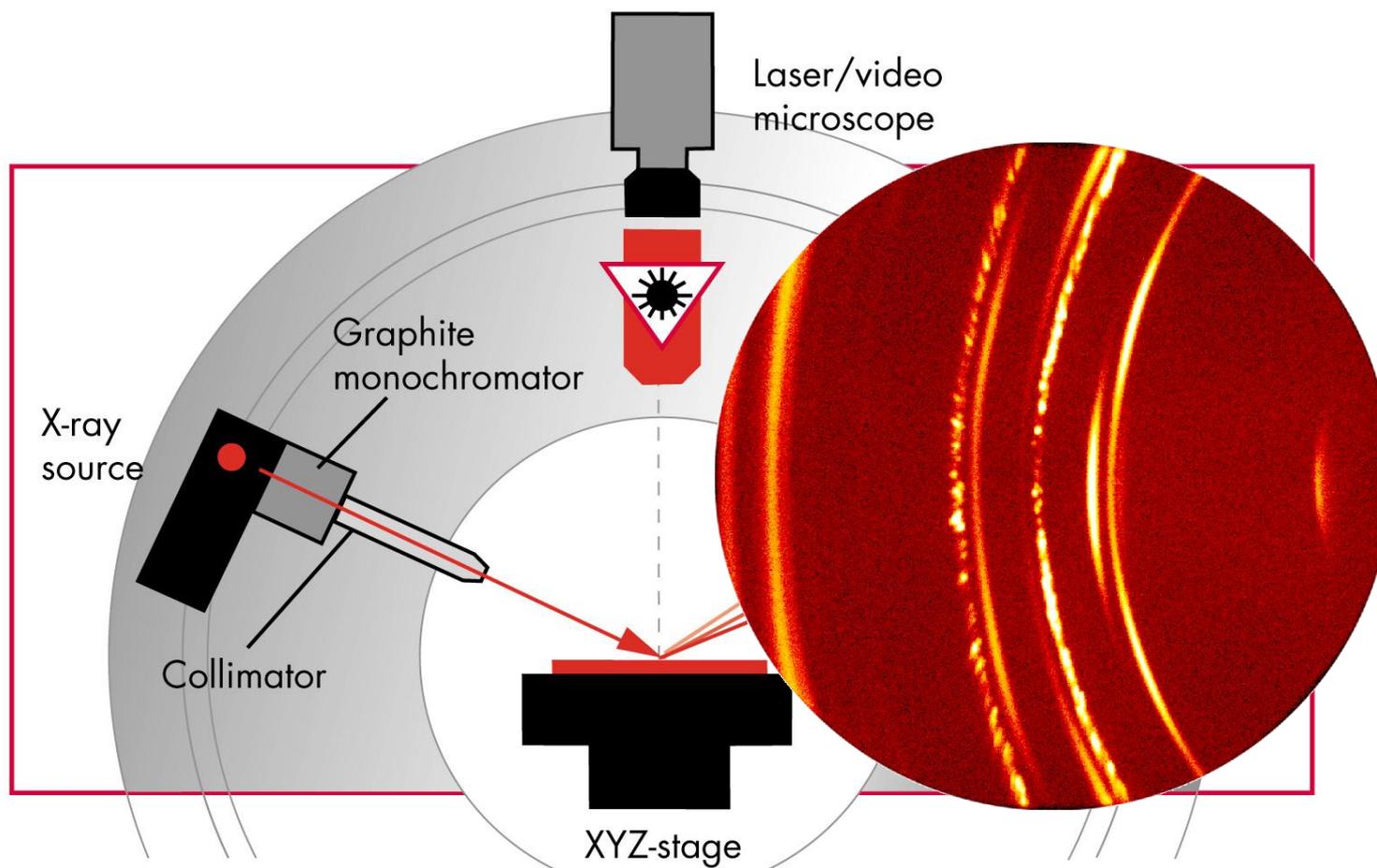

XRD²: Single Frame from Battery Anode Collected with Vantec-500 Detector

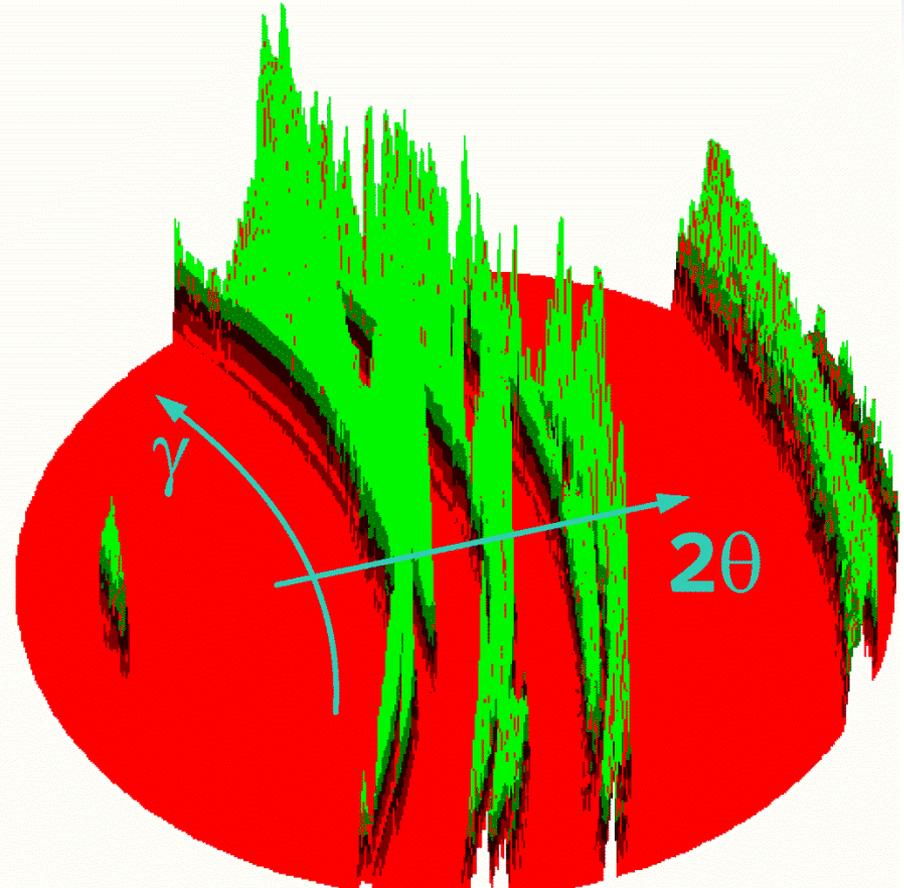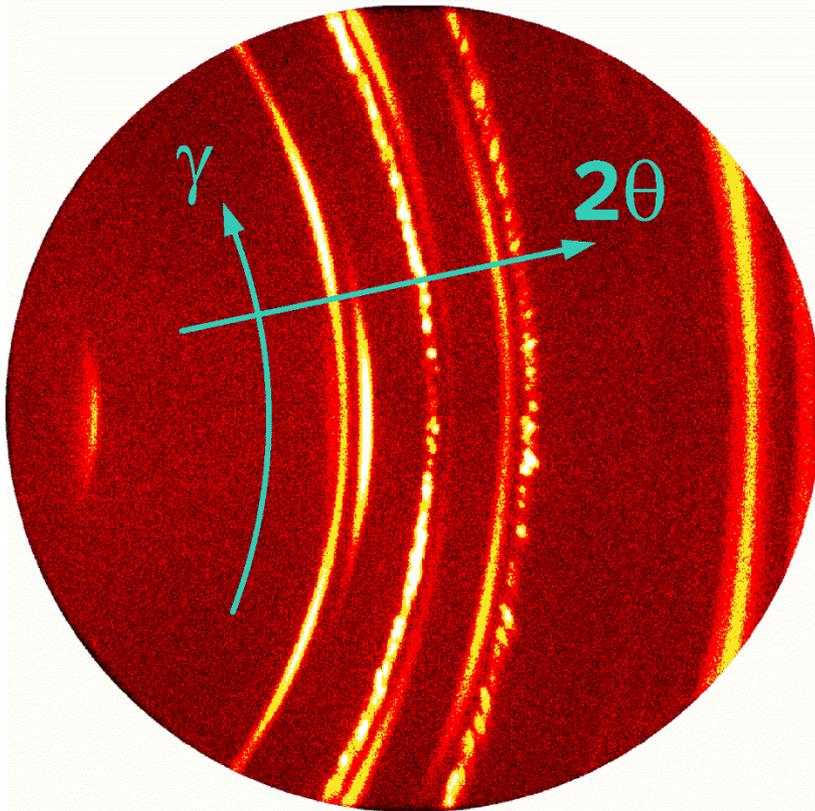



- 2θ coverage: 70° at 8 cm detector distance
- Contains information on phase, stress, texture and grain size

XRD²: 2D pattern in I on γ - 2θ coordinates



Powder

Texture

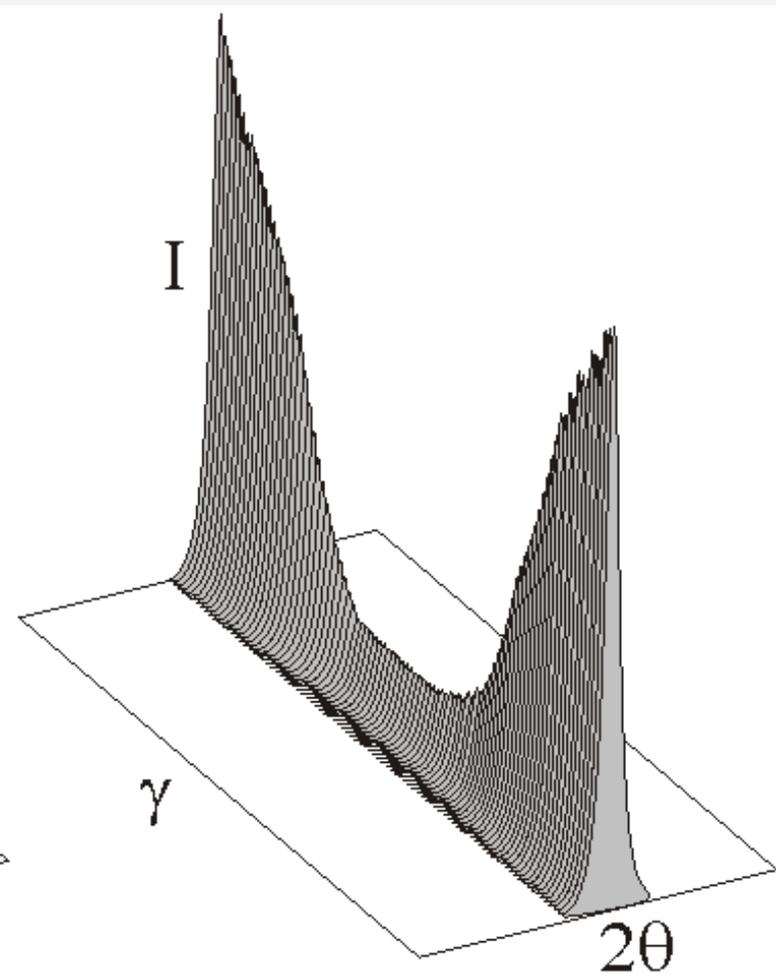
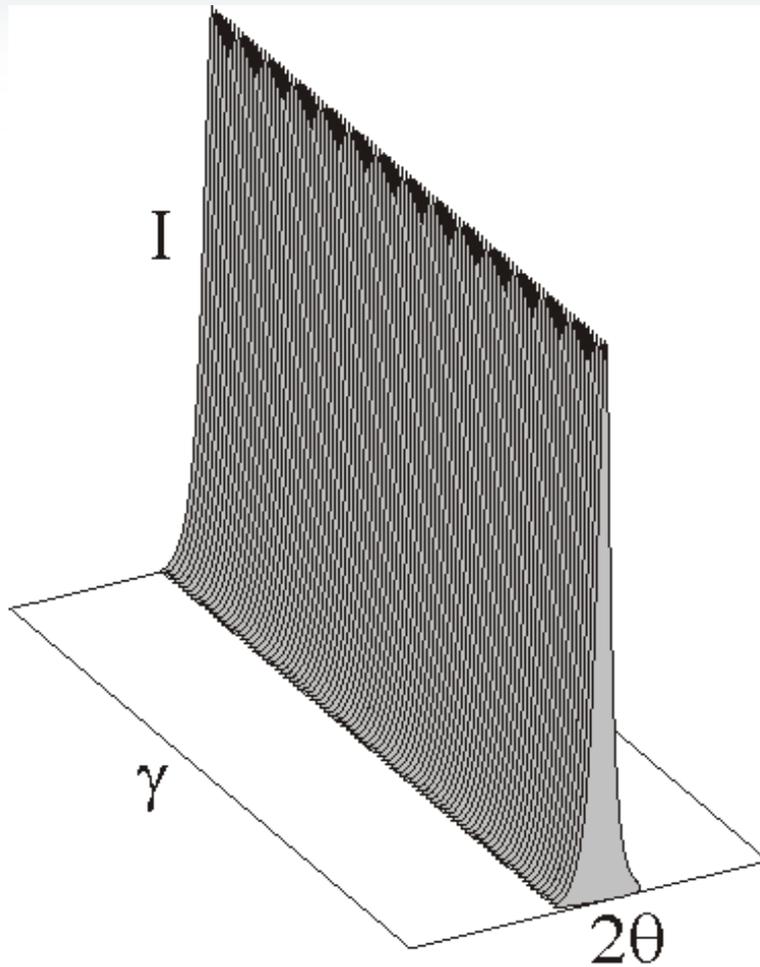


XRD²: 2D pattern in I on γ - 2θ coordinates



Stress

Large grains

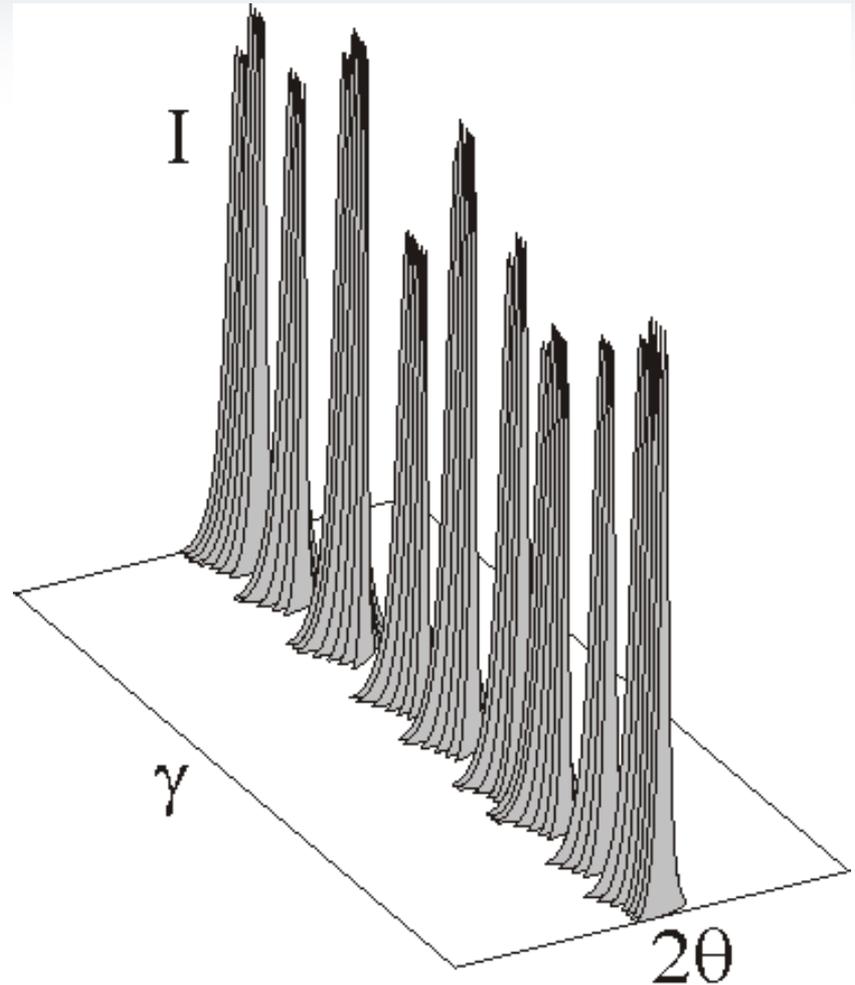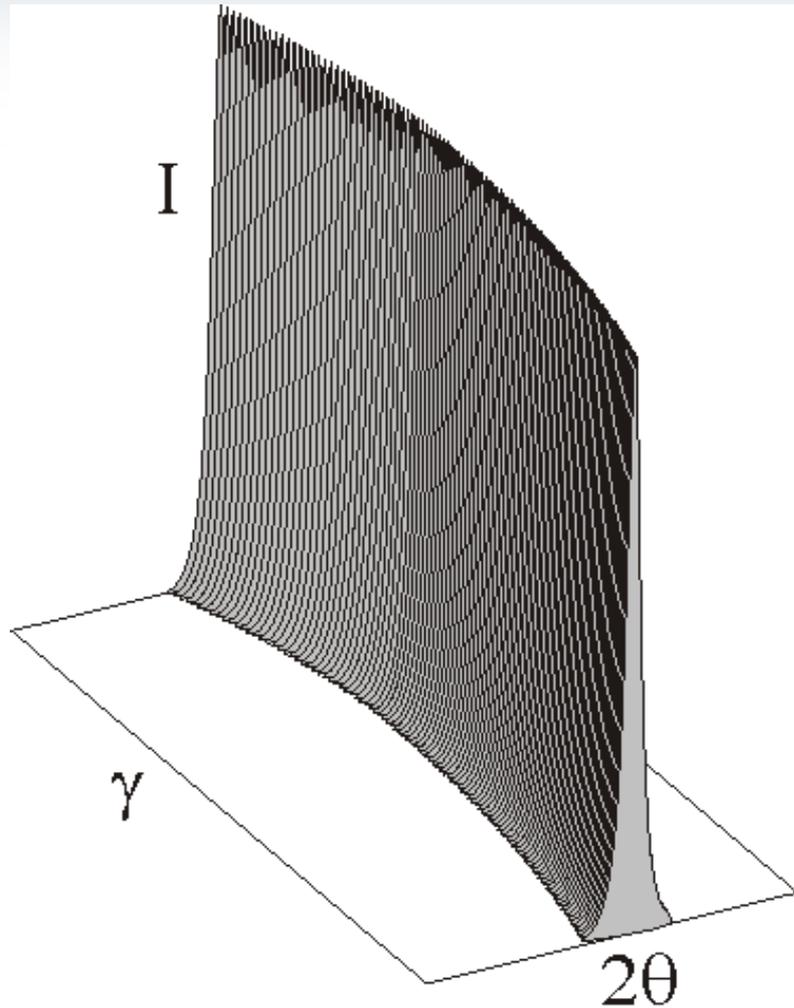

Fundamental Equations and Diffraction Vector Approach

XRD²: Diffraction Vector as a Function of 2θ and γ



Laue equations with 3 lattice axes:

$$\mathbf{a} \cdot (\mathbf{s} - \mathbf{s}_0) = h\lambda$$

$$\mathbf{b} \cdot (\mathbf{s} - \mathbf{s}_0) = k\lambda$$

$$\mathbf{c} \cdot (\mathbf{s} - \mathbf{s}_0) = l\lambda$$

The magnitude of the vector





$$\left| \frac{\mathbf{s} - \mathbf{s}_0}{\lambda} \right| = \frac{2 \sin \theta}{\lambda} = |\mathbf{H}_{hkl}| = \frac{1}{d_{hkl}}$$

Leads to Bragg law:

$$2d_{hkl} \sin \theta = \lambda$$


Diffraction condition in vector

$$\frac{\mathbf{s} - \mathbf{s}_0}{\lambda} = \mathbf{H}_{hkl}$$

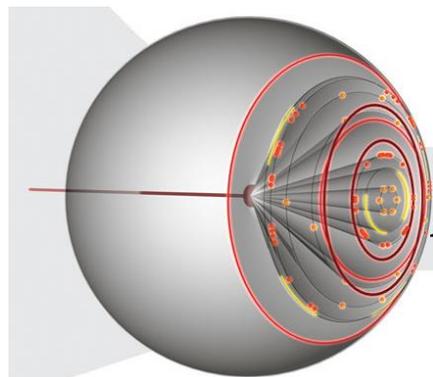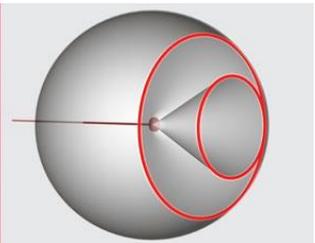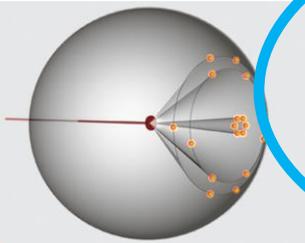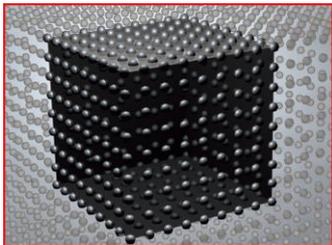

Introduce γ for XRD²:

$$\mathbf{H} = \frac{\mathbf{s} - \mathbf{s}_0}{\lambda} = \frac{1}{\lambda} \begin{bmatrix} \cos 2\theta - 1 \\ -\sin 2\theta \sin \gamma \\ -\sin 2\theta \cos \gamma \end{bmatrix}$$


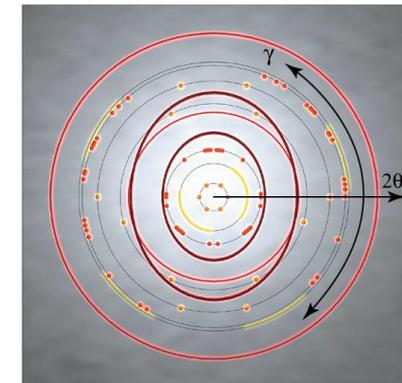


XRD²: Sample Space and Unit Diffraction Vector



The components of the unit vector \mathbf{h}_S in the sample coordinates $\mathbf{S}_1\mathbf{S}_2\mathbf{S}_3$ is then given by

$$\begin{bmatrix} h_1 \\ h_2 \\ h_3 \end{bmatrix} = \begin{bmatrix} a_{11} & a_{12} & a_{13} \\ a_{21} & a_{22} & a_{23} \\ a_{31} & a_{32} & a_{33} \end{bmatrix} \begin{bmatrix} h_x \\ h_y \\ h_z \end{bmatrix}$$

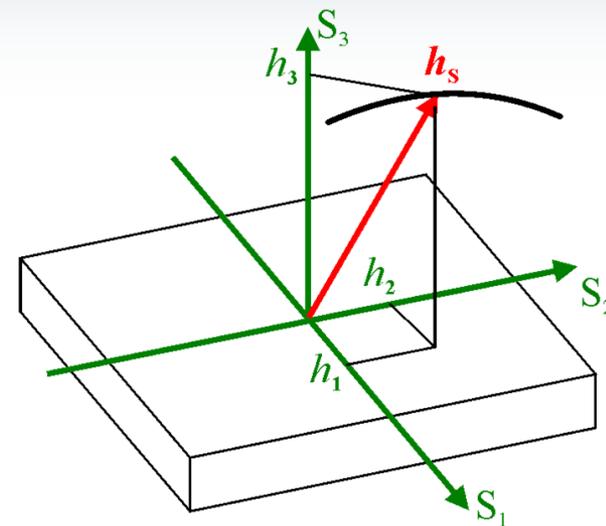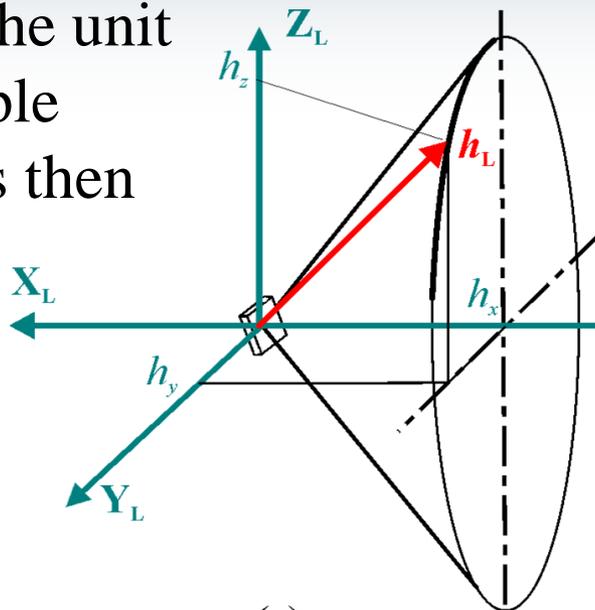

Or
in expanded form
for Eulerian
geometry:

$$\begin{aligned} h_1 &= \sin \theta (\sin \phi \sin \psi \sin \omega + \cos \phi \cos \omega) + \cos \theta \cos \gamma \sin \phi \cos \psi \\ &\quad - \cos \theta \sin \gamma (\sin \phi \sin \psi \cos \omega - \cos \phi \sin \omega) \\ h_2 &= -\sin \theta (\cos \phi \sin \psi \sin \omega - \sin \phi \cos \omega) - \cos \theta \cos \gamma \cos \phi \cos \psi \\ &\quad + \cos \theta \sin \gamma (\cos \phi \sin \psi \cos \omega + \sin \phi \sin \omega) \\ h_3 &= \sin \theta \cos \psi \sin \omega - \cos \theta \sin \gamma \cos \psi \cos \omega - \cos \theta \cos \gamma \sin \psi \end{aligned}$$

XRD²: Diffraction Vector as a Function of γ



Stress and strain:



Texture (pole-figure):

$$\alpha = \sin^{-1} |h_3|, \quad \beta = \pm \cos^{-1} \frac{h_1}{\sqrt{h_1^2 + h_2^2}}$$

$$\epsilon_{(\gamma, \omega, \psi, \phi)} = \epsilon_{ij} \cdot h_i \cdot h_j$$



Unit vector:

$$\mathbf{h}_L = \begin{bmatrix} h_x \\ h_y \\ h_z \end{bmatrix} \quad \mathbf{h}_s = \begin{bmatrix} h_1 \\ h_2 \\ h_3 \end{bmatrix}$$

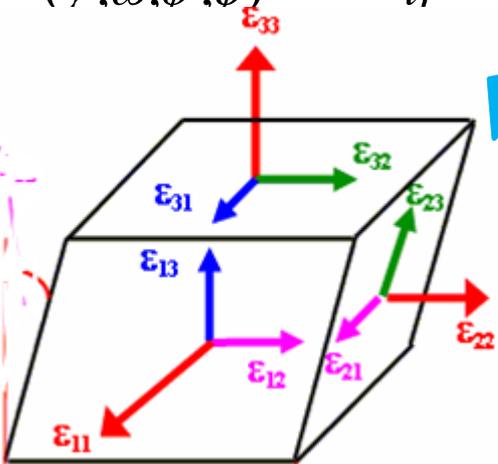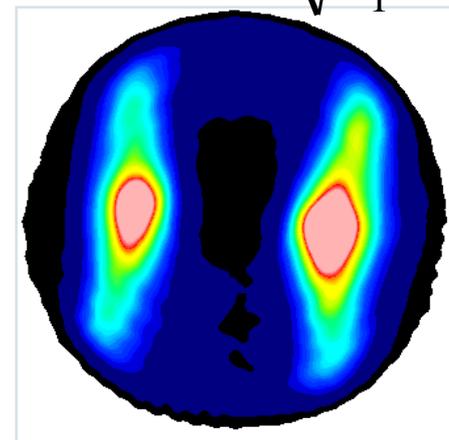

Crystal size:


$$N_s = p_{hkl} \cdot \frac{V f_i}{v_i} \cdot \frac{\Omega}{4\pi}$$

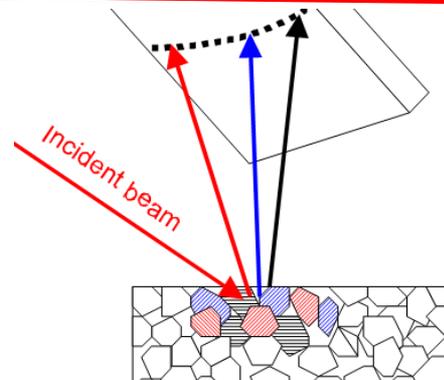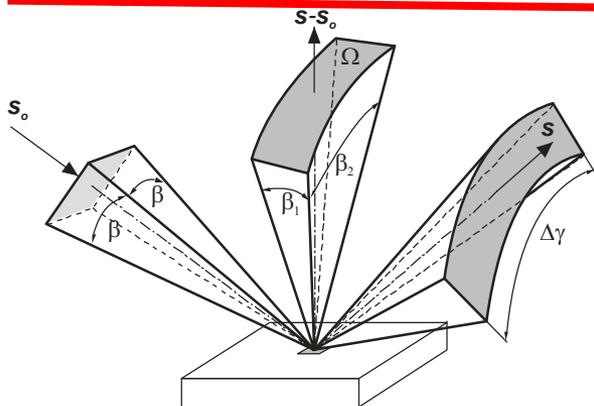

Sources & Optics

How to make brighter source I: Microfocus sources

- Brightness (**B**) is proportional to power loading (**p**)
- >99% power turns to heat and needs to be removed
- Power loading is higher for *smaller spot focus*

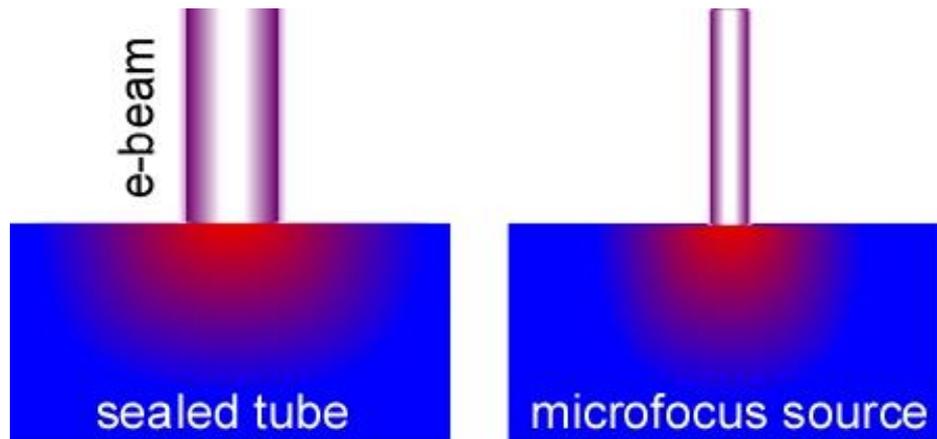

$$P_{\max} = \frac{2\kappa(T_m - T_0)}{r\sqrt{\pi \ln(2)}}$$

- | | |
|--|---|
| <ul style="list-style-type: none"> ■ Large spot ■ Quasi-one dimensional heat flow limits power loading | <ul style="list-style-type: none"> ■ Small spot ■ Two dimensional heat flow (more efficient cooling) ■ Relative performance improved by 10 times |
|--|---|



I μ S microfocus source



- Intensity 3×10^{10} X-rays/mm²-sec (Cu K α)
 - 8 times higher than conventional 5.4 kW rotating anode
- Typical lifetime >5 years
 - High reliability
 - 3 year warranty
 - >300 installed
- Air-cooled
- Available in Cr, Cu, Mo, Ag



$I_{\mu}S$ & VÅNTEC-2000 vs. Classic Set-up Corundum Comparison

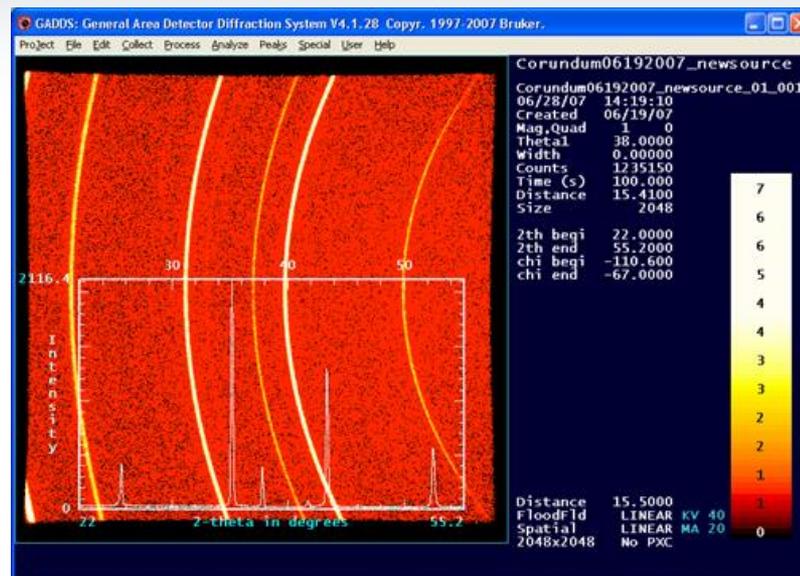
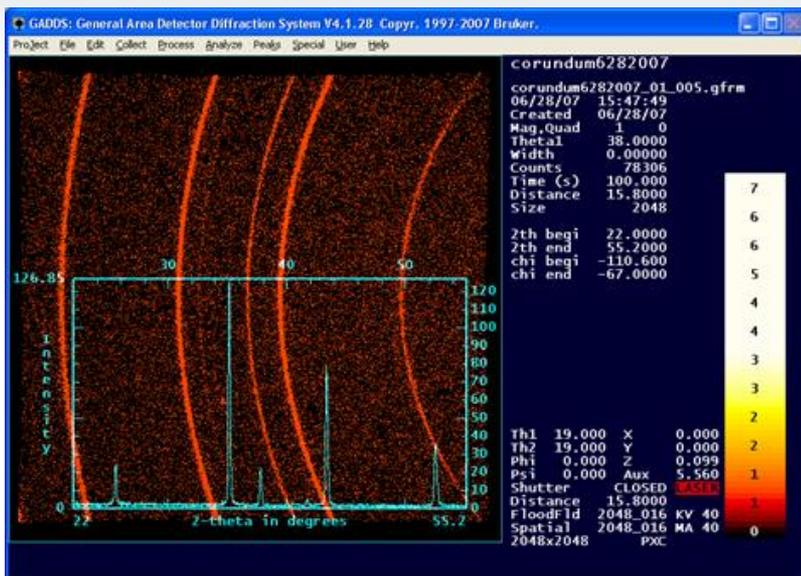


Sealed Tube with Göbel Mirror
45kV, 40mA, (1800 W)
0.3mm collimator
total counts: 78K


Microsource ($I_{\mu}S$)TM
45kV, 0.650mA, (30 W)
0.3mm collimator
total counts: 1235K

Intensity: 15.8x; Efficiency: 948x !


NEW: Liquid metal sources

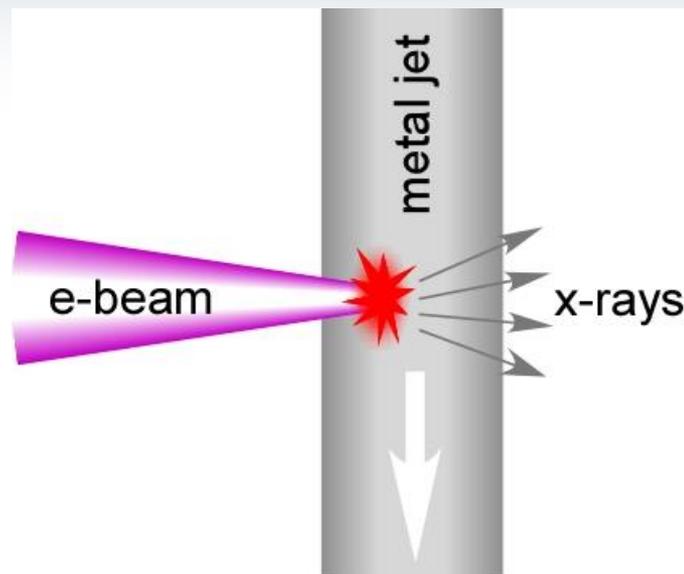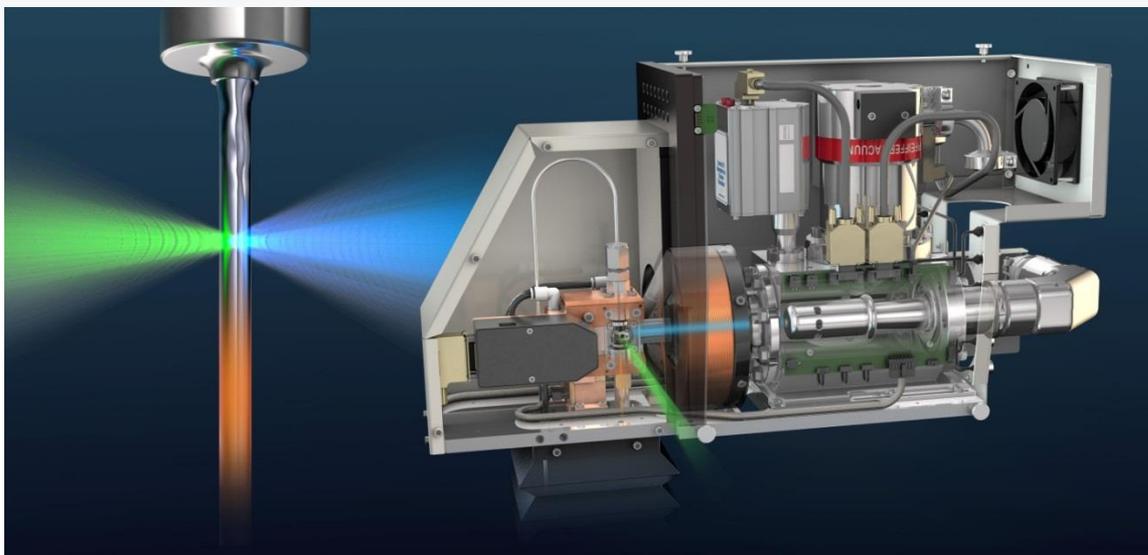

- High-speed liquid-metal-jet anode
- Anode is regenerative
- No longer limited by melting
- $>500 \text{ kW/mm}^2$ e-beam power density
 - Rotating anode limited to maximum 50 kW/mm^2

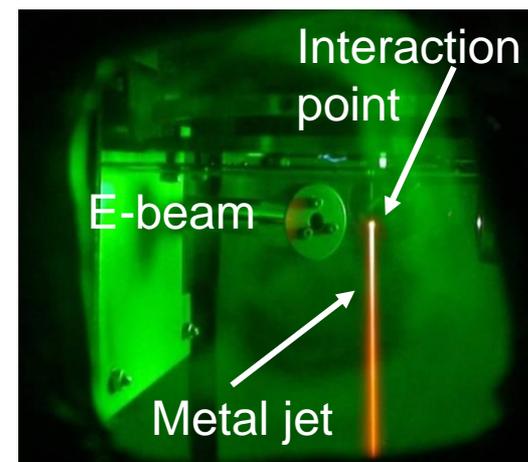

Source Spectrum and Brightness

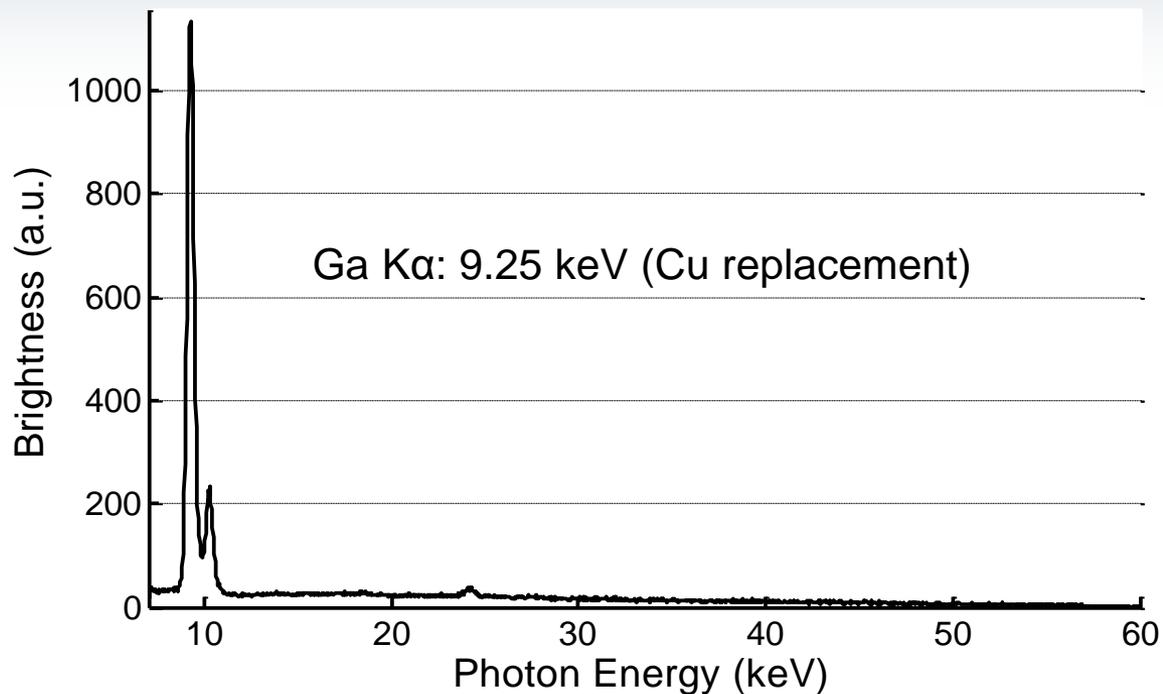


Spot size [μm , FWHM]	Voltage [kV]	Power [W]	Ga K α Brightness [Photons/(s \times mm ² \times mrad ² \times line)]
5	60	50	1.5×10^{11}
10	60	100	7.6×10^{10}
20	60	200	3.8×10^{10}

So, is it possible to put a synchrotron beamline on a table top?




- Yes, at least the equivalent of a typical present generation bending magnet beamline

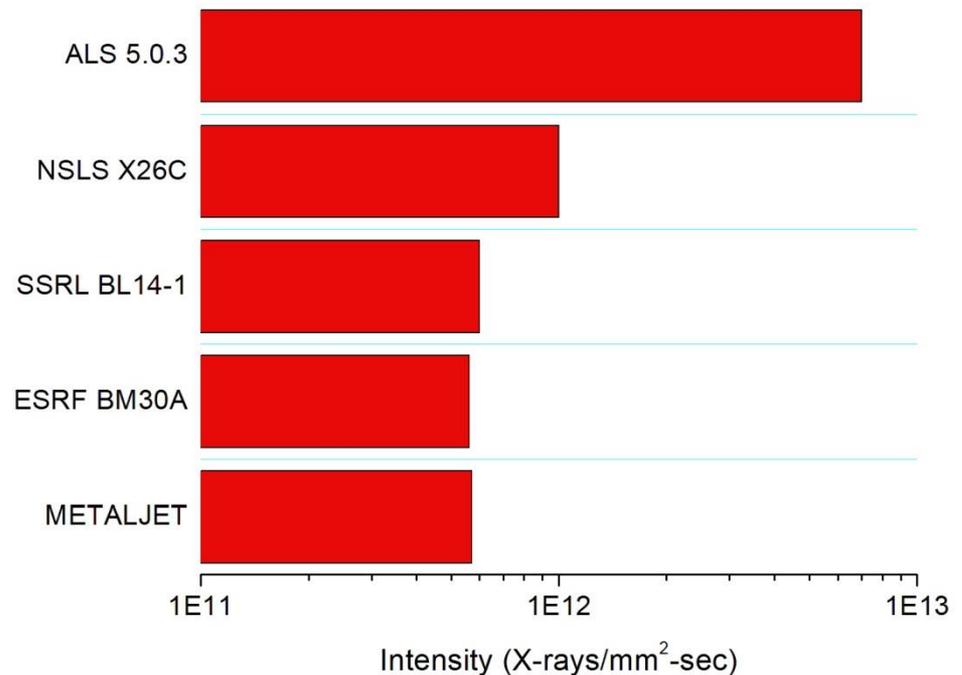

Detector

Choice of Detectors: 0D/1D/2D

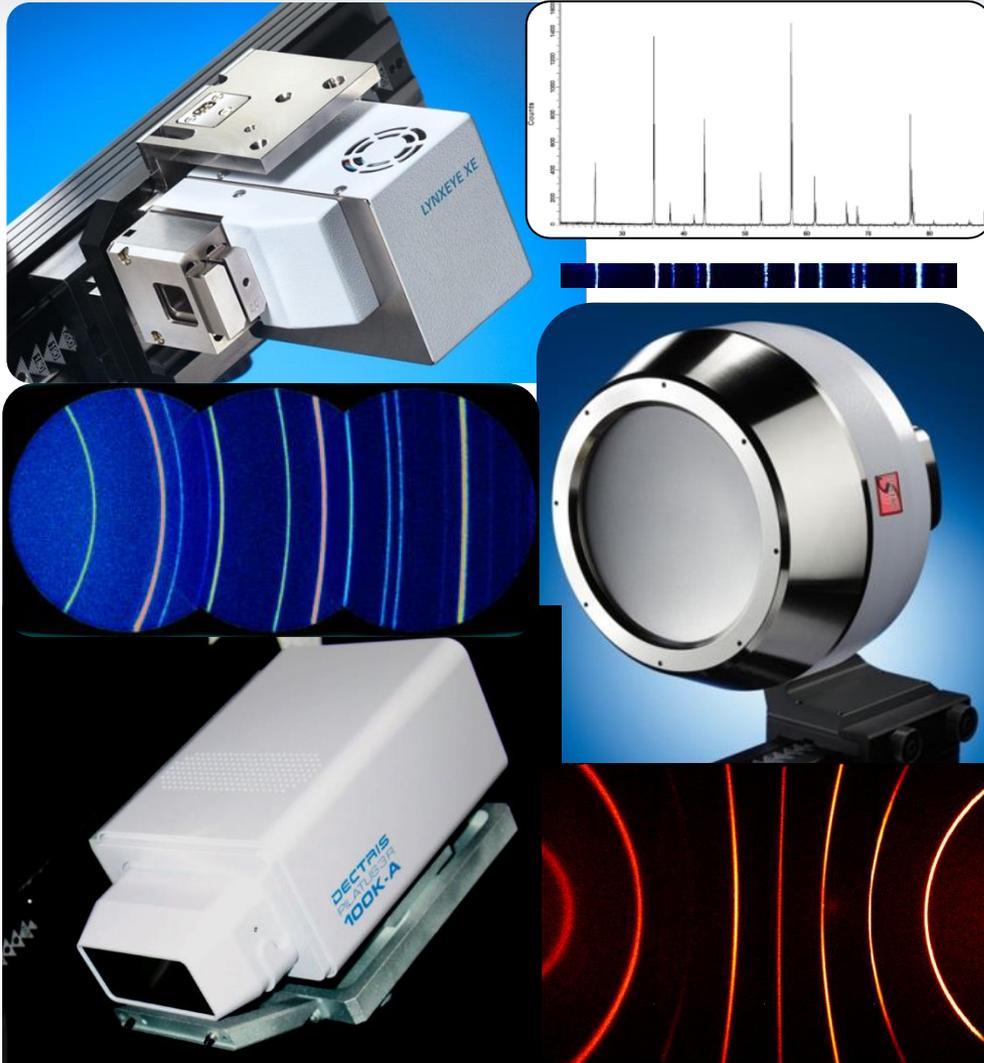

LYNXEYE XE

- Silicon strip technology
- Best Energy Discriminator of any 1D detector
- 0D, 1D and 2D modes
- Ideal detector for random powders and RSMs
- No Maintenance

VANTEC 500

- Detector with the Largest Active Area
- Radiation Hard
 - Can take the direct beam and strong reflections
- No Maintenance

PILATUS3 R 100K-A

- Hybrid Photon Counting (HPC) technology
- High count rate capability
- Sensitivity: Co, Cu + hard radiation (Mo, Ag)
- Active area: 83,8 x 33,5 (2.807) mm²No Maintenance
- Pixel size: 172 μ m (195 x 487 pixel)

VÅNTEC-500 – Outperforms all previous gaseous detectors.





- Large active area: 140 mm in dia.
- Frame size:
 - 2048 x 2048 pixels
 - 1024 x 1024 pixels
 - 512 x 512 pixels
- Pixel size:
 - 68 μm x 68 μm
 - 136 μm x 136 μm
 - 272 μm x 272 μm
- High sensitivity: 80% DQE for Cu
- High max linear count rate:
 - 0.9 Mcps – global;
 - 160 kcps/reflection -local
- Low background noise:
 - $<10^{-5}$ cps/pix
- Maintenance-free:
 - no re-gassing



D8 DISCOVER with PILATUS3 R 100K-A

2D HPC Technology for Your Lab Instrument

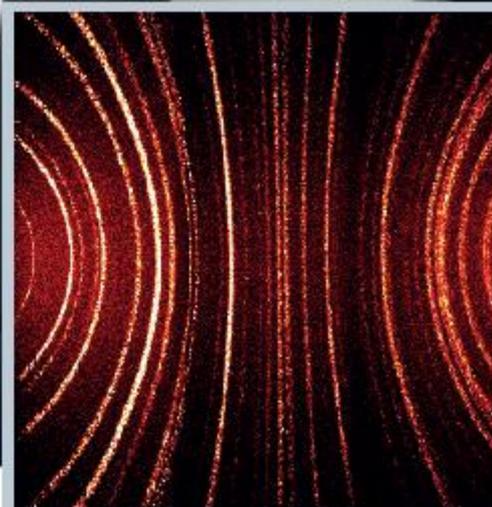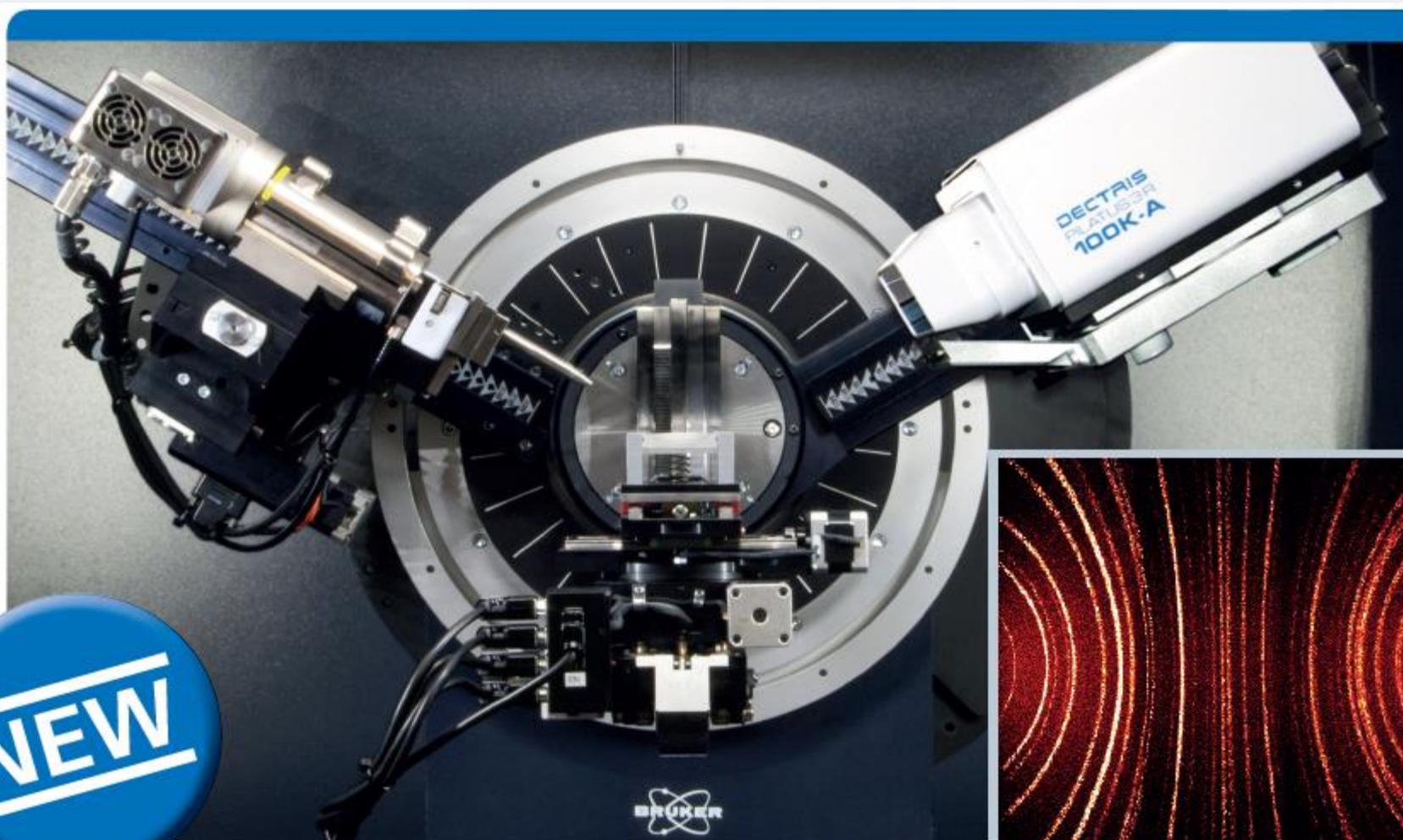

XRD² : Detector Orientation: γ -optimized vs. 2θ optimized





XRD² : Choice of Detectors: Active Area and Orientation



PILATUS3 R 100K-A

VANTEC-500

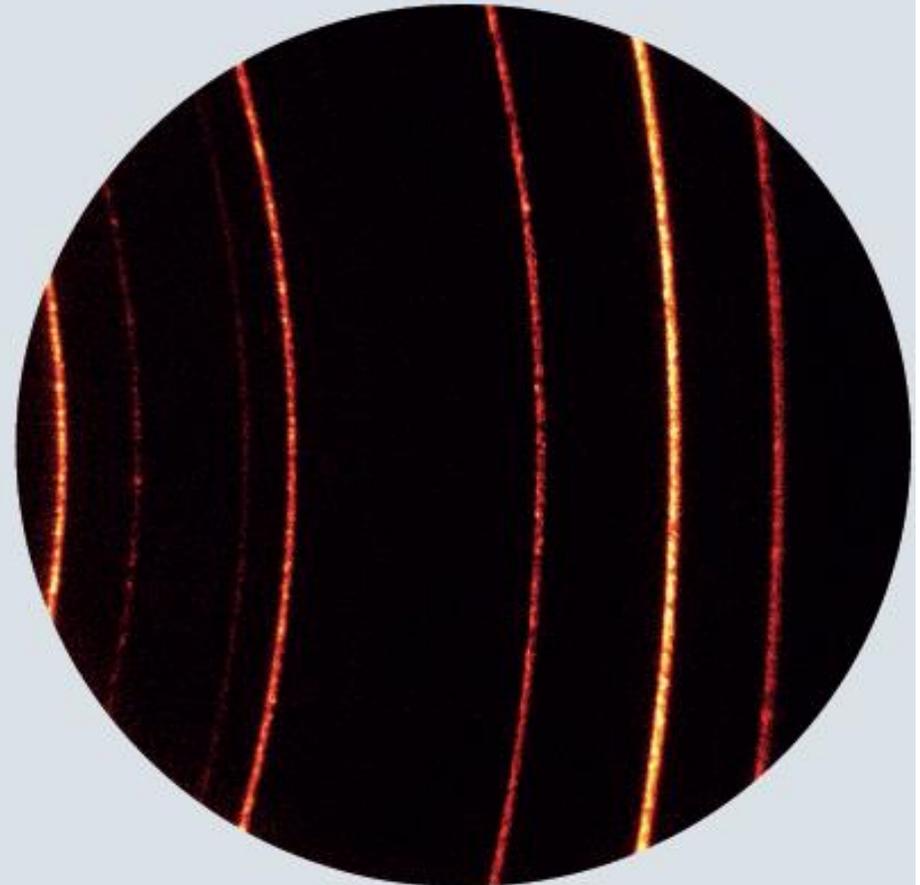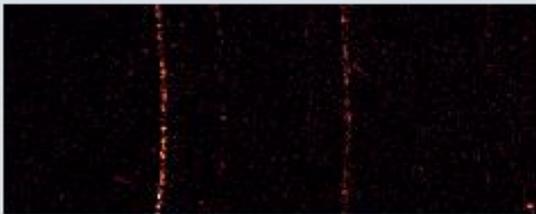

NIST 1976 (corundum) data collected at
sample-to-detector distance of 20 cm.

Left: PILATUS3 R 100K-A (2Theta-optimized Mode)


Middle: PILATUS3 R 100K-A (Gamma-optimized Mode)

Right: VANTEC-500

XRD² : Detector Orientation and Distance: frames collected with Aspirin




γ optimized – 0° Mode

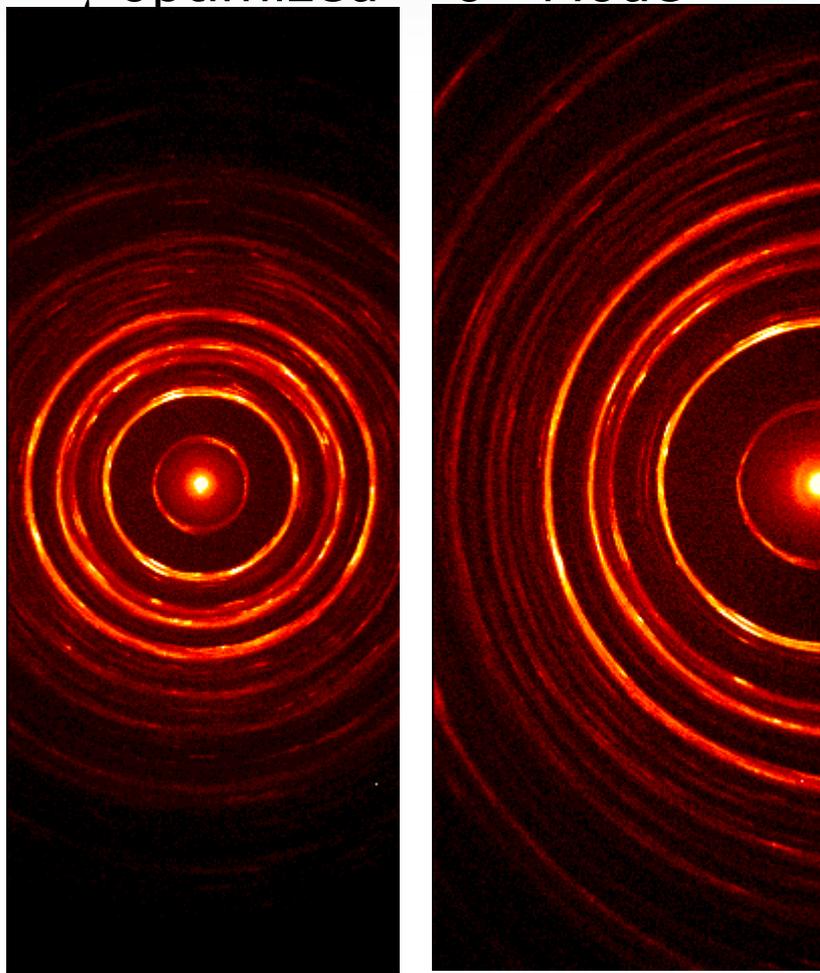


30 mm

50 mm & offset 19deg

2 θ optimized – 90° Mode

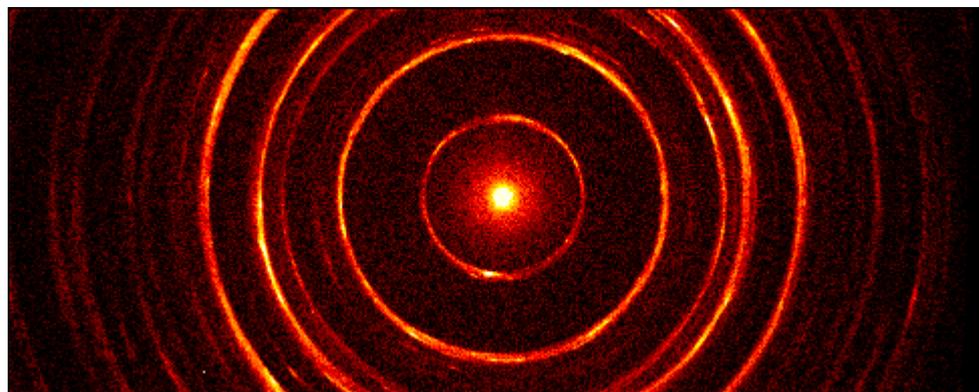

50 mm

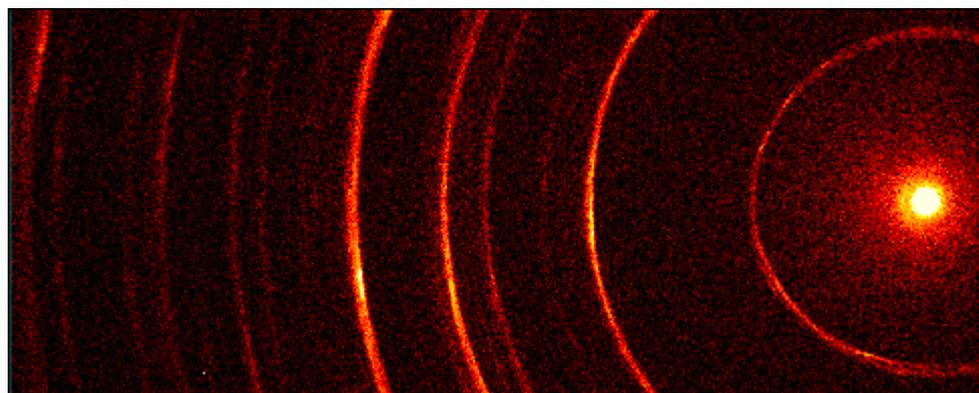

10 mm & 20 deg off

Sample Stage & System Configuration

XRD² : Sample Stages for Various Application



Centric Eulerian Cradle

- Most versatile stage on the market
- Ψ , ϕ , X, Y and Z are always mounted

Universal Motion Concept Stages


- Sits in front of the goniometer
 - Allows more weight and travel
- UMC 1516
 - Ψ , ϕ , X, Y and Z
- UMC 150 HTS
 - Designed for Reflection and Transmission
 - 96 well plates
 - X, Y and Z

Accessory Attachments

- Attach directly to the XY table
- Standard Powder Adapter (included)
- Dome Temperature stages
- Capillary Attachment
- Wafer Chucks

Horizontal System

Horizontal $th-2th$, CEC

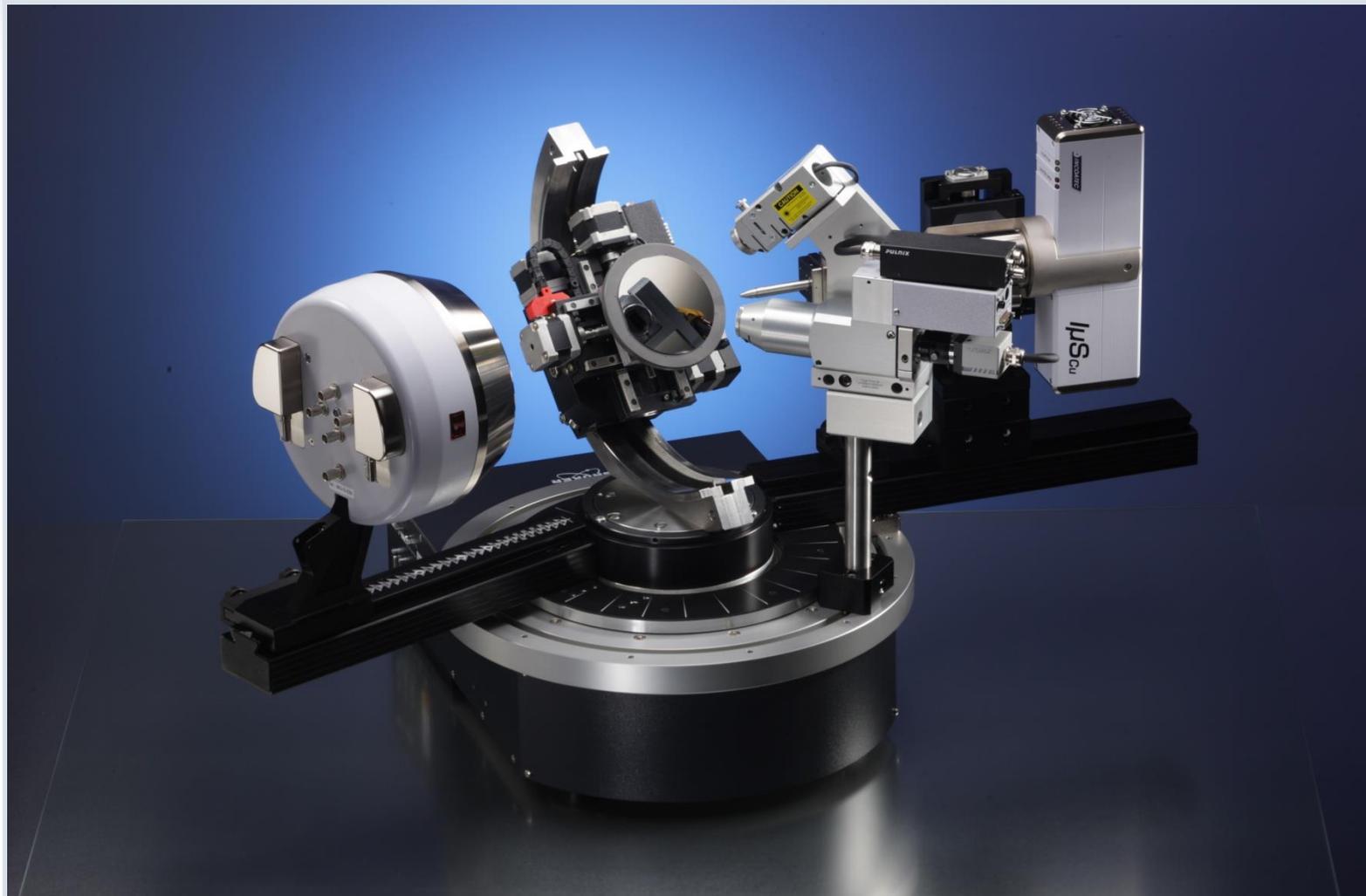

No barrier between 0D/1D/2D

Vertical theta-theta, CEC for microdiffraction/stress/texture




High-throughput Screening (HTS)

Vertical theta-theta, Reflection/Transmission




D8 DISCOVER High Throughput Screening

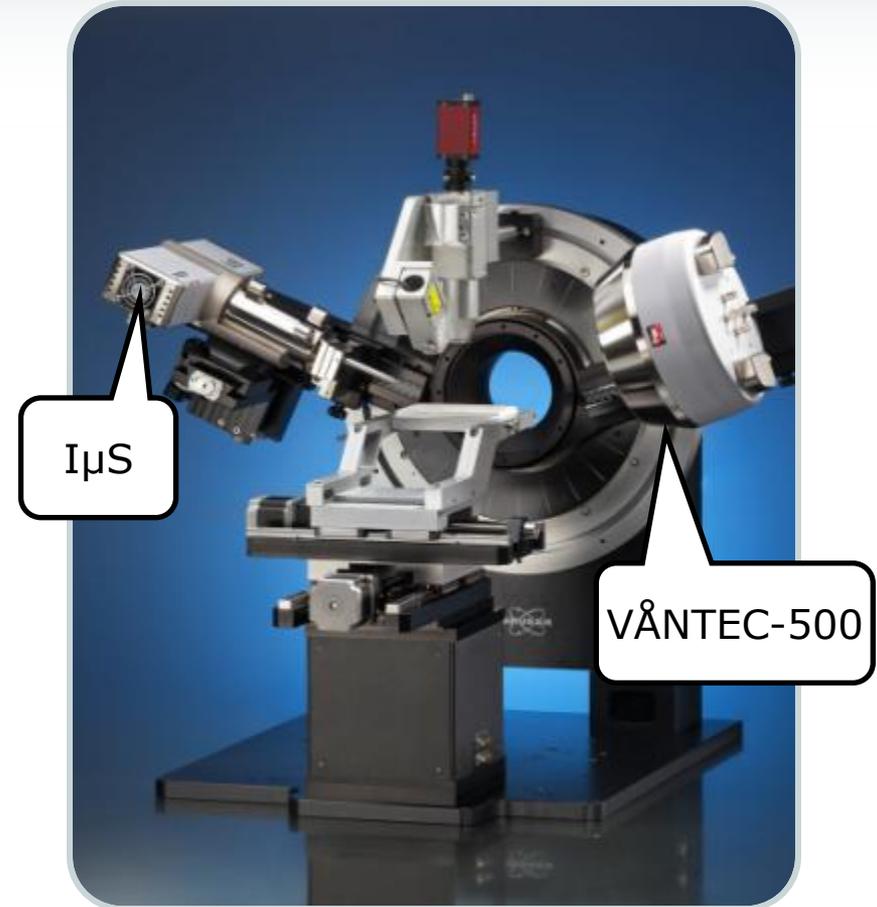

Phase Identification

XRD²: Single Frame Covering All

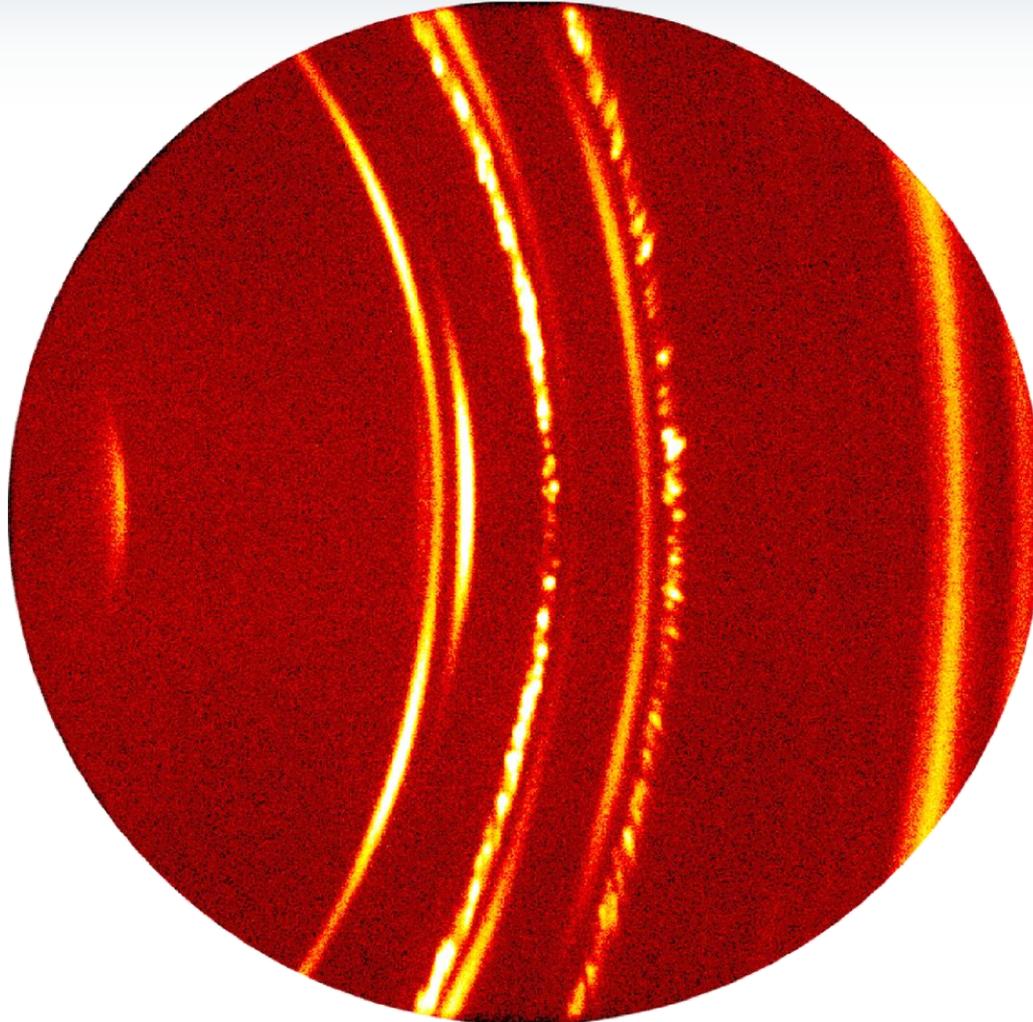



- Multilayer battery anode.
- 2θ coverage: 70° at 8 cm detector distance
- A single frame showing information on phase, stress, texture and grain size
- 2D detector is essential for In-situ measurement

Ibuprofen Integrated to 1-D data

t	Dist	Beam
15 s	15cm	300 μm



Ibuprofen - 15s

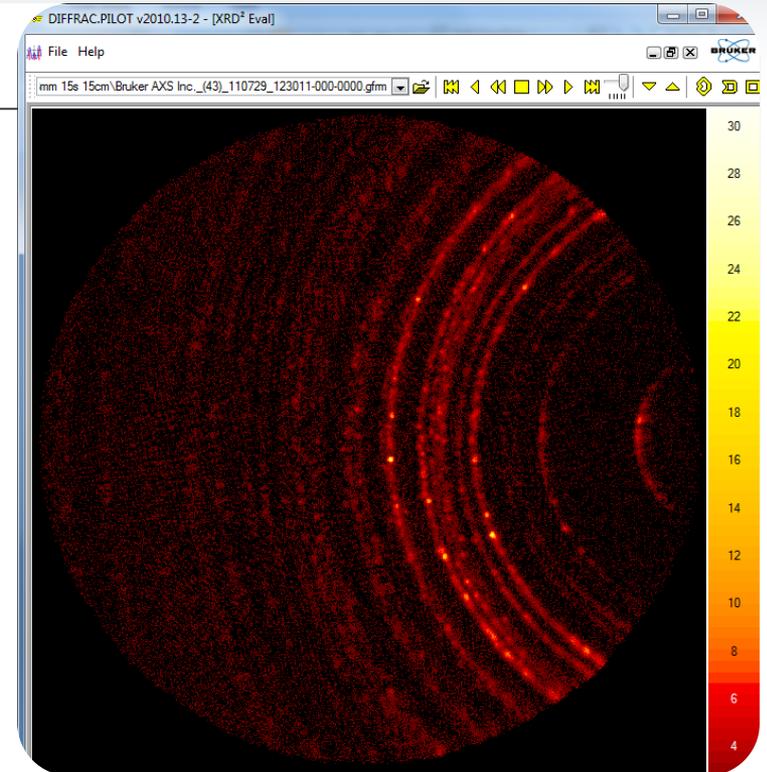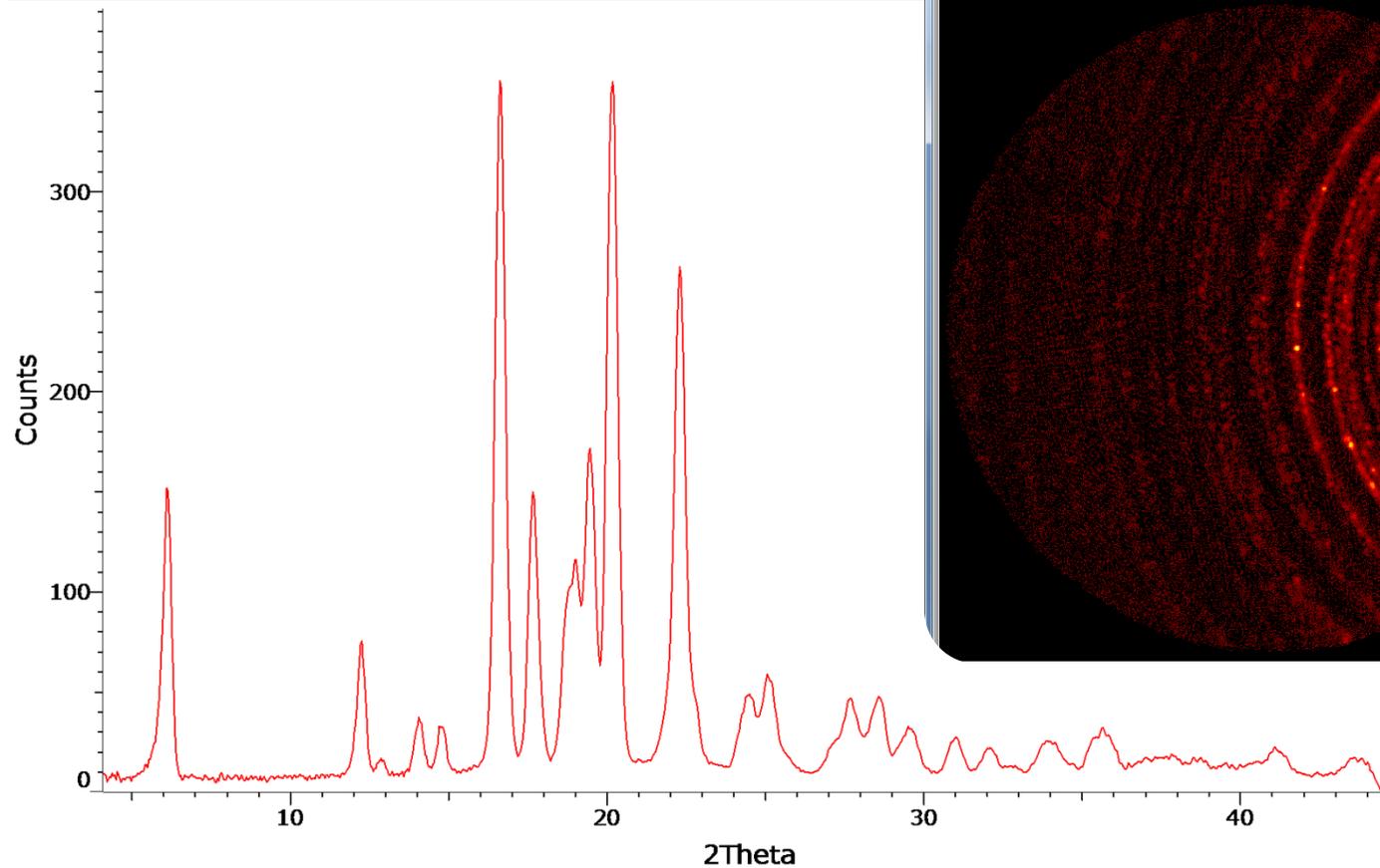




XRD²: Data Collection:




Acetaminophen powder

5 second data collection

30 second data collection

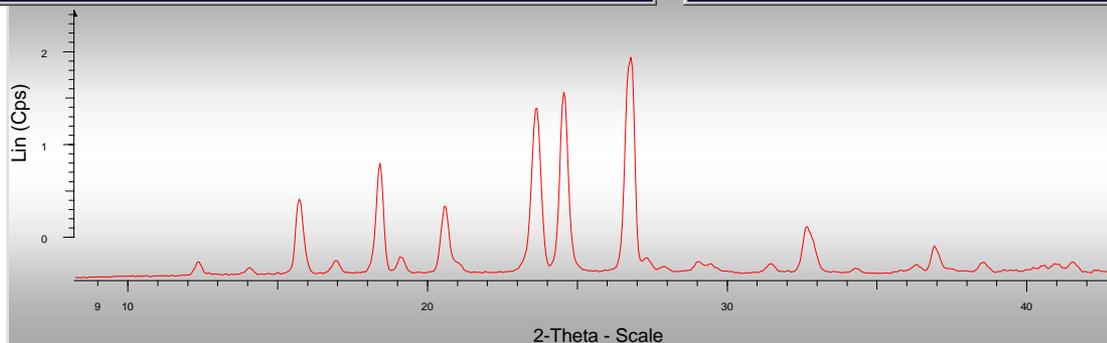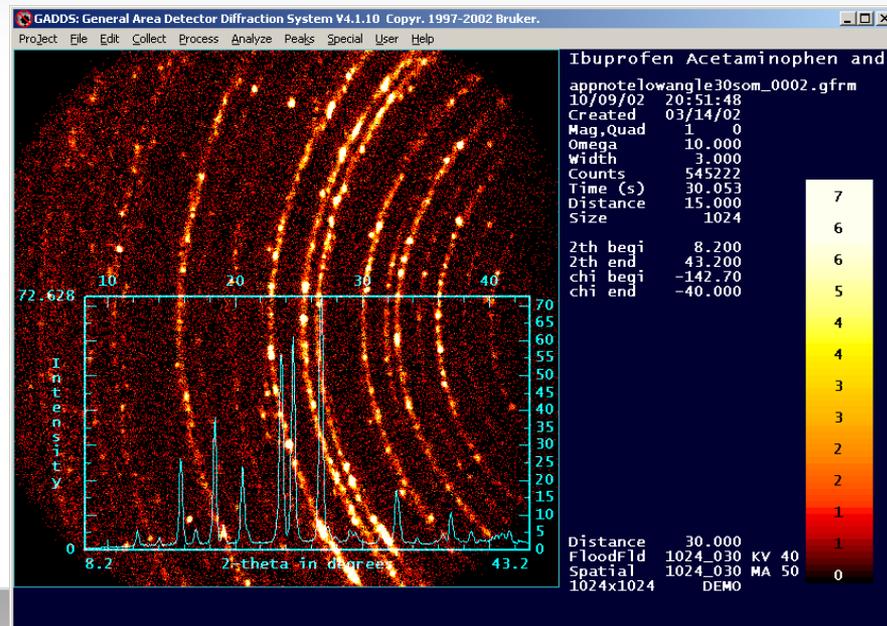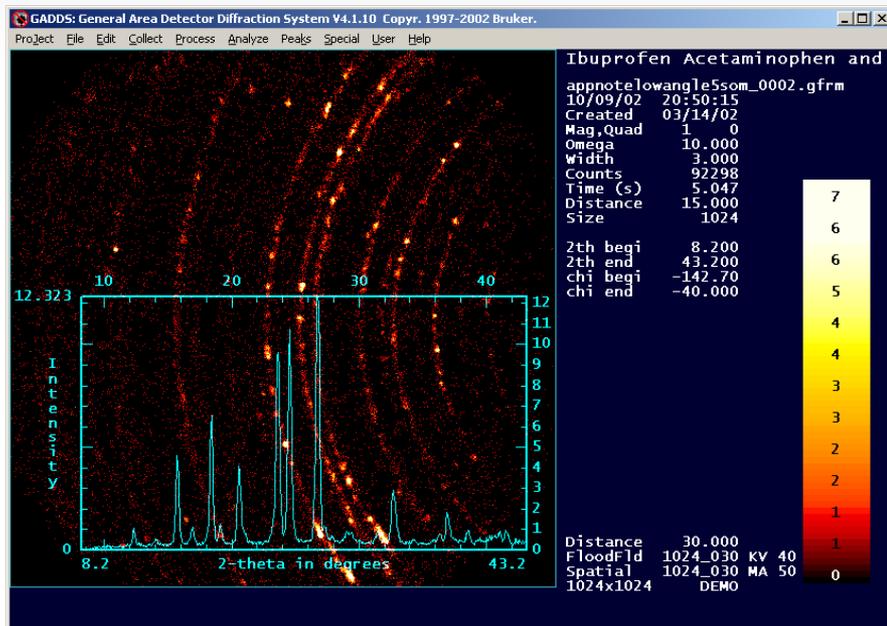


XRD²: Frame Merge and Integration

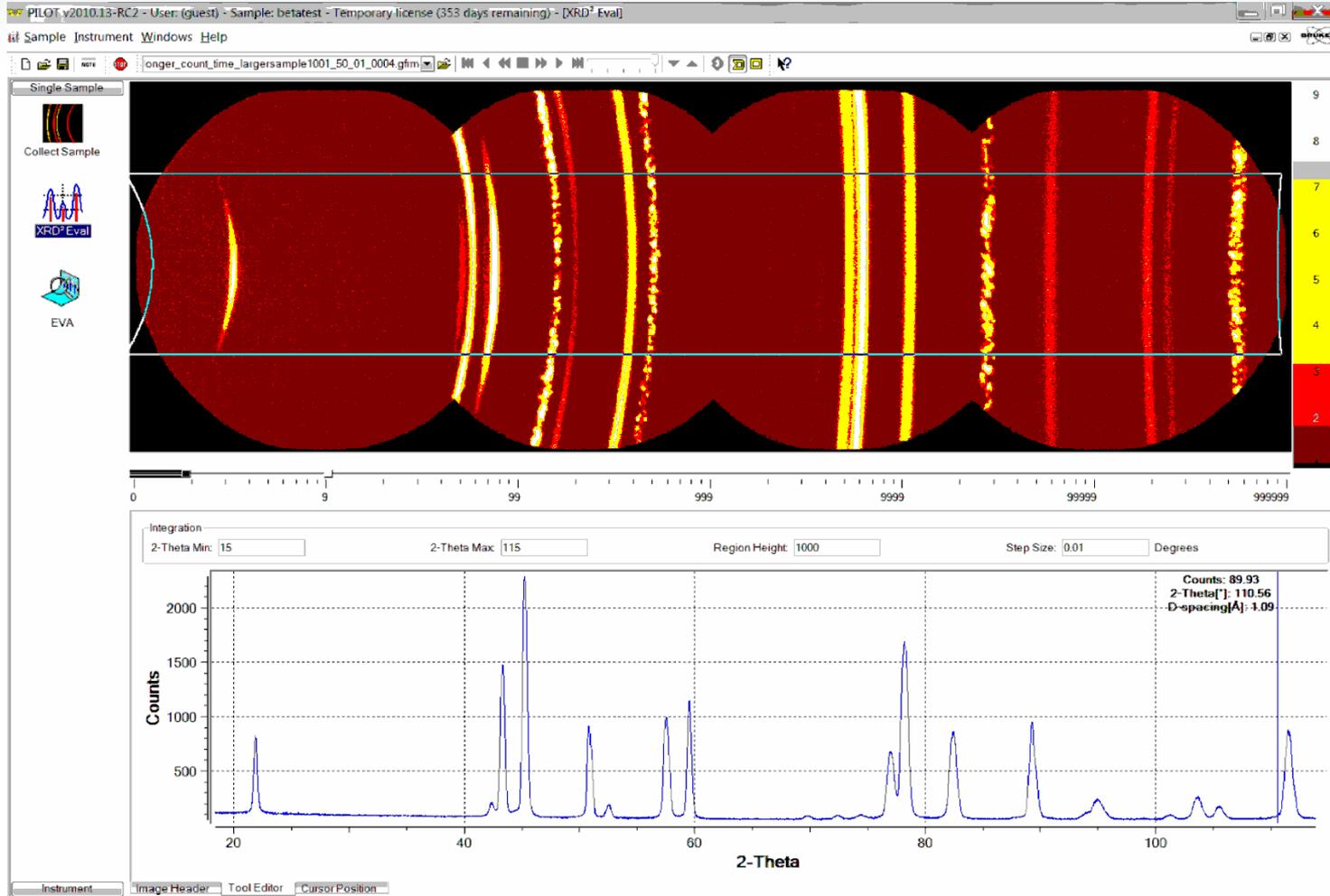

- 4 frames at 20 cm
- Merged frames 2 θ coverage: 100°
- Integrated profile for phase ID search/match

XRD²: Mapping: API Distribution in a Pill


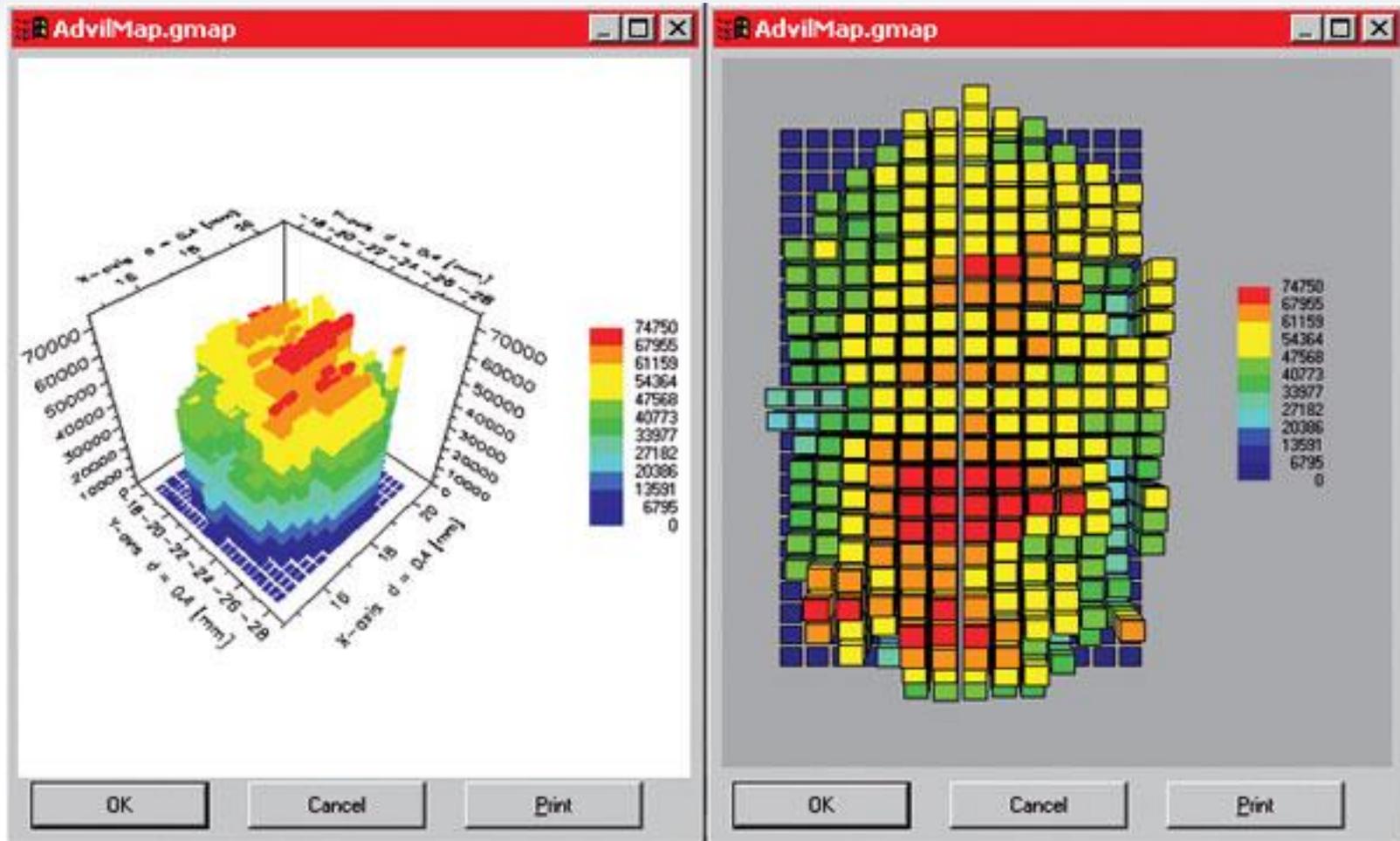


XRD²: High throughput screening

Laser/video sample alignment



Easy and accurate sample positioning without touching the sample surface

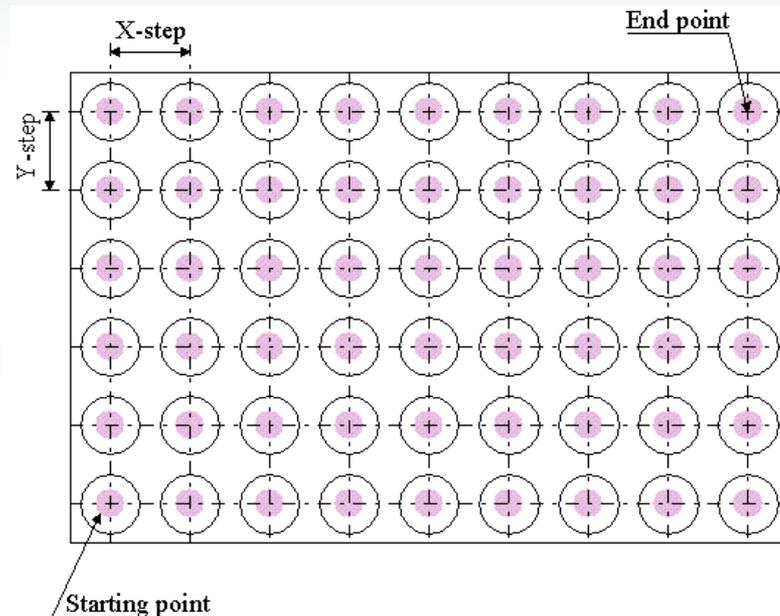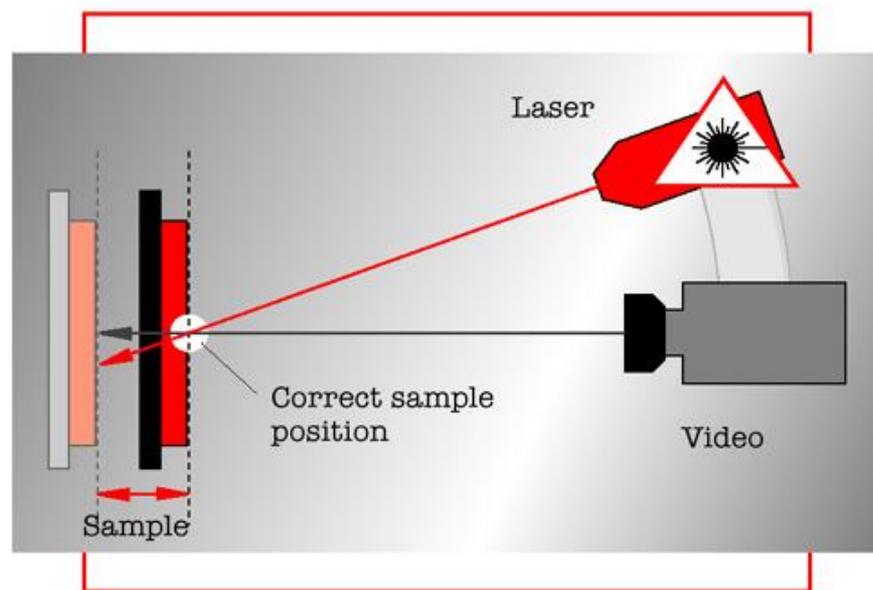

Video image of each material library spot can be automatically stored during data scan

Particle/Crystal Size

XRD²: Crystal Size by γ profile analysis: Organic glass for food & drugs




Food

Bio-preservation



In sugar glasses


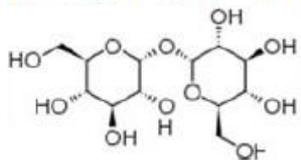


In amber





Annealing of GC crystals

LM

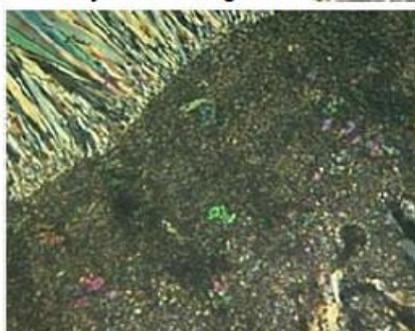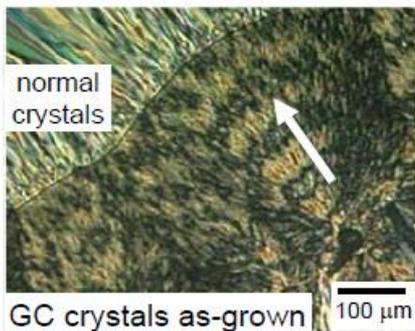


50 min/313 K

AFM

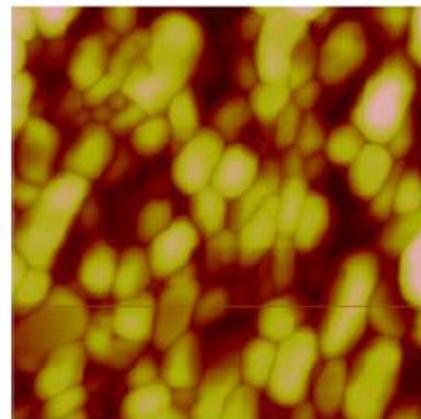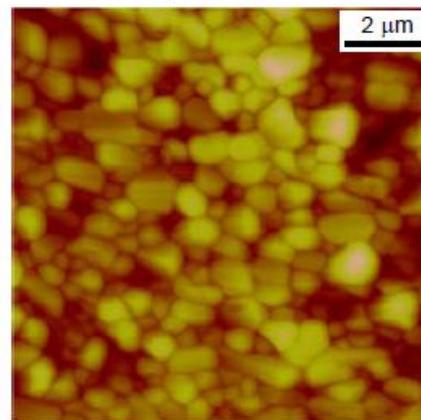


30 min/313 K


2D XRD

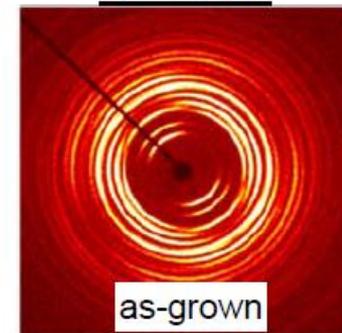

as-grown

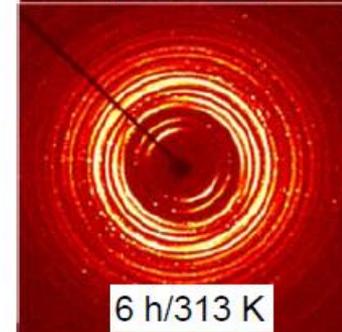

6 h/313 K

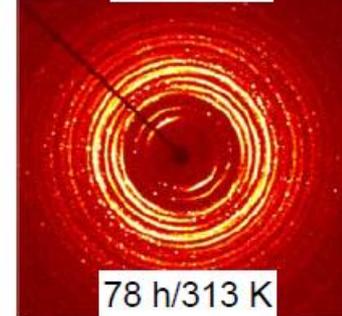


78 h/313 K


*Courtesy of Prof. Lian Yu, U. of Wisconsin -
Madison

XRD²: Data Collection:

Acetaminophen powder

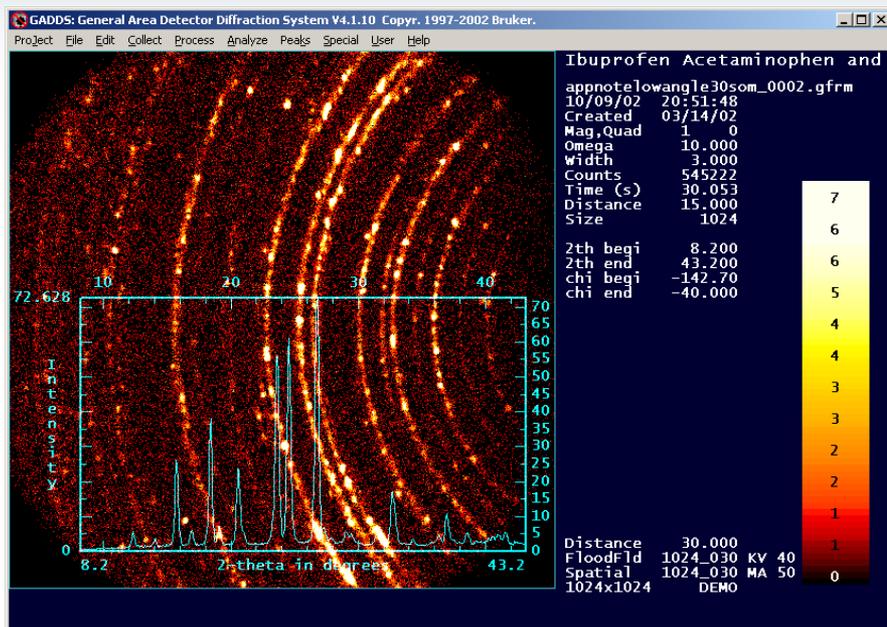

- The spotty diffraction ring is due to the large crystallites compared to the sampling volume (beam size).
- The number of spots on the ring is determined by crystallite size, instrumental window (γ -range), multiplicity of the crystal plane, and effective diffraction volume.
- The size of jelly beans and candy bin determines how many you can fill.

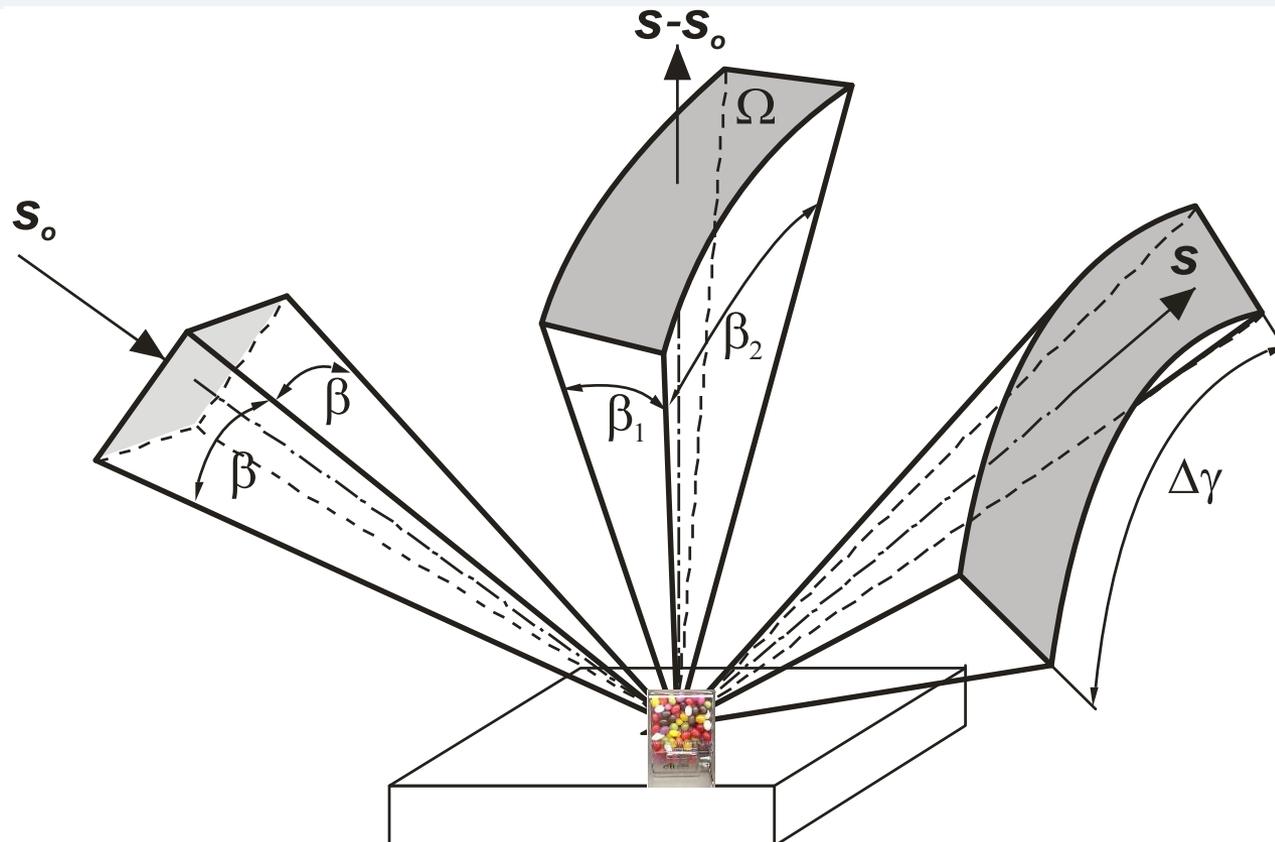

For XRD², the instrumental window Ω is given by

$$\Omega = \beta_1 \beta_2 = 2\beta \arcsin[\cos \theta \sin(\Delta\gamma / 2)]$$

XRD²: Particle size measurement by γ profile analysis: **BRUKER**



For XRD² in reflection mode, the crystallite size is given by

$$d = k \left\{ \frac{p_{hkl} b^2 \arcsin[\cos \theta \sin(\Delta\gamma / 2)]}{2\mu N_s} \right\}^{1/3}$$

where μ is the linear absorption coefficient

For transmission mode with the incident beam perpendicular to the sample surface, the crystallite size is given by

$$d = k \left\{ \frac{p_{hkl} b^2 t \arcsin[\cos \theta \sin(\Delta\gamma / 2)]}{N_s} \right\}^{1/3}$$

where t is the sample thickness.

k is the instrument calibration factor or can be calculated from:


$$k = \left(\frac{3\beta}{4\pi} \right)^{1/3} \text{ if the instrument broadening in } 2\theta \text{ direction is known.}$$

XRD²: Particle Size Analysis





2θ profile analysis


γ profile analysis

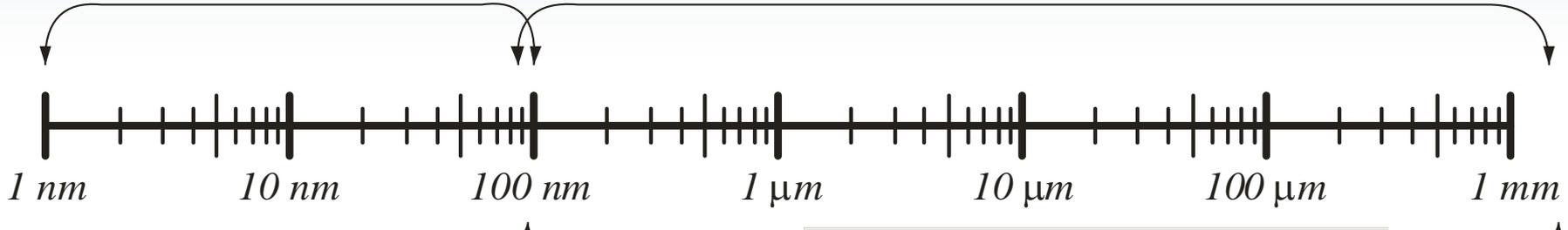

Scherrer equation:

$$t = \frac{C\lambda}{B \cos \theta}$$

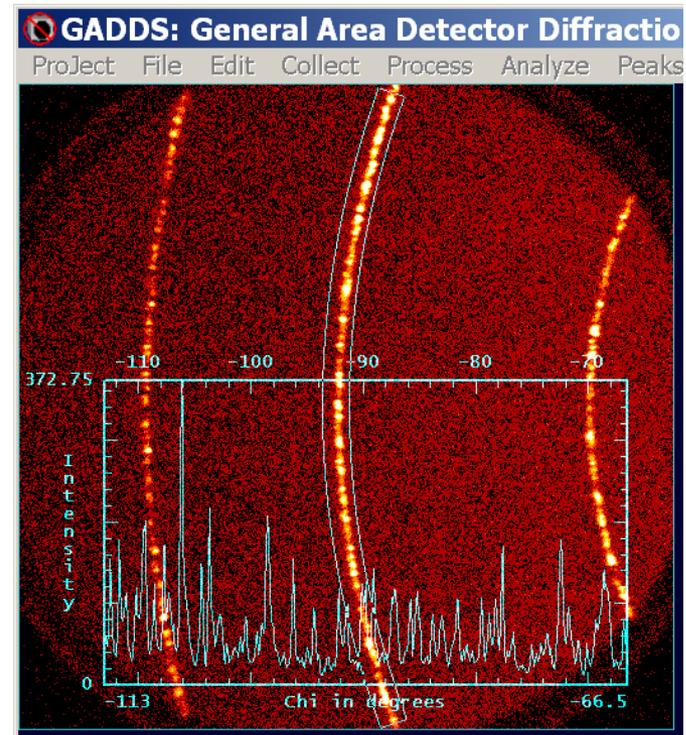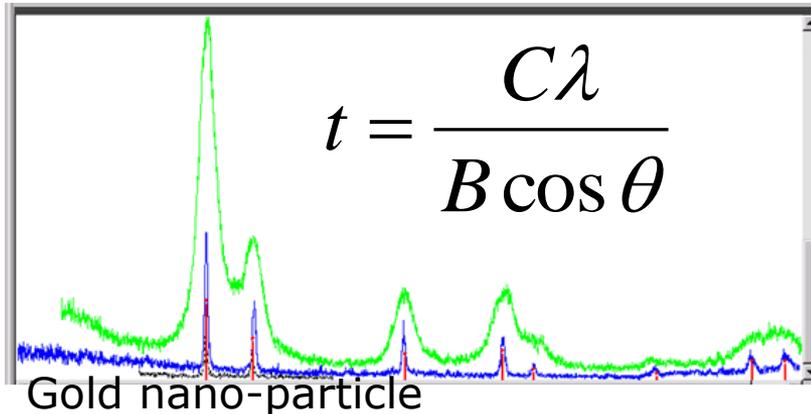

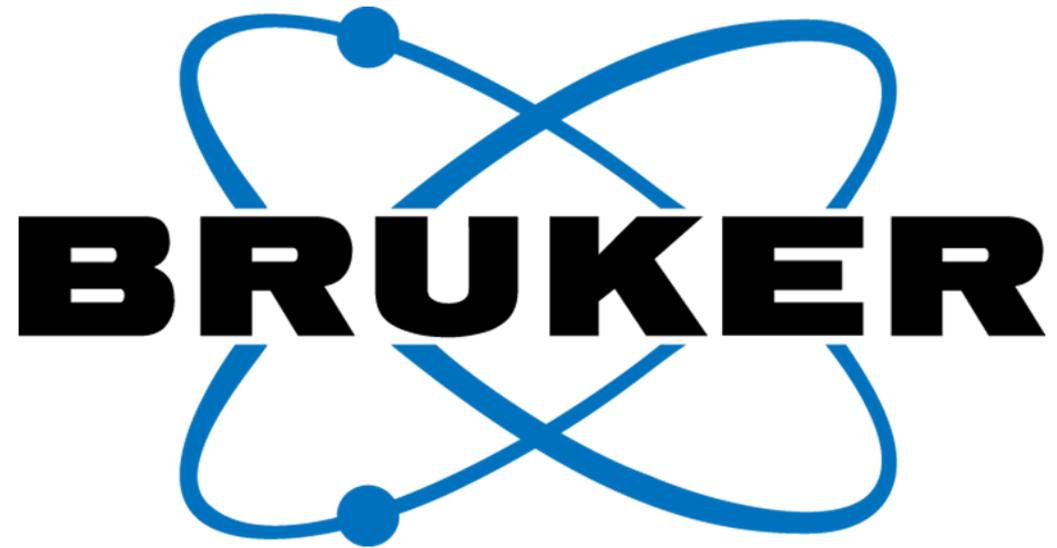

www.bruker.com